

## Product Release Notes

**Product:** PONEMAH Physiology Platform  
**Model:** P3 Plus  
**Version:** 4.70-SP1 (Service Pack 1 for version 4.70)  
**Build:** J03263 (CD Build)  
**Date:** October, 2008

Product Release Notes for PONEMAH Physiology Platform (P3 Plus) Version 4.70-SP1, indicate revisions made to the P3 Plus core application since release of P3 Plus Version 4.70. For information regarding changes to the software from previous versions, please refer to the Release Notes folder located on the Version 4.70 CD. Product Release Notes indicate only revisions to application contents that are part of CD Part #J02910 – Build Version 4.70.

Product Release Notes for P3 Plus do not include information regarding revisions of the P3 Plus Analysis Modules. Information regarding revisions of the P3 Plus Analysis Modules is indicated in separate Product Release Notes.

**Notice for organizations that must comply with FDA's Good Laboratory Practices (GLP) and 21 CFR Part 11 Electronic Records; Electronic Signatures:** P3 Plus Versions may contain **Preview Features**. These **Preview Features** are listed in the Product Release Notes table under the column, "Type of Change". A **Preview Feature** indicates that enhancements have been made to P3 Plus, but have not been validated. Instead, Data Sciences International (DSI) has opted to delay complete validation until receiving comments from customers regarding use of these features. Further validation of these features will be performed in later releases of P3 Plus. There may be additional **Preview Features** that had been documented in previously released versions that are not documented here. These features are not available unless manually enabled by the user. If documentation is needed regarding these features, please contact the Technical Support Group at DSI.

Reference #	Type of Change	Key: N = New Feature; E = Enhancement; F = Fix
		Description
VivoMetrics		
N/A	E	<p>Added the ability to graphically view the three individual accelerometer channels (x, y, and z) collected by the Vivometrics hardware.</p> <p>Note: when opening a previously configured protocol using the VivoMetrics interface, a message will be displayed informing the user that additional channels have been added. The protocol will need to be modified to account for the three new channels.</p>

**DSI Corporate Headquarters**

119 14<sup>th</sup> Street NW  
St. Paul, MN 55112  
P: +1 (651) 481.7400 • F: +1 (651) 481.7404  
[www.datasci.com](http://www.datasci.com)

**Ponemah Operations**

5525 Cloverleaf Parkway  
Valley View, OH 44125  
P: +1 (216) 328.7000 • F: +1 (216) 328.7400  
[www.ponemah.com](http://www.ponemah.com)