

INSTALLATION NOTES

MX2 FIRMWARE 1.0.11

Prior to installing, it is recommended to review the MX2 Firmware 1.0.11 Release Notes. The following are the instruction to install the firmware update.


ACCESSING MX2 FIRMWARE

The website is the primary means for delivering firmware. If a CD is desired, one can be requested by contacting DSI Customer Service (CService@datasci.com).


1. Navigate to www.datasci.com
2. Browse to **Resources | DSI Gateway Portal**.
3. Login using your **User Name** and **Password**. If you have forgotten or do not have a User Name and Password, the ability to request a new password or register at the site is available.
4. Select **Software Updates**.
5. Click the **MX2 1.0.11.zip** to download the contents to a desired location.
6. Extract the files by **right-clicking** the .zip file and choosing **Extract**.

Installing MX2 Firmware


1. Launch **Ponemah**.
2. Select the **Hardware | MX2 Diagnostics...**
*Note: If using Ponemah v5.20, select **Hardware | Diagnostics...***

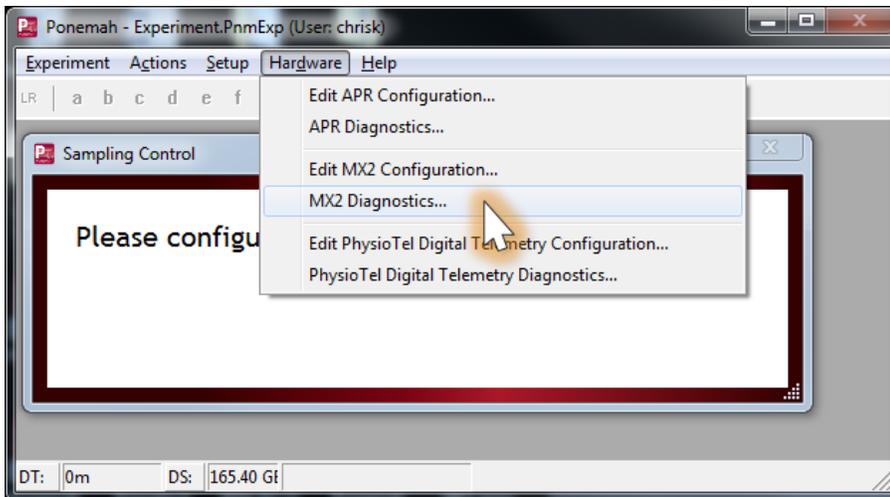

- Choose the MX2 from the *Name – Model Name (Serial Number)* dropdown.
Note: Select the Refresh List button if the MX2 does not appear in the dropdown list.

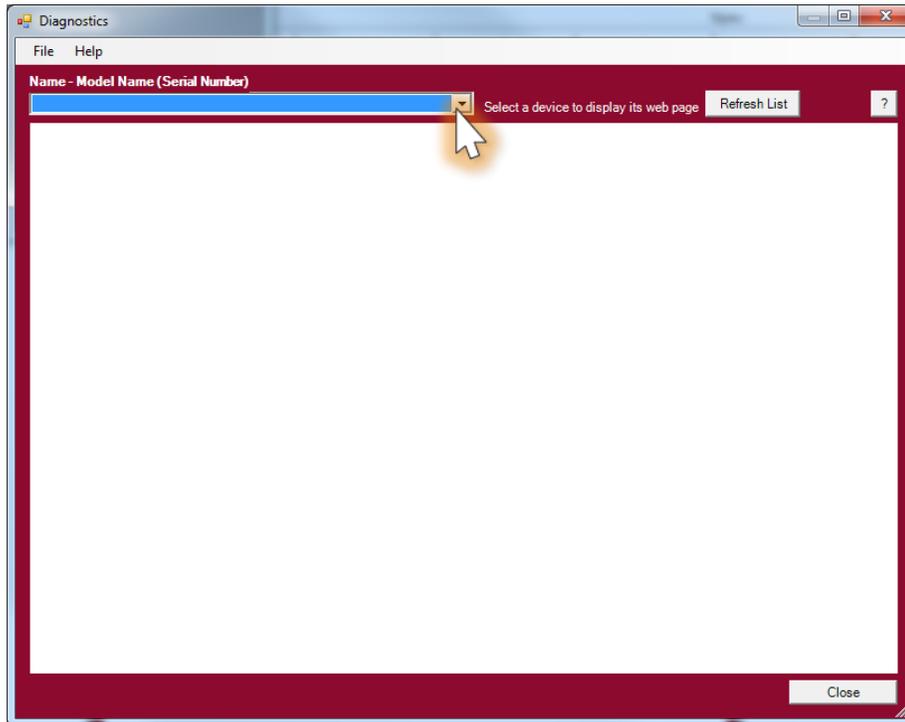


- Select the *Upload MX2 Firmware* page from the list view on the left side of the diagnostic page.

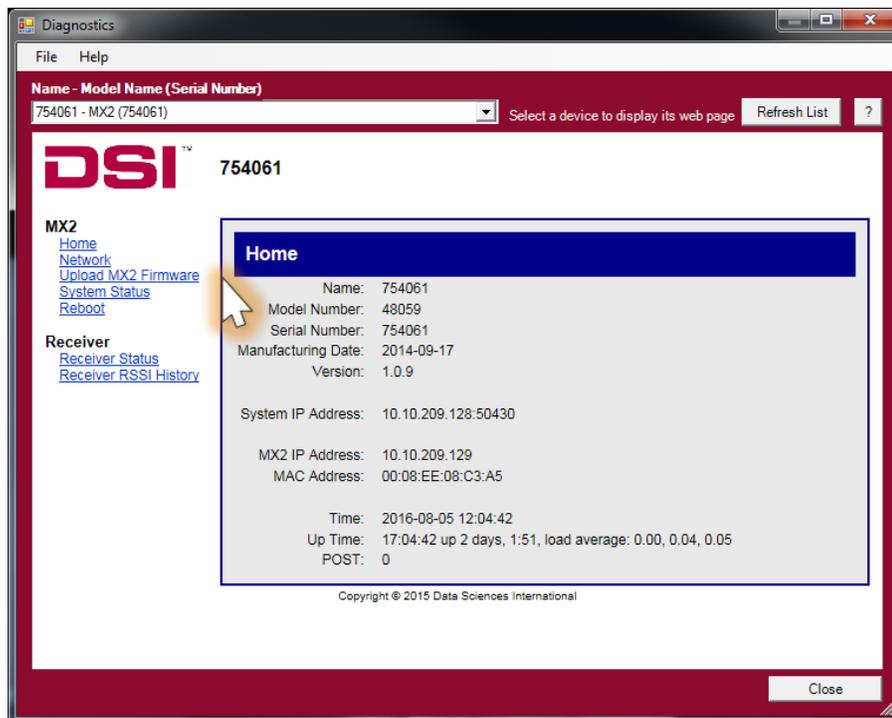


5. Please complete steps a. – d. below twice.


- a. Select the **Browse...** button.

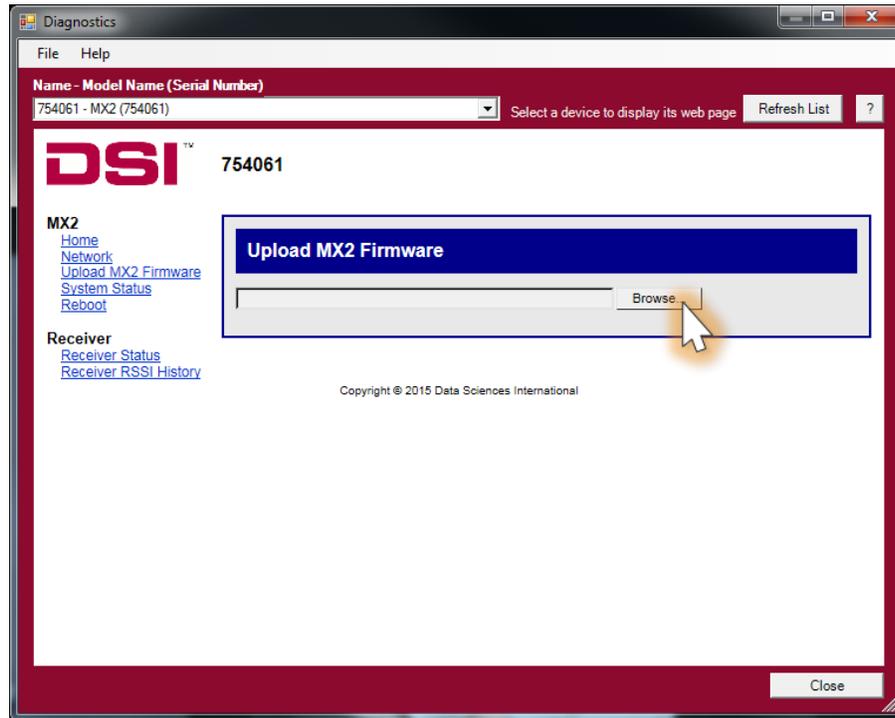

- b. Select the *MX2 1.0.11* file obtained during the .zip extract step above. Once uploaded, the following message will appear: *Validated. Upgrade will be applied during reboot.*

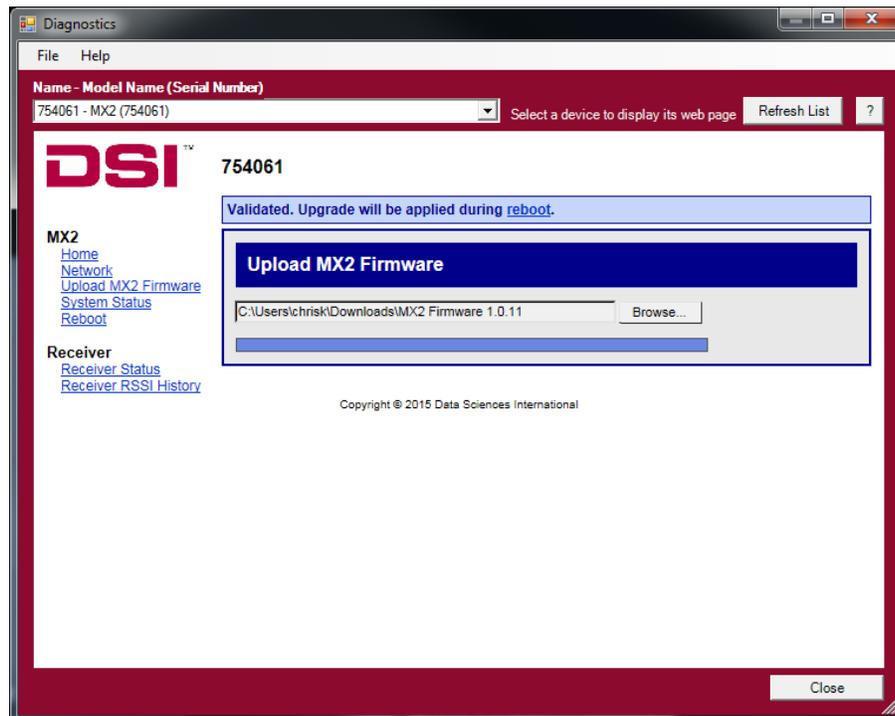

- c. Select the **reboot** link within the message to navigate to the *Reboot* diagnostics page.

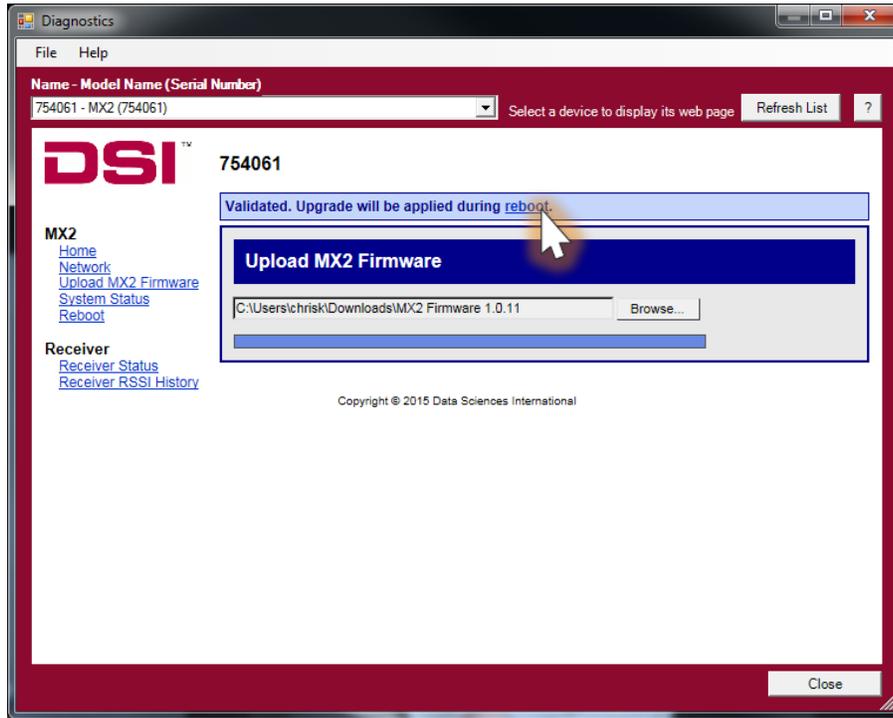

- d. Click the **Reboot** button.

