

# FINDING THE IP ADDRESS OF A COMMUNICATION LINK CONTROLLER (CLC) OR MX2

There may be times when it is helpful to know the IP Address of the CLC or MX2. There are two easy ways to locate the IP Address of the CLC or MX2. The instructions below outline the 2 methods for locating the IP Address.

## PONEMAH DIAGNOSTICS

1. With the CLC connected to the system, launch **Ponemah**.
2. Select **Hardware | PhysioTel Digital Diagnostics...**

If using the MX2, select **MX2 Diagnostics...**

*Note:* If using Ponemah v5.20, select Hardware | Diagnostics...In the main Ponemah Window, go to Hardware | Diagnostics.

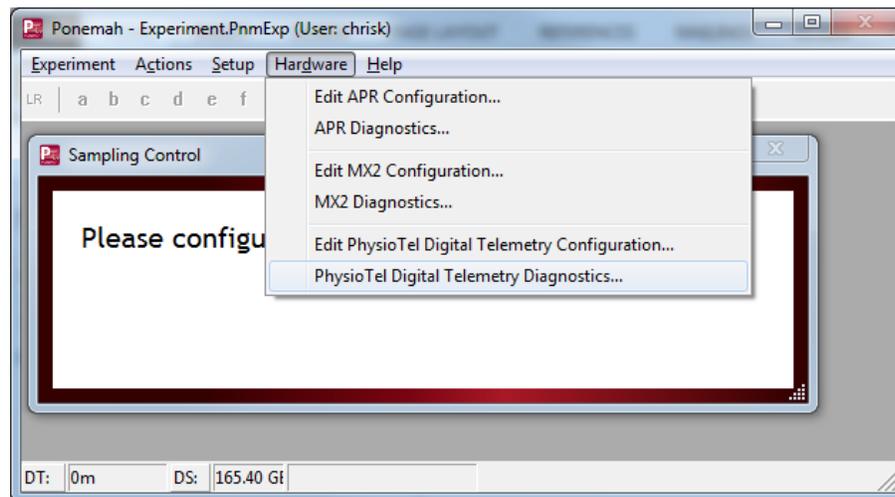


Figure 1 - Diagnostics Pulldown

3. Select the CLC you would like the IP address for from the dropdown list.
4. Located the **CLC IP Address** for your CLC's IP Address. See **Figure 2**.  
If using the MX2, locate the **MX2 IP Address** within the *Home* page.

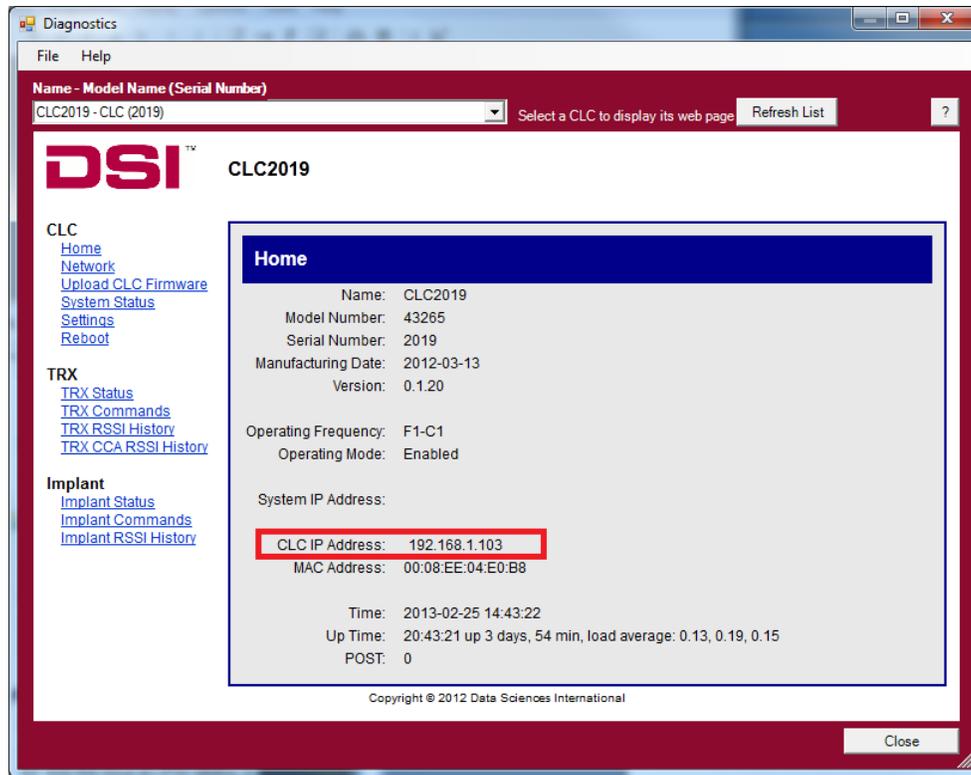


Figure 2 - CLC IP Address on Diagnostic Page

## COMPUTER NETWORK LOCATIONS

1. In Windows 7, click on the **Windows Orb**  (the Start button) and go to **Computer**.
2. On the left-hand menu pane of Computer, scroll down and select **'Network'** from the list.
3. In the right-hand pane, under **Other Devices**, the CLCs and MX2s connected to the network can be seen.
4. **Right-click** on the CLC or MX2 you would like the IP address for and select **Properties**. See **Figure 3** for an example.

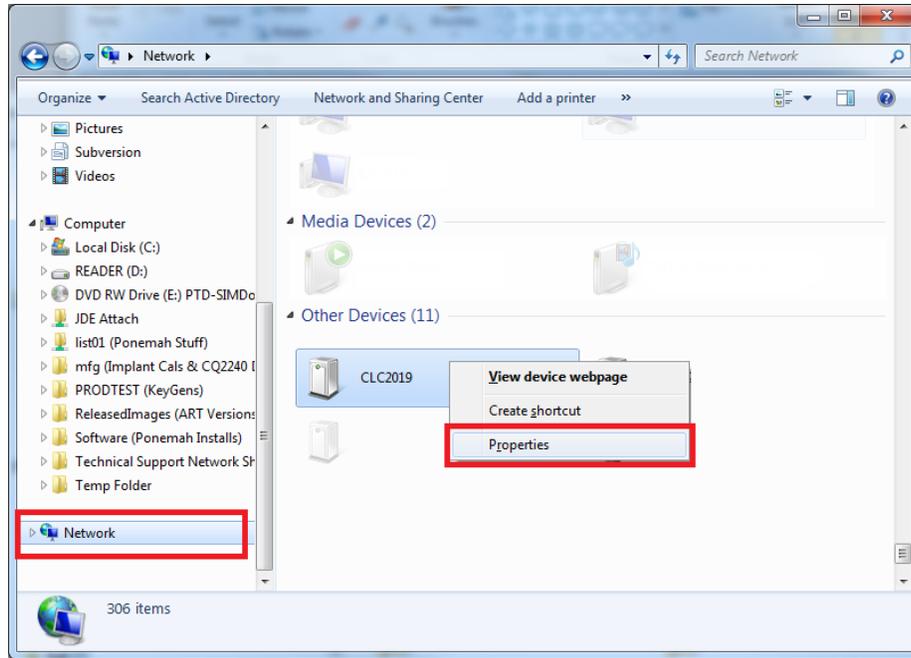


Figure 3 - Network Locations and Properties

5. Under **Properties**, look at the bottom of the window next to **IP Address**, this is the IP Address of your CLC or MX2.

*Note:* It may take up to 2 minutes for Network to display all the discovered connected devices.

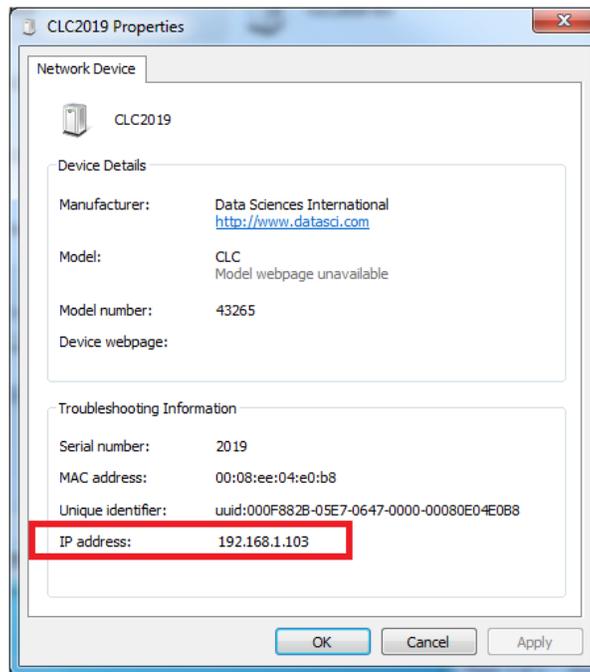


Figure 4 - CLC IP Address in Properties

