

# PHYSIOTEL DIGITAL RF ASSESSMENT PROCEDURE USING PONEMAH


The RSSI Threshold for the PhysioTel™ Digital Implants and Transceivers (TRX) can be adjusted in the Communication Link Controller (CLC) Diagnostic page in Ponemah. Changes to the RSSI Threshold should only be made if advised by DSI. The guide below explains how to perform a basic RF Assessment to determine the noise level in the room. This information should be provided to DSI Technical Support.

## PERFORMING AN RF ASSESSMENT


Make sure the CLC is set to the frequency you want to scan and that TRXs are connected. Turn off any implants that are set to the same frequency you are checking.

1. With the CLC connected to the system, launch **Ponemah**.
2. Select **Hardware | PhysioTel Digital Diagnostics...**


*Note:* If using *Ponemah v5.20*, select **Hardware | Diagnostics...**

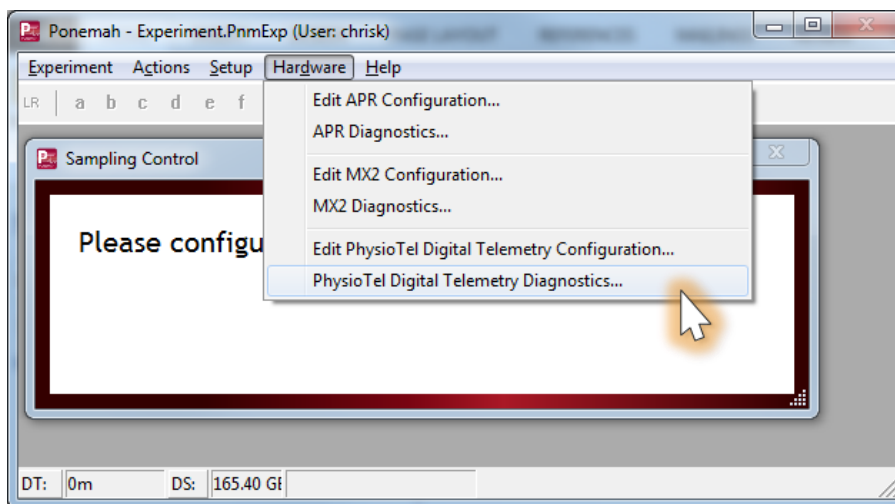


- Choose the CLC from the *Name – Model Name (Serial Number)* dropdown.  
*Note:* Select the Refresh List button if the CLC does not appear in the dropdown list.

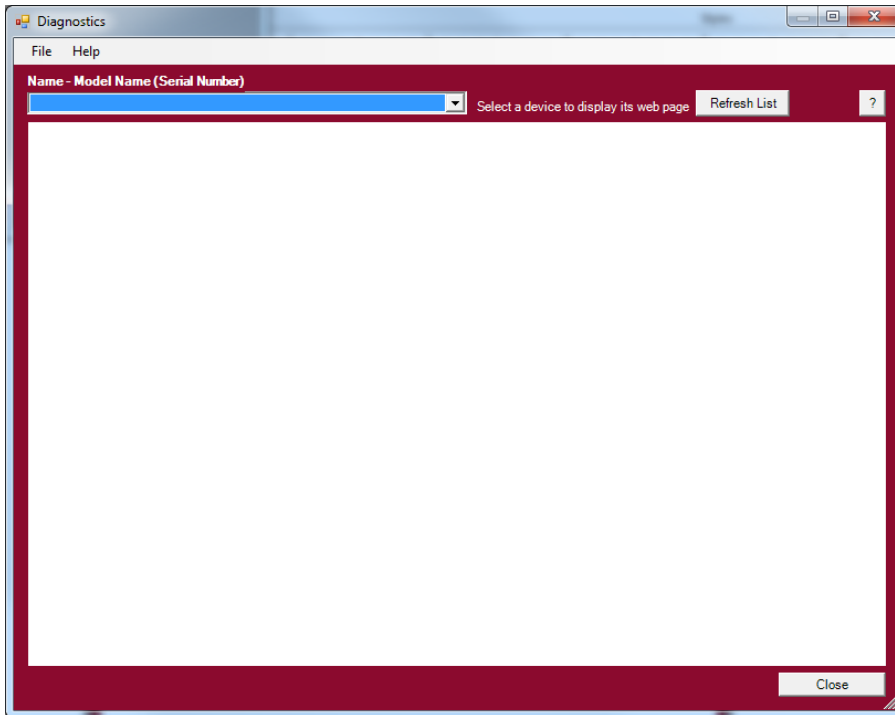


- Select the *Settings* link from the list view on the left side of diagnostics page.

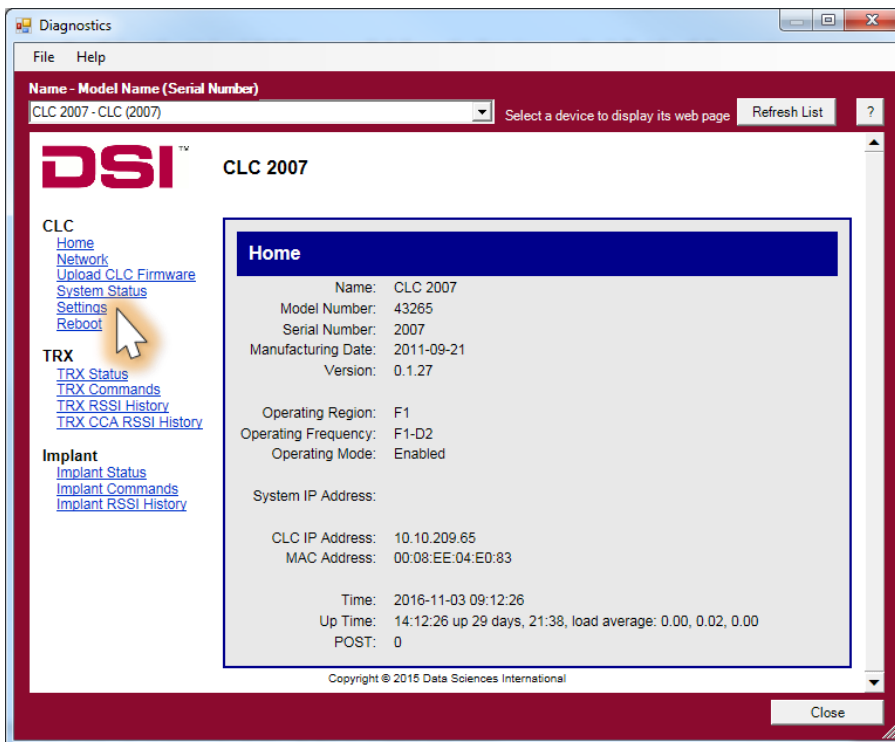


- Use the *RFMode* dropdown to select **Assessment**.

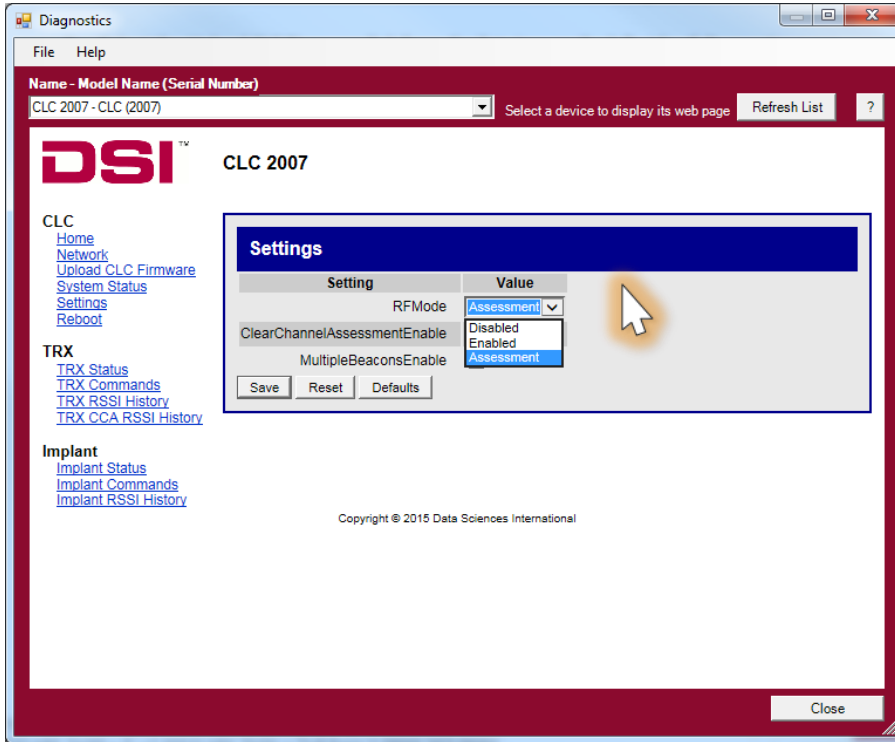

- Select the **Save** button.
- Select the *TRX RSSI History* link.

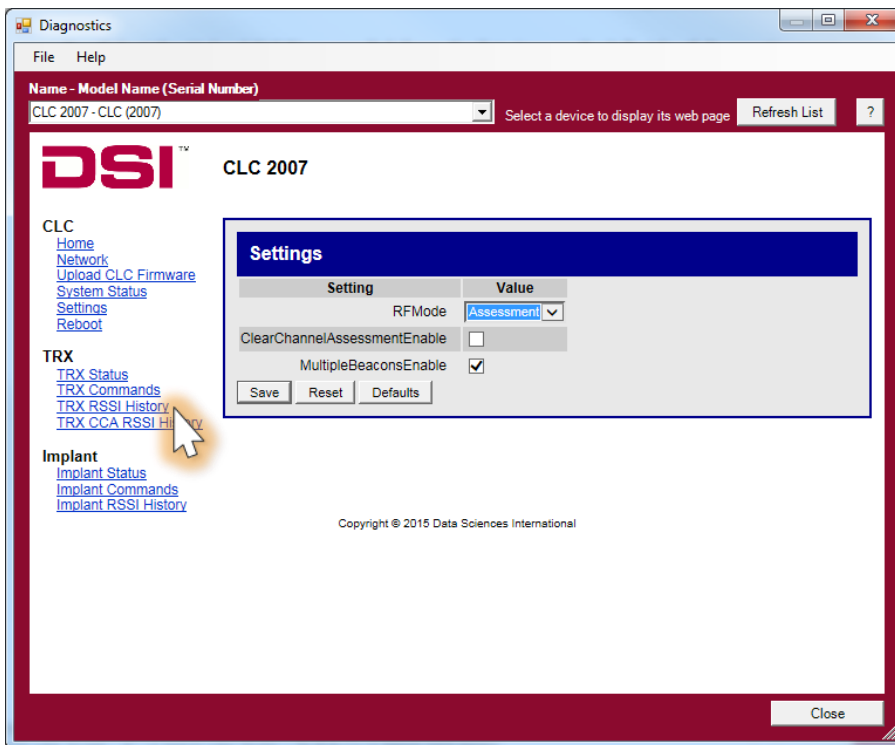


- At the top of the TRX RSSI History page, set **Auto refresh page in...** to **10** seconds.

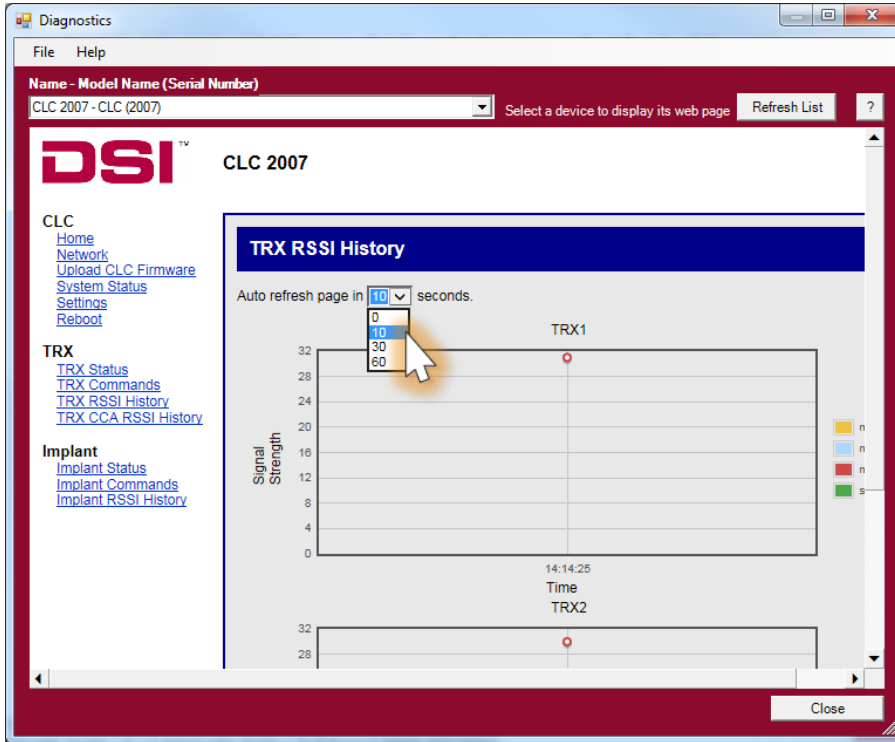


- Wait about 60 seconds for the graph below to populate with data related to your noise environment.
- Take a screenshot of the first graph on the page. Send the screenshot by e-mail to DSI Technical Support at [support@datasci.com](mailto:support@datasci.com). They will evaluate the graph and send further instructions.

11. Click the *Settings* link.

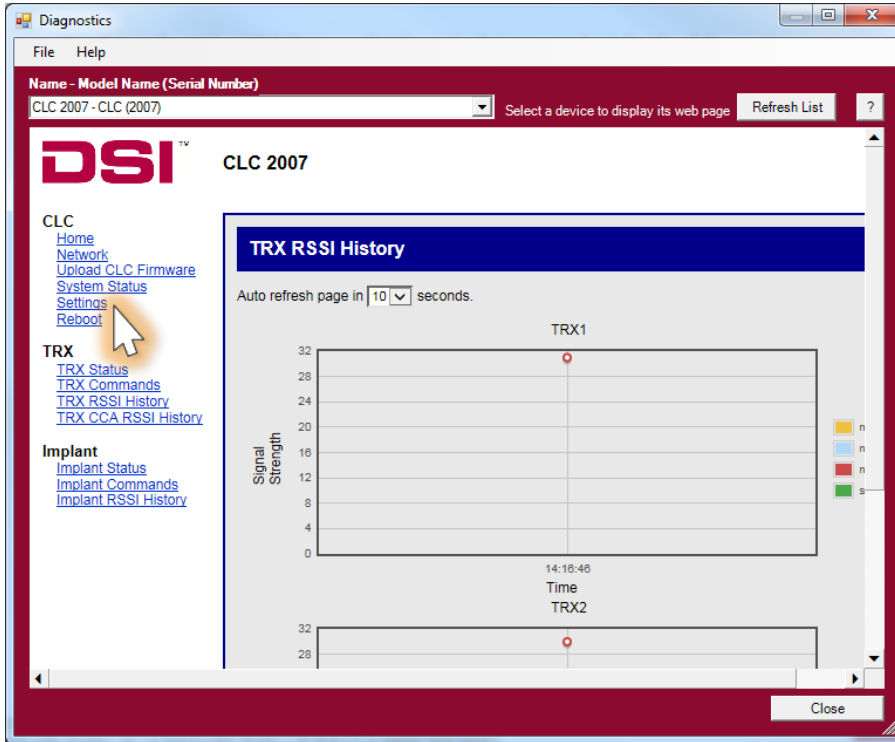

12. Use the *RFMode* dropdown to select **Enabled**.

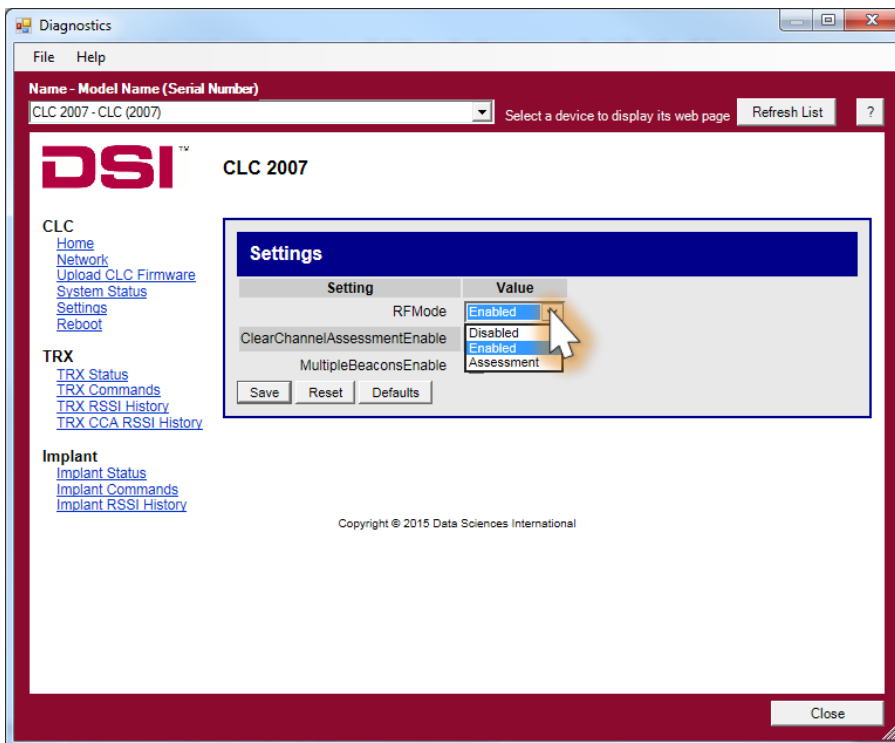


13. Select **Save**.

**TIP:** A screenshot can easily be taken in Windows 7 using the ‘snip’ tool. Click the Windows ‘Start’ Button. Type **snip** and the snip program should appear at the top of the list. Using this tool, you can select your assessment graph with a box, and then save it to your hard drive.

**EVALUATING THE GRAPH YOURSELF (NOT RECOMMENDED)**

1. Examine the first graph on the TRX RSSI History page. Look at the line created by the red dots; this refers to the Mean Noise Floor in your room.
2. Record the rough value of the Mean Noise Floor on the graph. In the example below, the Mean Noise Floor is about 9.

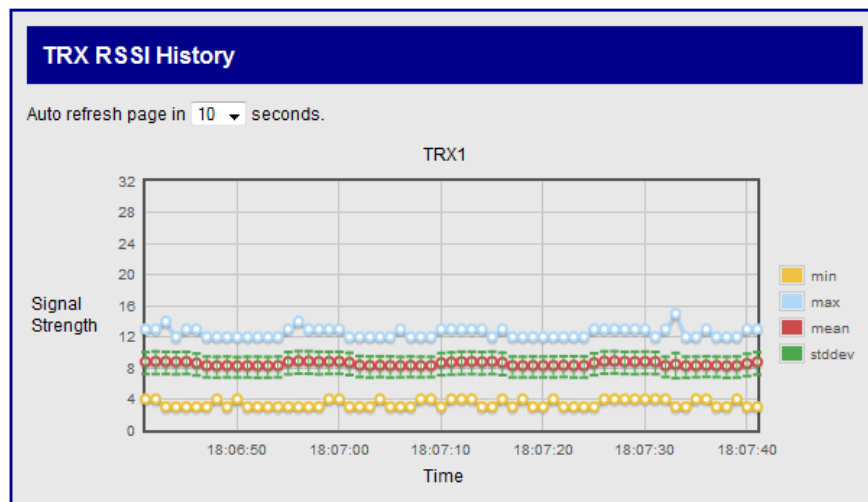

3. Next, record the width of the standard deviation (green hashes). In the example above, the deviation is about 2 (1 level above the mean and 1 level below it).
4. Calculate the necessary RSSI Threshold for the TRX and Implant as follows:

$$TRX\ RSSI\ Threshold = Mean\ Noise\ Floor + 3 * Std\ Dev$$

$$Implant\ RSSI\ Threshold = TRX\ RSSI\ Threshold - 3$$

Example:

$$TRX\ RSSI\ Threshold = (9) + 3 * (2)$$

$$TRX\ RSSI\ Threshold = 15$$

$$Implant\ RSSI\ Threshold = 15 - 3$$

$$Implant\ RSSI\ Threshold = 12$$

5. Last, these values will need to be converted from decimal to hexadecimal. This can be done in Windows by using the calculator function. Open the Windows Calculator.
6. In the calculator, click **View | Programmer**. Type in the TRX RSSI Threshold value from above.
7. Click the bubble button next to **Hex** on the left side of the calculator. The number should change to the HEX equivalent. In the example above, **15** becomes **F**.

8. Record this number or letter. If the calculator only gives a single digit or number, record a leading zero. In the example above, the output **F** should be written **0F** instead.
9. Refer to the technical note *Adjusting the RSSI Threshold for a Transceiver (TRX) and an Implant* for instructions on where to input the number.