

# TRX FIRMWARE INSTALL INSTRUCTIONS


Prior to installing, it is recommended to review the TRX Firmware Release Notes. The following are the instructions to install the firmware update.

## ACCESSING TRX FIRMWARE


The firmware file is typically provided on a CD from DSI. If the firmware was e-mailed by DSI, save the attached firmware file to a location on your hard drive. Make a note of this location for **Step 8** below.

## INSTALLING TRX FIRMWARE


1. Launch **Ponemah**.
2. Select the **Hardware | PhysioTel Digital Telemetry Diagnostics...**  
*Note: If using Ponemah v5.20, select **Hardware | Diagnostics...***

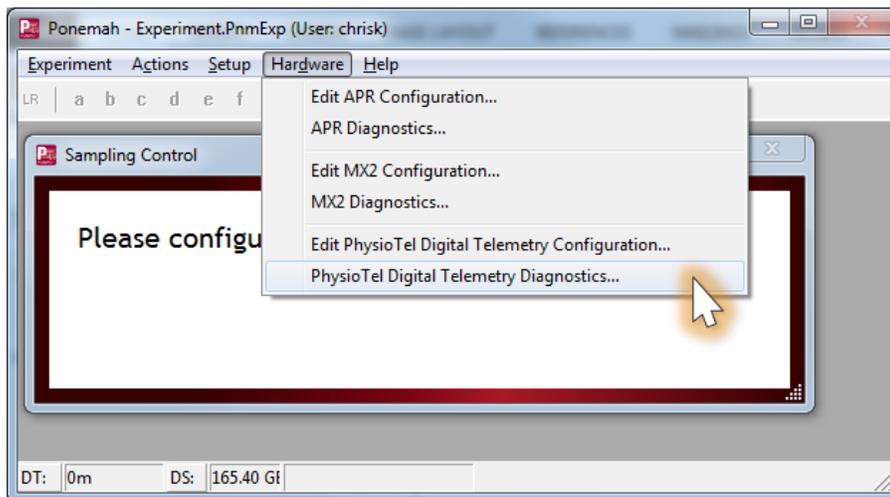


- Choose the CLC from the *Name – Model Name (Serial Number)* dropdown.  
*Note:* Select the Refresh List button if the CLC does not appear in the dropdown list.

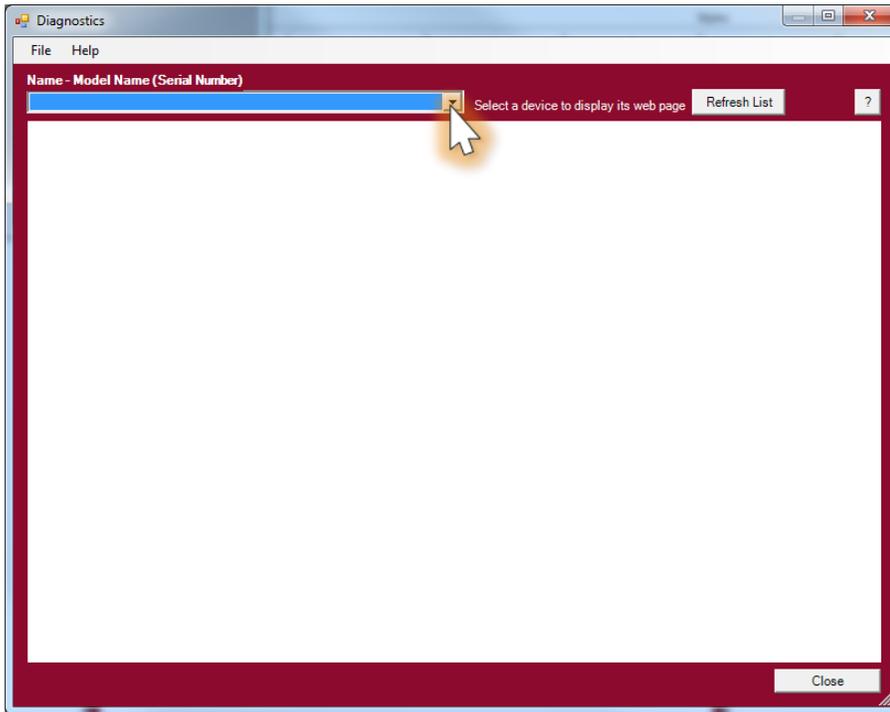


- Select the *TRX Commands* link from the list view on the left side of the diagnostic page.

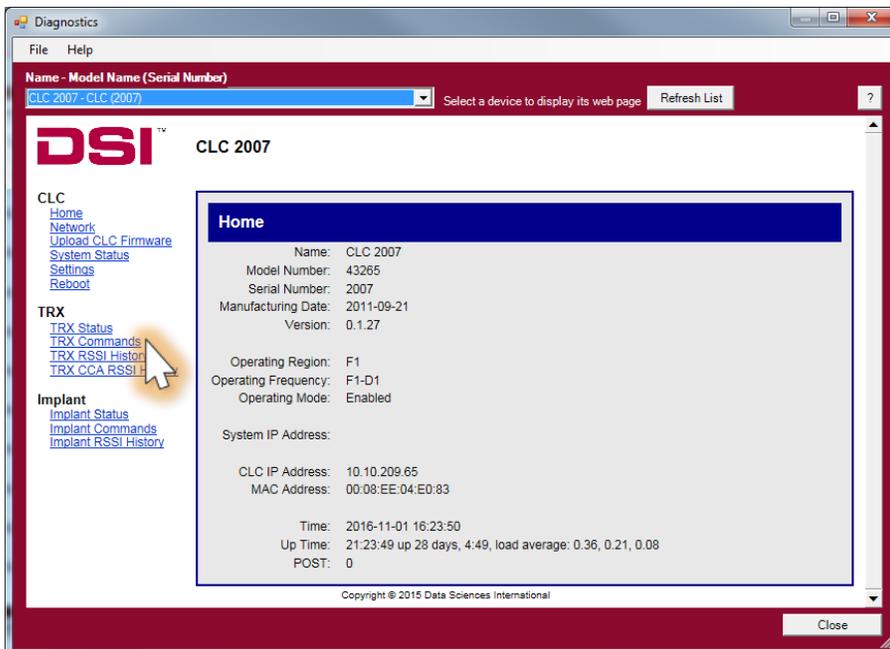


- Select the TRX number for the firmware update from the **TRX No.**  
*Note:* The TRX number corresponds to the CLC jack that the TRX is connected to.

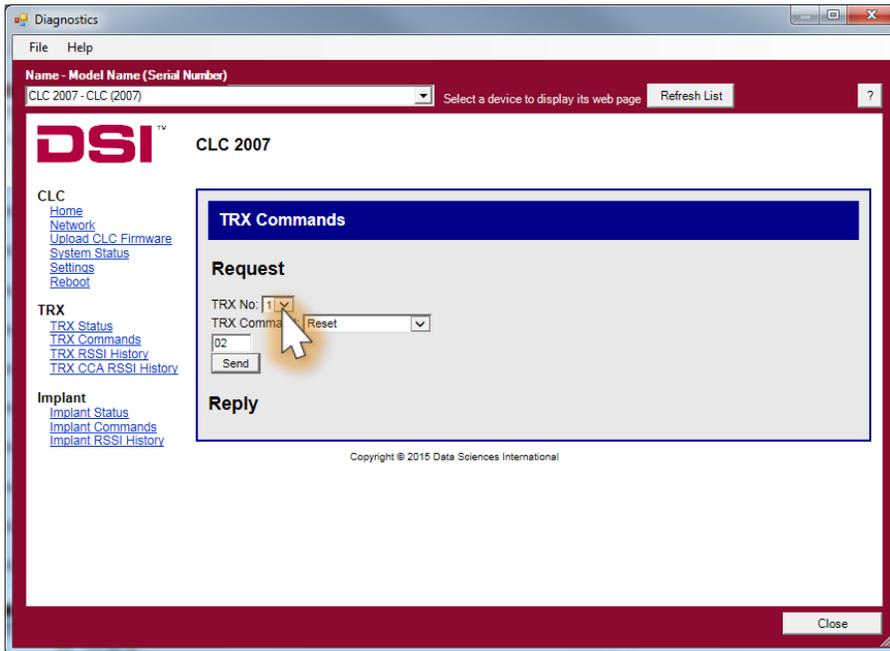


- Select **Upload Firmware** from the *TRX Command* dropdown.

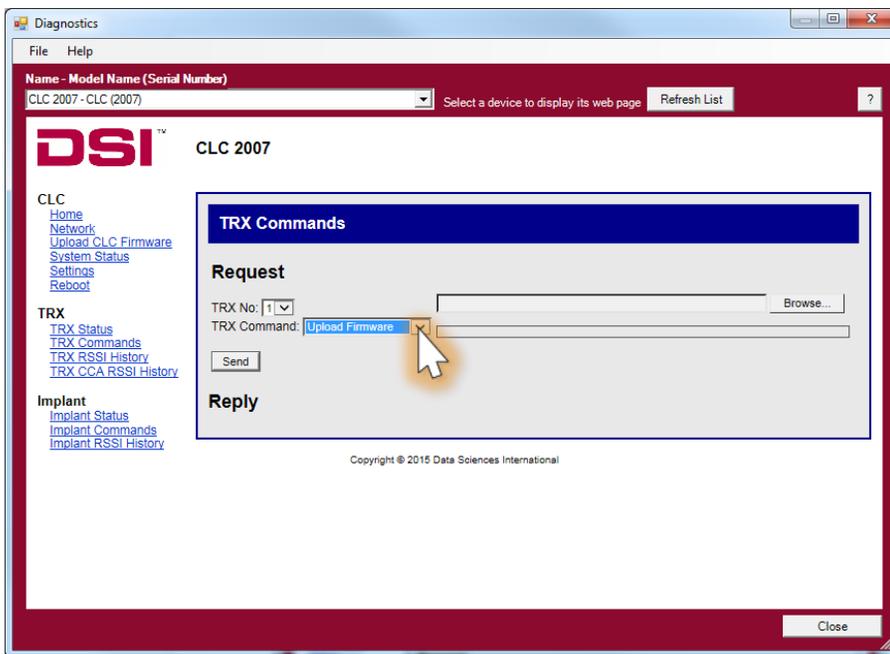

- Click the **Browse** button and navigate to the firmware file that was provided.

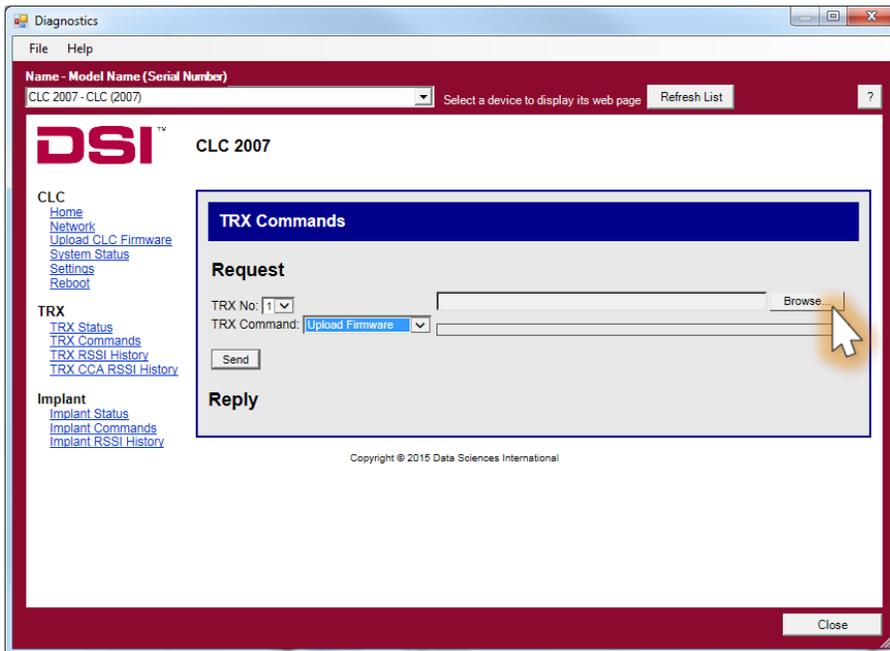

- Navigate to the file location for the TRX firmware, select it, and click OK.
- Click **Send**.

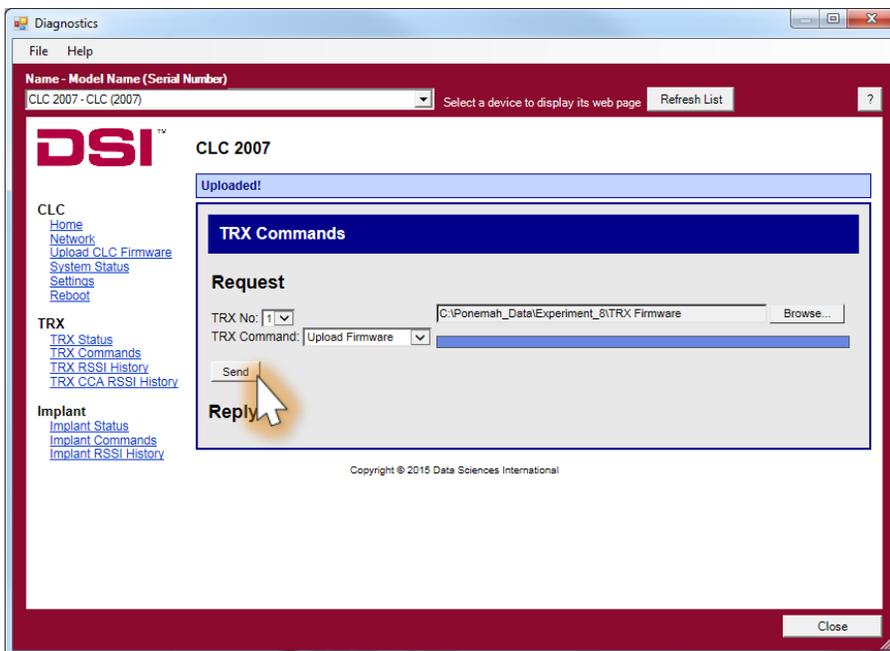

- After the firmware has been updated, verify the firmware revision by clicking on the **TRX Status** link from the left menu.