

PRODUCT RELEASE NOTES

Product: Ponemah™ Physiology Platform

Model: Ponemah Client

Version: 5.51

Build: 10622.2

Date: August, 2023

Product Release Notes for Ponemah Physiology Platform version 5.51 indicate revisions made to the Ponemah core application and analysis modules since version 5.50. For information regarding changes to the software from previous versions, please refer to the Release Notes folder located on install key or in the manuals and release notes section of the DSI support website.

Compatibility

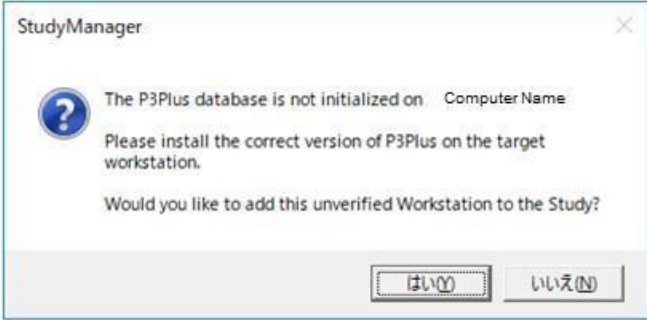
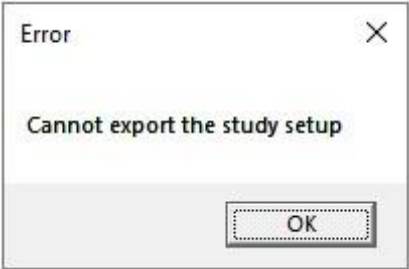
Ponemah v5.51 is compatible with the 64-bit version Windows 10 version 1607 and later. It is also compatible with the 64-bit version of Microsoft Office 2007 or greater.

Note: Microsoft .NET framework 3.5.1 and 4.7.2 required.

Fixes

Reference #	Ponemah Version	Description	Disposition
60218	5.51	Fixed an issue that could cause signals from implants of the model L11 and L21 to have inappropriate calibration ranges when swapping implants on an existing protocol. This results in a display of unrealistic ranges.	Fix
60975	5.51	Fixed an issue that could cause values for Custom Derived Parameters in the list view not to update when making changes to the analysis template used. Changing the logging rate would cause the CDP values to update.	Fix
60979	5.51	Remedied an issue that would cause the remaining battery life of a DSI implantable device not being displayed in the Implant Status dialog.	Fix

Known Issues

Reference #	Description
45933	<p>When using Japanese Language Pack with Windows 10 Operating System, the following error may be displayed when attempting to add a Workstation to Study:</p>  <p>This is caused by special characters being used in the data path name. The recommended workaround is to change the Windows display language to English.</p>
46259	<p>When trying to concatenate multiple RAW files created by achieving RAW file limit to cause file incrementation, data break lines may be seen at the rollover point. Additionally, a time overlap error may be displayed.</p>
47231	<p>If previous version of Ponemah have been installed on the workstation and the C:\Ponemah_Data\ folder includes subfolders with customized permissions configured for P3Users, then installing Ponemah v5.4x or v5.50 will result in removal of the customer permissions.</p>
48407	<p>When attempting to export a Study setup or print LOGs, an Error message may be displayed indicating that the function cannot be executed. E.g.</p>  <p>To correct this issue, navigate to C:\Ponemah\Utilities\ and run the AccessDatabaseEngine_x64.exe.</p>
48408	<p>When using PhysioTel Digital or MX2 configuration with HD implants, the Battery Voltage channel must be enabled for Battery Days Remaining (Digital)/Battery Time ON (HD) to display properly in Derived Parameter List Views and Implant Status dialogs.</p>

