

Certain hotkeys have been made available within Review to help facilitate quicker execution of certain functions from the keyboard and/or mouse without having to click additional icons or access menu options. Please note that they hotkeys are currently available only within Ponemah Review and cannot yet be augmented by the user. Please note: these hotkeys are only compatible with Ponemah v5.20-SP8 and later. For available hotkeys in v6.30, please refer to the v6.30 User Manual Data Review section.

Key Combination	Description of Action
Zoom/Graphical Navigation	
<Alt> + <S>	Synchronize Primary, Trend, and Scatter graphs with associated Derived List View. When using this hotkey from the Scatter graph, only the Derived List View is synchronized. To sync the other graphs, double-click the selected row.
<+> (on number pad)	Expand graph
<-> (on number pad)	Compress graph
<Ctrl> + <+> (on number pad)	Expand graph by factor
<Ctrl> + <-> (on number pad)	Compress graph by factor
<Shift> + Mouse wheel	Expand/Compress Scroll Wheel up – zoom in (expand time span) Scroll Wheel down – zoom out (compress time span)
<Ctrl> + Mouse wheel	Expand/Compress graph by a factor Scroll Wheel up – zoom in (expand time span) Scroll Wheel down – zoom out (compress time span)
Mouse Wheel	Scroll waveform window left/right
Page Up	Scrolls waveform window to the left (increase in time) by screen time
Page Down	Scrolls waveform window to the right (decrease in time) by screen time
<Shift> + <F9>	Auto scale waveform for selected pane within graph page
<Ctrl> + <F9>	Auto scale waveforms for all panes within graph page
<Ctrl> + <1> through <5>	Change time scale to predefined entries defined within Advanced settings of Options Application Configuration: Review.Graph.Times n
<Ctrl> + <6> through <9>	Scale Y axis of all graph panes within the graph page to one of the 4 defined entries within Advanced settings of Options Application Configuration Review.Graphs.High.Scales n

Key Combination	Description of Action
	Review.Graphs.Low.Scalen
<Shift> + <6> through <9>	Scale the Y axis of the selected graph pane to one of the 4 defined entries (see above)
<Ctrl> + <Z>	Undo zoom on waveform (only one level)
<↑>	Move up one trace area – activate trace area
<↓>	Move cursor down one trace area – activate trace area
<←>	Move cursor left
<→>	Move cursor right
<Ctrl> + <←>	Move cursor the left by a factor
<Ctrl> + <→>	Move cursor the right by a factor
<Shift> + <F6>	Moves to the previous Parser Segment
<Shift> + <F7>	Moves to the next Parser Segment
Events	
<Shift> + <a> through <j>	Add Event to currently selected Group
<Shift> + <Ctrl> + <a> through <j>	Add Event to all Groups
 on Event	Delete selected Event
Mark Placement/Positioning	
<Tab>	Move to next validation mark
<Shift> + <Tab>	Move to previous validation mark
<Shift> + <x>	Add Bad Data Mark
<Shift> + <F2>	Move to next Bad Data Mark
<Shift> + <F3>	Move to previous Bad Data Mark
<Shift> + <F4>	Find next change (only if Diff Marks are on)
<Shift> + <F5>	Find previous change (only if Diff Marks are on)
<Ctrl> + <←> or <→> on Validation Mark, Event, or Bad Data Mark	Move selected mark left or right
<Alt> + <v>	Executes reanalysis for the visual region of the <i>selected</i> channel within a Primary graph.
 on Validation Mark	Delete selected mark Note: If on main Cycle mark (e.g. R-mark) deletes complete cycle
 on Bad Data Mark	Delete selected Bad Data Mark
<Alt> + <r>	Deletes all Bad Data Marks within the selected channel and automatically executes a reanalysis of the channel

Key Combination	Description of Action
<Alt> + 	Adds Bad Data Marks across the visible region of data within Primary and Trend graphs for the currently <i>selected</i> channel
<Alt> + <z>	Adds Bad Data Marks across the visible region of data within Primary and Trend graphs for <i>all groups and their associated channels</i>
ECGPRO	
<Ctrl> + <n>	Add New Template Cycle
<Shift> + <u>	Delete Unmatched Cycles on a Channel (User will be prompted for confirmation)
Data Insights	
Mouse Wheel	Scroll waveform window left/right
<Ctrl> + Mouse wheel	Expand/Compress graph by a factor Scroll Wheel up – zoom in (expand time span) Scroll Wheel down – zoom out (compress time span)
Left-click-and-drag	Zoom in to selected area
<Enter> or Double left-click	Sync to Match Result within Primary graph, Trend graph, Scatter Graph and Logged Line of Derived List View.
<Ctrl> + <r>	Reject selected Match Result(s)
<Ctrl> + 	Add Bad Data Marks around selected Match Result(s)
<Ctrl> + <d>	Delete cycle validation marks from selected Match Result(s)