



SEND 3.1: Standard for Exchange of Nonclinical Data



Data Sciences International

**119 14TH STREET NW
SUITE 100**

ST. PAUL, MN 55112 USA

(651)481-7400 • 1(800)262-9687

www.datasci.com

PN: MU00285-001 REV01

OVERVIEW

This manual highlights how DSI's SEND 3.1 can be used to produce output needed for data submission to the FDA. This document will provide an overview of the system components and basic operation.

Table of Contents

Welcome	3
Background	3
SEND Setup from PPP3 Setup	4
SEND dialog	4
Groups	7
Data Reduction	8
Using Study Protocol Option and SEND	13
SEND Configuration When Using Study	13
SEND Manager	16
SEND Viewer and Output	18
SEND Export To Excel	19
SEND Viewer	20
Send Export to CSV	25
Synchronize SEND Data	27
Synchronize SEND Data	27
Synchronize SEND Study Data	29
High Level Process Flow for Generating SEND Output	29
Troubleshooting	30
Messages Posted Due to Incorrect SEND Configuration	31
Technical Support	34
Appendix 1	35

WELCOME

The structure of this manual was designed to sequentially guide you through setup of the various SEND components and ultimately generating the standardized output needed for data submission. This manual references functionality that may not be fully detailed here. It may be necessary to reference additional Ponemah manuals for detailed feature information. This information can be found under the Help menu within Ponemah.

WHAT YOU WILL BE LEARNING

1. Determine what functions are needed in order to generate SEND information
2. Understand how to configure SEND specific features
3. Generate SEND output

BACKGROUND

SEND (Standard for Exchange of Nonclinical Data) is a standardized method to exchange data between organizations and for data submission to the FDA or other regulatory bodies for non-clinical data. Detailed information on formatting, standards, terminology and other information can be found at <https://www.cdisc.org/>.

In order to produce output using the SEND module, users must ensure the following:

- SEND Output or SEND SQL Server Output has been enabled in license file
- Derived Parameters selected
 - Not all derived parameters generated by Ponemah are currently defined by CDISC. Derived parameters will only be present in SEND output if those parameters defined by CDISC have been enabled in the Channel Input Setup menu.
 - See Appendix 1 for a list of currently accepted parameters
- Data reduction enabled for subjects
 -
 - The currently defined calculations are Average (Avg), Maximum (Max), and Minimum (Min)
- SEND.XML file is created and listed in the Ponemah directory or the location where the Ponemah software has been installed (automatically created with software installation)
 - The SEND.XML file provides a mapping of the terminology between the Ponemah derived nomenclature and the expected SEND controlled terminology. This file can be modified by the user to include additional Ponemah parameters. Additionally, as parameters are updated by CDISC, this file can be easily modified to include the new information.

This manual assumes users have read and understand the core components of the Ponemah software application. In some instances, SEND is built upon existing menus or functions in the software. If additional information is needed on functionality outside of the SEND application, please review the manuals located under the Help pull-down menu in the Ponemah application.

SEND SETUP FROM PPP3 SETUP

In order to create SEND output, users must have enabled Data Reduction, selected derived parameters in the Channel Input Setup menu that correspond to output designated by CDISC, and performed a Save of the data while in Review (Save Marks Section, Save Derived Data, or Save Run).

SEND output can be generated within our outside of the Study Protocol Option (Study). This section will describe how to set up components that are used both with and without Study. Specific use with Study will be described in a later section.

SEND DIALOG

SEND can be configured in idle mode and saved as part of the protocol or modified after collection in Review. To modify SEND settings, select P3 Setup from the Setup pull-down menu and click on SEND in the tree-view menu on the left-hand side of the dialog. By default, SEND will be disabled. In order to edit, select "Enable SEND Output". Once enabled, fields will become editable.

Definitions for the functions displayed in the SEND dialog are listed below.

PPP3 Setup - SEND

PPP3 Setup SEND

Channel Input Setup
Template Setup
Groups
Events
Experimental Protocol Header
Data Reduction Setup
Variability Analysis
Graph Setup
Settings
SEND

Enable SEND Output

Enable Excel Output

Study Name: Study Name Edit Study Name

Study Day: 1 Edit Study Day

Workup: Edit Workup

Output Server: Test connection...

Laboratory Name:

Append Subject ID to Study Name Subject Separator:

Update SEND database during Acquisition

Consciousness State: CONSCIOUS

Position: UNCONSTRAINED

EG Method: 12 LEAD STANDARD

CV Method: INTRAVASCULAR

RE Method: HEAD OUT PLETHYSMOGRAPH

OK Cancel Apply

Figure 1. – SEND dialog from the PPP3 Setup menu.

1 – Enable SEND Output

Toggles the ability to modify the editable fields displayed in Figure 1 and controls whether or not SEND output is turned on or off.

2 – Enable Excel Output

SEND output can be ported to Microsoft Excel along with the derived, log, and subject information currently created when saving a Marks section, Derived Data, or a Run (within Study). If enabled, a new worksheet for each SEND domain will be created within the Excel workbook that contains SEND information along with the expected Derived, Data Reduction, and Experimental Log information. Refer to section SEND Export To Excel for details on the output generated when enabling this feature.

If this option is not checked, SEND output can be generated by creating a CSV file using SEND Viewer (see SEND Viewer below) only with the SEND SQL Output Server option enabled in the license file. The SEND Output option only allows output to Excel.

3 – Study Name

Allows the user to create a unique study name for the SEND output. If using the Study Protocol Option, this information will be automatically pulled by the software. However, the option to modify the Study name is possible by checking the box next to the edit field. This may be needed to further define a set of data or update an incorrect Study name that may have been created.

4 – Study Day

The day of collection may be entered here. Note that CDISC does not recognize Day 0. If using the first day of the study collection, a “1” should be entered. If using the Study Protocol Option, this information will be automatically pulled by the software. However, the option to modify the Study Day is possible by checking the box next to the edit field. This may be needed to further define a set of data or update an incorrect Study entry.

5 – Workups

Workups can be manually defined, or If using the Study option, this information can be automatically pulled by the software. If pulled from Study, users can modify the workup name used in Study by checking the box next to the edit field.

6 – Output Server

Permits defining a dedicated server or workstation as the primary repository for SEND data. Additionally, the SEND instance name, as it is installed with the application, must be appended to the server/workstation name. Example naming convention would be “server name\P3Plus_V2”. P3Plus_V2 is the SQL Instance used by Ponemah for SEND related data.

NOTE: It is important to note that this functionality is only available if the SEND SQL Server Option has been enabled in the license file. If only SEND Output is enabled, this feature will not be available.

7 – Laboratory Name

Allows a unique identifier to be manually entered which will populate the EGNAM variable in the SEND output for the EG Domain

8 – Append Study ID to Subject Name

Will append the Subject ID to the Study Name in the SEND output. By default, the separator is “_” but can be changed. This modifies USUBJID in the SEND output for the domain.

9 – Update SEND database during acquisition

Generation of SEND data will typically be generated during Review after post processing the data. However, there may be instances where post analysis of the data is not needed such as in short duration, Tox applications. Enabling this checkbox will allow data generated during acquisition to be used by SEND.

10 – Consciousness State

Pull-down field allows selection of CDISC controlled terminology that represents the state of the animal, such as conscious or unconscious. These definitions are managed by the SEND.XML file.

11 – Position

Presents a pull-down menu allowing the selection of CDISC controlled terminology that represents the subjects position from unconstrained to various restrained states. These definitions are managed by the SEND.XML file.

12 – EG Method

Allows the selection of the CDISC controlled terminology that represents the ECG lead presentation used for collection. These definitions are managed by the SEND.XML file.

13 – CV Method

Allows the selection of CDISC controlled terminology for definition of pressure measurements. These definitions are managed by the SEND.XML file.

14 – RE Method

Permits selection of CDISC controlled terminology such as Head Out Plethysmograph, Whole Body Plethysmograph, or Invasive to describe the respiratory system being used. These definitions are managed by the SEND.XML file.

See Technical Support for details, if modification to the SEND.XML file is needed.

GROUPS

The Groups dialog allows the user to control the Group name and dose which is associated with a given Subject. In order to edit the Study Subject Name or Study Dose, click in the edit field under the appropriate column. Figure 2 shows subjects one and two (Group A and Group B) with the names and doses modified.

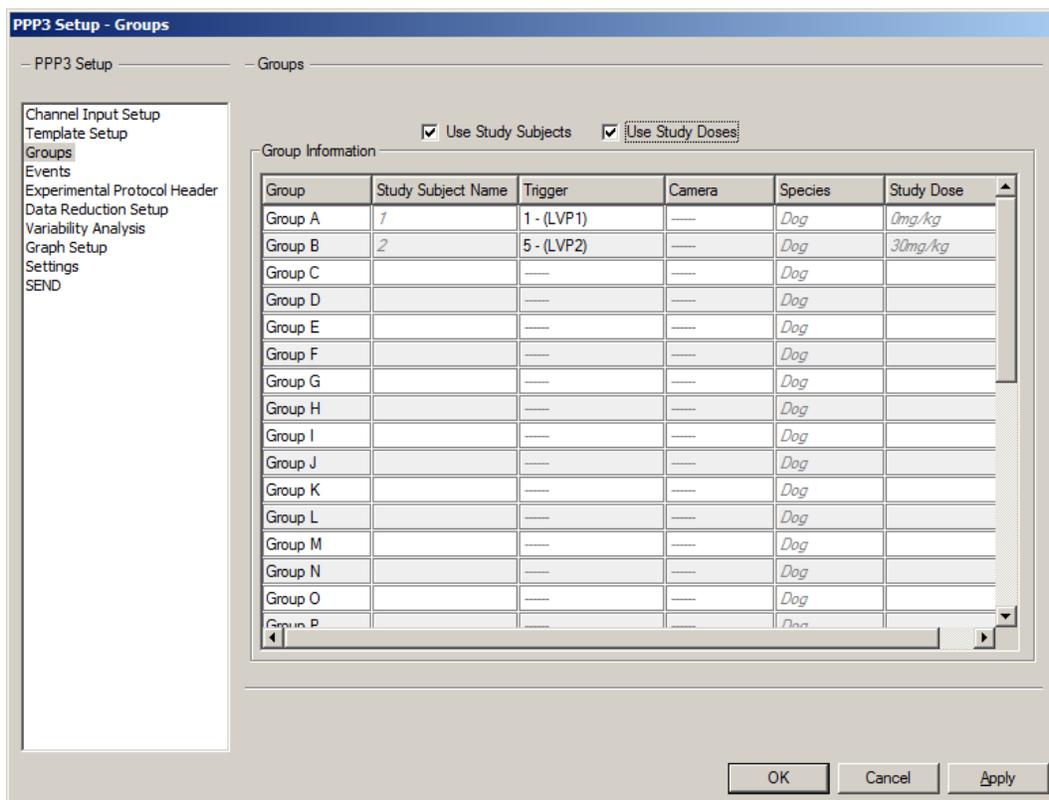


Figure 2 – Groups dialog from the PPP3 Setup menu.

Providing a unique name in the Groups dialog will propagate the subject information to SEND.

Below shows example SEND output where Group A Study Subject Name was changed in the Groups tab (Figure 2). This identifier, ID 123, is shown under USUBJID in Figure 3.

Two check boxes are available at the top of the Groups dialog, Use Study Subjects and Use Study Doses. These two features are related to data generated using the Study Protocol Option. If enabled, Subject and Dose information entered in Study will be automatically pulled into the Groups tab. When these features are enabled, manually editing Study Subject Name and Study Dose will not be permitted. However, it may be necessary to update or correct information that was entered incorrectly into the Study Protocol Option. If needed, uncheck these boxes and enter the appropriate information manually. When these boxes are unchecked, the column headers revert back to the default values of Name and Dose. If not using the Study Protocol Option, these features do not need to be enabled.

1 – Use Study Subjects

Subject information defined within the Study Protocol Option can be automatically pulled into SEND output by selecting this checkbox

2 – Use Study Doses

Dose information entered within the Study Protocol Option will be pulled into SEND by selecting this checkbox

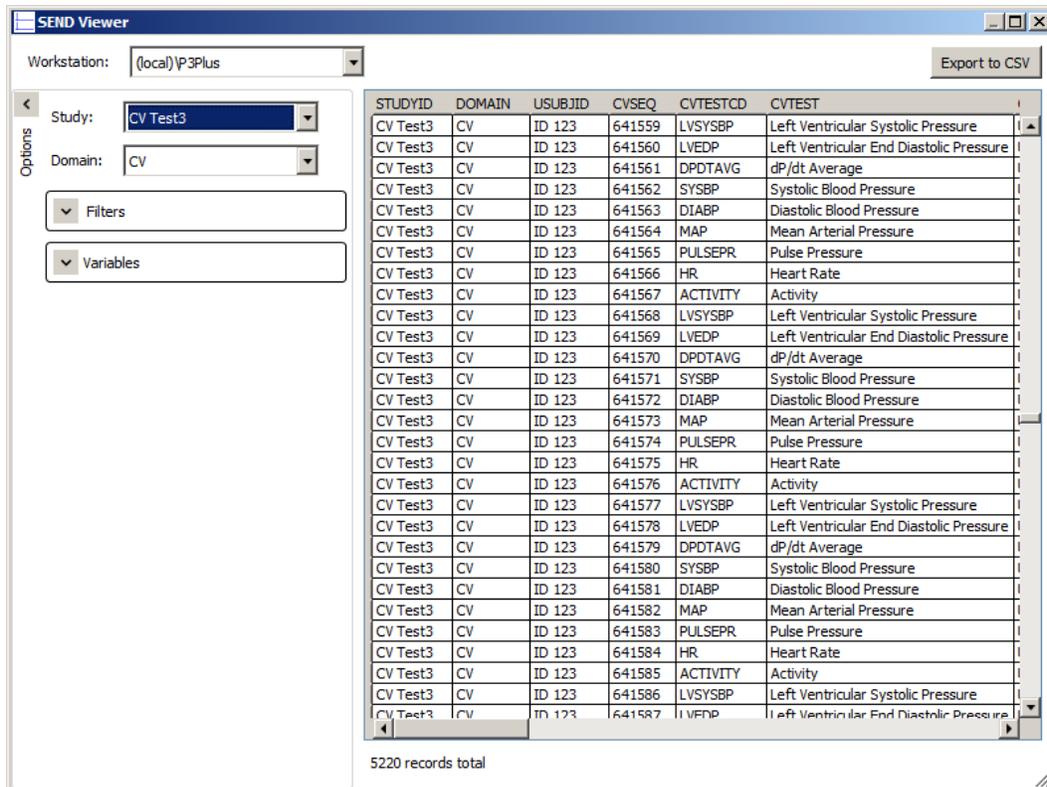


Figure 3. –Example SEND Viewer dialog showing Group information pulled into SEND (USUBJID).

DATA REDUCTION

Data Reduction must be enabled for all desired subjects in order to generate SEND output. Only parameters defined by CDISC will be available in the SEND output.

With SEND enabled, users can no longer create user specified labels as that will be handled automatically to conform to SEND requirements. When SEND has been enabled, Data Reduction will display “SEND compatibility enabled” at the top of the dialog to notify users that they will be complying with SEND requirements and some functions will not be able to be modified.

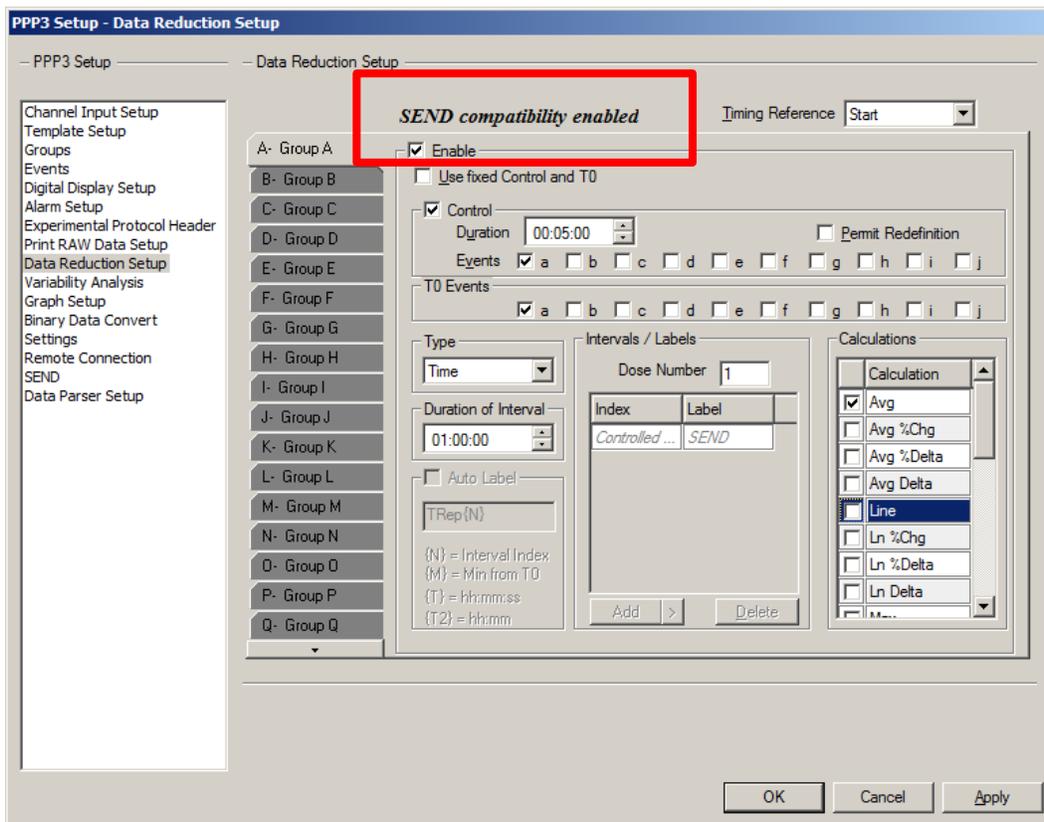


Figure 4. – Data Reduction dialog showing SEND enabled.

Several features have been added to the Data Reduction dialog to provide details on the data generated for SEND output.

1 – Timing Reference

The Timing Reference selection determines the point on each data reduction interval that will be used to report timing information for the interval (Figure 5). This is used to define Data Reduction labels along with TPT and ELTM.

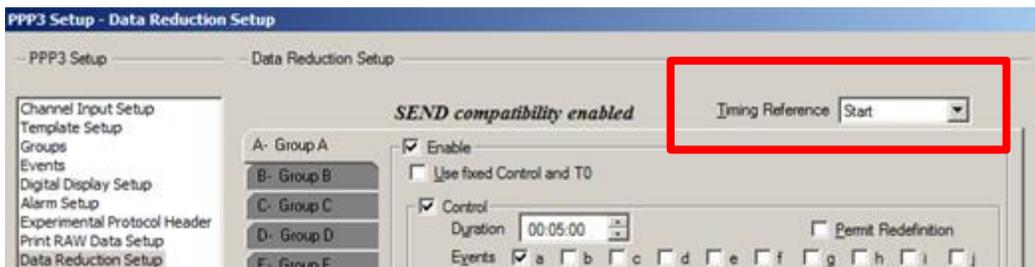


Figure 5. – Data Reduction showing Timing Reference selection.

Three selections are available which include Start, Middle, and End. The selection chosen will update the Data Reduction label to display the time information at the start of the logging interval, the middle of the interval, or the end of the interval as shown in Figures 6 and 7.

Figure 6 shows labels starting at 0s and incrementing by 10 minutes (10m, 20m, 30m) as dictated by the “Duration of Interval” chosen by the user.

The screenshot shows a software window titled "Ponemah - SampleECGData.RVW (User: mjb) - [Data Reduction: a]". The interface includes a menu bar (Functions, Setup, SEND, Data Parser, Options, Window) and a toolbar. A table displays data reduction parameters. The 'Label' column is highlighted with a red box. The table contains the following data:

Label	Start Time	End Time	Duration	1: Sys	1: LVEDP	1:+dP/dt	2: Sys	2: Dia	2: Mean	2: PH	2: HR
T0	00:00:00										
0s Postdose	00:00:00	00:10:00	0000:10:00								
Avg				118.40	5.5708	3310.2	117.87	81.248	99.849	36.621	74.423
10m Postdose	00:10:00	00:20:00	0000:10:00								
Avg				111.20	4.7871	3023.4	112.41	76.446	94.795	35.967	60.68
20m Postdose	00:20:00	00:30:00	0000:10:00								
Avg				115.28	5.6705	3043.0	117.20	78.698	98.12	38.50	64.35
30m Postdose	00:30:00	00:40:00	0000:10:00								
Avg				117.80	6.0854	3062.5	119.69	81.78	100.54	37.906	69.436
40m Postdose	00:40:00	00:50:00	0000:10:00								
Avg				127.26	7.4468	3525.0	128.53	89.39	109.08	39.142	86.194
50m Postdose	00:50:00	01:00:00	0000:10:00								
Avg				121.04	6.2823	3302.2	121.49	82.506	102.14	38.985	74.726
1h Postdose	01:00:00	01:10:00	0000:10:00								
Avg				143.49	12.489	4083.5	142.69	98.744	120.67	43.951	106.27
1h 10m Postd...	01:10:00	01:20:00	0000:10:00								
Avg				119.46	7.1452	3384.1	120.36	83.135	102.49	37.225	76.565
1h 20m Postd...	01:20:00	01:30:00	0000:10:00								

Figure 6. – Data Reduction label showing “Start” selected for the Timing Reference.

Figure 7 shows the same Start Time, End Time, Duration and averages for the data, but the label has been modified to display the “Middle” selection. Since the duration of the averaging interval is 10 minutes, selecting Middle will update the label information to specify the middle of that logging interval. Figure 7 shows those updated labels starting at 5m and incrementing every 10 minutes (5m, 15m, 25m, etc.).

Ponemah - SampleECGData.RVW (User: mjb) - [Data Reduction: a]

Functions Setup SEND Data Parser Options Window

Label	Start Time	End Time	Duration	1: Sys	1: LVEDP	1:+dP/dt	2: Sys	2: Dia	2: Mean	2: PH	2: HR
T0	00:00:00	00:00:00	0000:10:00								
5m Postdose	00:00:00	00:10:00	0000:10:00	118.40	5.5708	3310.2	117.87	81.248	99.849	36.621	74.423
Avg											
15m Postdose	00:10:00	00:20:00	0000:10:00	111.20	4.7871	3023.4	112.41	76.446	94.795	35.967	60.68
Avg											
25m Postdose	00:20:00	00:30:00	0000:10:00	115.28	5.6705	3043.0	117.20	78.698	98.12	38.50	64.35
Avg											
35m Postdose	00:30:00	00:40:00	0000:10:00	117.80	6.0854	3062.5	119.69	81.78	100.54	37.906	69.436
Avg											
45m Postdose	00:40:00	00:50:00	0000:10:00	127.26	7.4468	3525.0	128.53	89.39	109.08	39.142	86.194
Avg											
55m Postdose	00:50:00	01:00:00	0000:10:00	121.04	6.2823	3302.2	121.49	82.506	102.14	38.985	74.726
Avg											
1h 5m Postdose	01:00:00	01:10:00	0000:10:00	143.49	12.489	4083.5	142.69	98.744	120.67	43.951	106.27
Avg											
1h 15m Postd...	01:10:00	01:20:00	0000:10:00	119.46	7.1452	3384.1	120.36	83.135	102.49	37.225	76.565
Avg											
1h 25m Postd...	01:20:00	01:30:00	0000:10:00								

Current Group Name Change Reference Group Name Change Raw File C:\Ponemah_Data\SampleECGData\SampleECGData

Figure 7. – Data Reduction label showing “Middle” selected for the Timing Reference.

2 – Dose Number

Dose Number is manually edited and refers to the dose that will be referenced in the SEND output. The dose number listed here will be placed in the SEND output under TPT. By default, the value is “1” but can be modified to represent the desired dose if multiple dosing events occur in a single day. If SEND is not enabled in the SEND dialog, this field will not be editable.

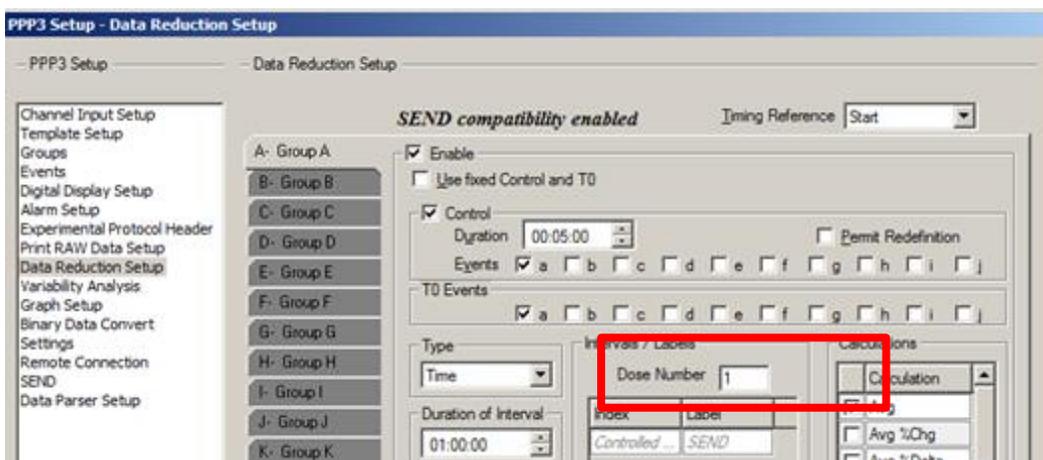


Figure 8. – Data Reduction dialog showing Dose Number edit field.

3 – Used Fixed Control and T0

The Use Fixed Control and T0 function allows the designation of the start of the Data Reduction without the use of an event. Entering in date and time into the T0 Time edit fields will trigger the start of Data Reduction. When using Time as the Data Reduction “Type”, all labels will be controlled by Ponemah for SEND compliance.

A time of 9:00AM was entered into the T0 Time field. In the list view dialog, the “0s Postdose” label is entered and controlled by Ponemah and the first logged line of data is at 9:00AM. All subsequent intervals are based on the Duration of Interval configured.

The screenshot shows the 'PPP3 Setup - Data Reduction Setup' dialog box. The 'SEND compatibility enabled' checkbox is checked. Under 'Use fixed Control and T0', the 'Control' checkbox is unchecked, and the 'T0' section is active. The 'T0 Time' is set to 5/17/2007 at 9:00:00 AM. The 'Type' is set to 'Time' and the 'Duration of Interval' is 00:10:00. The 'Intervals / Labels' table shows a single entry: Index 'Controlled ...' and Label 'SEND'. The 'Calculations' list includes 'Avg', 'Avg %Chg', 'Avg %Delta', 'Avg Delta', 'Line', 'Ln %Chg', 'Ln %Delta', and 'Ln Delta'. The 'OK', 'Cancel', and 'Apply' buttons are visible at the bottom.

Below the dialog box is a screenshot of the 'Ponemah - CV Pro.RVW' application window. The 'Data Reduction: a' table is displayed, showing the results of the data reduction process. The table has columns for Label, Start Time, End Time, Duration, and various data points (1: Sys, 1: LVEDP, 1: HR, 1: +dP/dt, 2: Sys, 2: Dia, 2: Mean). The first row is '0s Postdose' starting at 09:00:00 on 05/17/2007. Subsequent rows show 'Avg' values for 10m, 20m, 30m, 40m, and 50m post-dose intervals, followed by 'Avg' values for 1h, 1h 10m, and 1h 20m post-dose intervals.

Label	Start Time	End Time	Duration	1: Sys	1: LVEDP	1: HR	1: +dP/dt	2: Sys	2: Dia	2: Mean
T0	0001:01:12 09:00:00 05/17/2007									
0s Postdose	0001:01:12 09:00:00 05/17/2007	0001:11:12 09:10:00 05/17/2007	0000:10:00							
Avg				143.64	12.826	106.97	4080.8	142.74	99.035	120.88
10m Post...	0001:11:12 09:10:00 05/17/2007	0001:21:12 09:20:00 05/17/2007	0000:10:00	119.18	6.5803	77.269	3435.4	119.90	82.338	101.84
Avg				116.34	3.4051	74.239	3672.5	115.41	75.577	95.801
20m Post...	0001:21:12 09:20:00 05/17/2007	0001:31:12 09:30:00 05/17/2007	0000:10:00							
Avg				119.50	3.1541	79.70	3804.6	117.94	75.991	96.84
30m Post...	0001:31:12 09:30:00 05/17/2007	0001:41:12 09:40:00 05/17/2007	0000:10:00							
Avg				118.41	6.4415	66.17	3441.6	118.18	79.706	98.018
40m Post...	0001:41:12 09:40:00 05/17/2007	0001:51:12 09:50:00 05/17/2007	0000:10:00							
Avg				121.06	5.8841	73.593	3605.2	119.84	79.766	99.597
50m Post...	0001:51:12 09:50:00 05/17/2007	0002:01:12 10:00:00 05/17/2007	0000:10:00							
Avg				137.34	5.8307	96.607	4829.7	136.31	89.754	112.78
1h Postdose	0002:01:12 10:00:00 05/17/2007	0002:11:12 10:10:00 05/17/2007	0000:10:00							
Avg				137.10	5.3037	91.027	4460.5	136.10	91.636	113.74
1h 10m Po...	0002:11:12 10:10:00 05/17/2007	0002:21:12 10:20:00 05/17/2007	0000:10:00							
Avg										

Figure 9. – Data Reduction with Use fixed Control and T0 enabled.

Parser Segments can also be configured and used for Data Reduction. When using Parser Segments, the ability to add user defined labels is permitted. Figure 10 shows the Add button expanded under the Intervals/Labels field. All labels manually entered here will be fixed in the Data Reduction list view and any additional reduction line labels will be handled by Ponemah.

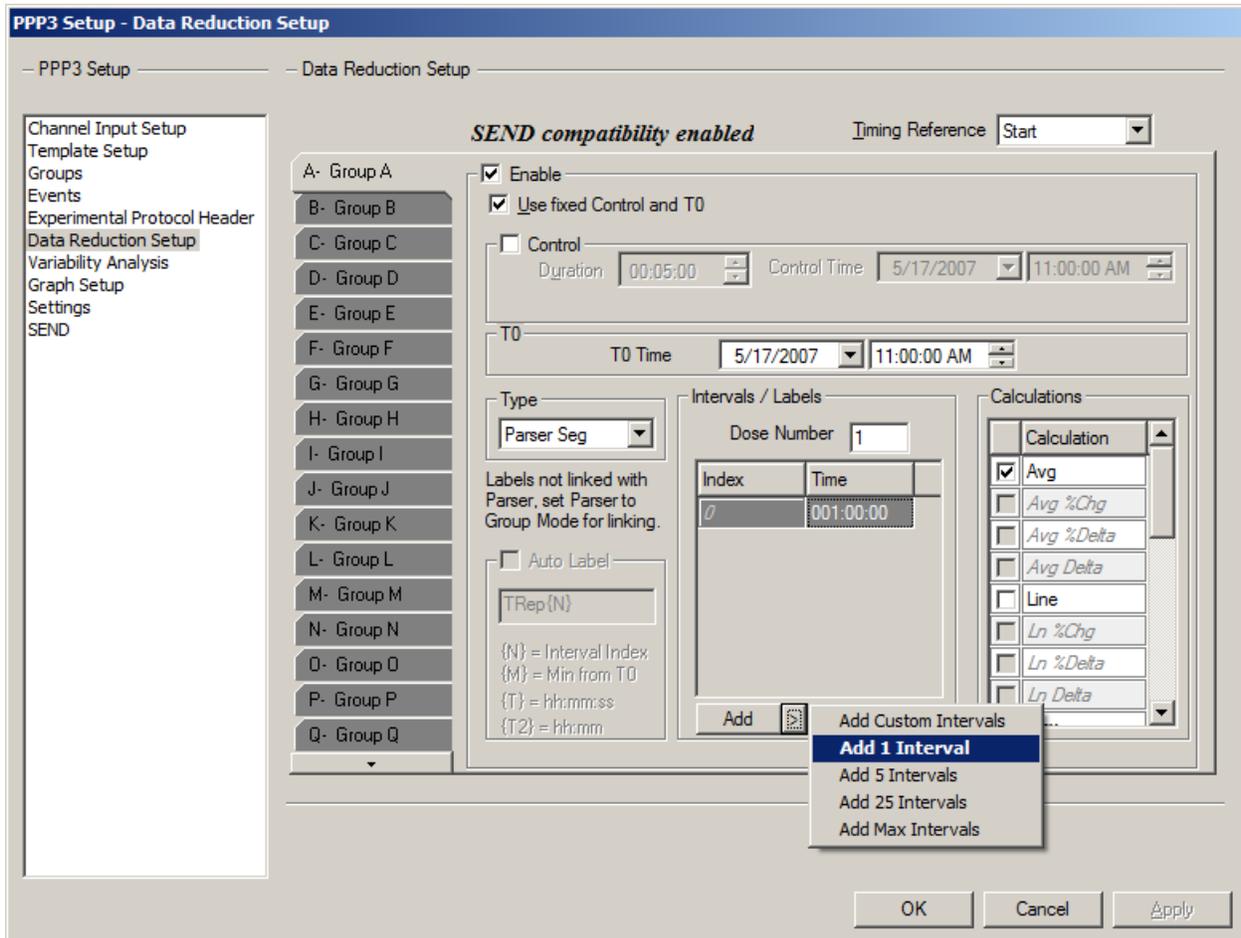


Figure 10. – I know I will add something here.

USING STUDY PROTOCOL OPTION AND SEND

The following outlines the differences in SEND options when using the Study Protocol Option. Refer to previous sections outlining the setup required for the SEND, Groups and Data Reduction dialogs to ensure requirements for generating SEND output have been configured.

SEND CONFIGURATION WHEN USING STUDY

If using the Study Protocol Option, the study configuration has been previously configured and saved as part of the Study prior to any collections. This includes Data Reduction, Subject IDs and other information. When using the Study Protocol Option, information can be automatically pulled into the SEND dialog and other menus. However, it

may be necessary to modify information prior to generating SEND output to correct entry errors or further define a set of SEND generated data.

Functions specific to the Study Protocol Option in the SEND dialog include the ability to automatically load the Study Name, Study Day, and Workup. In order for the Study information to be displayed in the SEND dialog, you must have accessed a Review file within the Run Study dialog. With the Review file open, navigating to the SEND dialog will display the information for the Study that is currently accessed. Additionally, a Review file created within Study can be opened outside of Study with access to this information.

When the checkboxes have been disabled as in Figure 11, the information created using the Study option will be pulled into the SEND dialog. However, no modification of these fields will be allowed. Figure 11 shows CVStudy and Study Day 1 pulled from the Study configuration. Workup has no information since Workups were not configured as part of this Study.

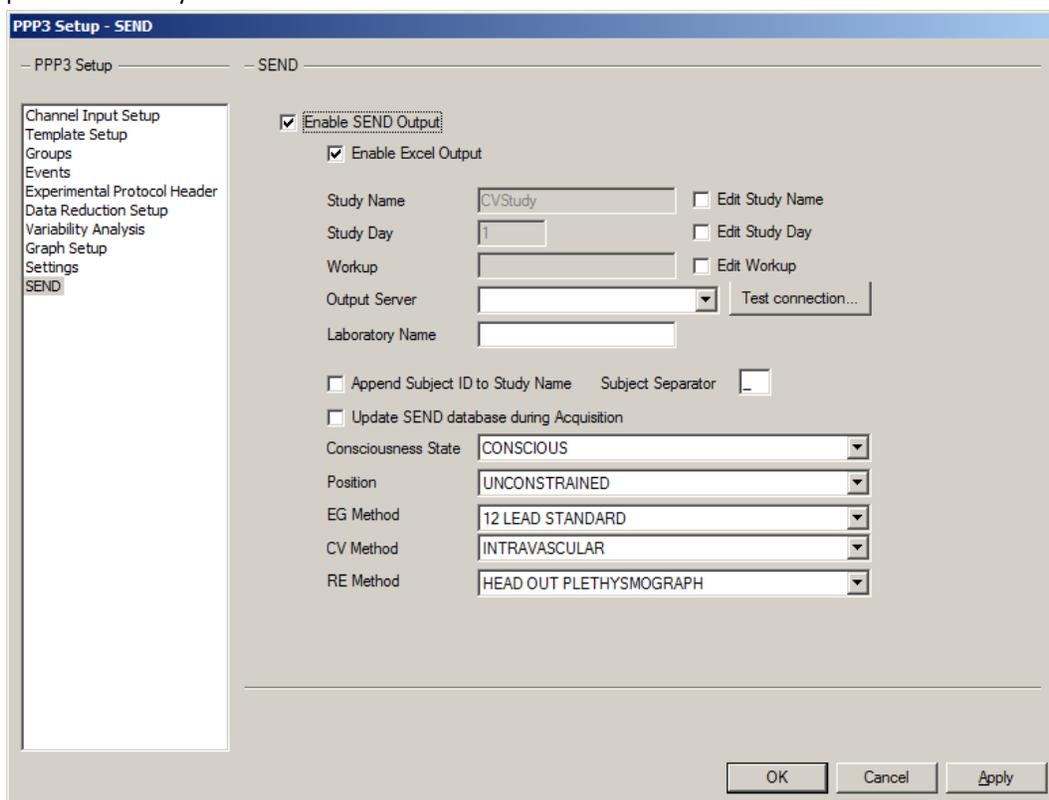


Figure 11. – Study Name, Study Day, and Workup pulled from the Study Protocol Option

To modify Study information, place a check in the Edit Study Name, Edit Study Day, or Edit Workup checkbox. If the check marks beside these options are enabled as shown in Figure 12, these fields may be manually modified from their original values from when the Study was created.

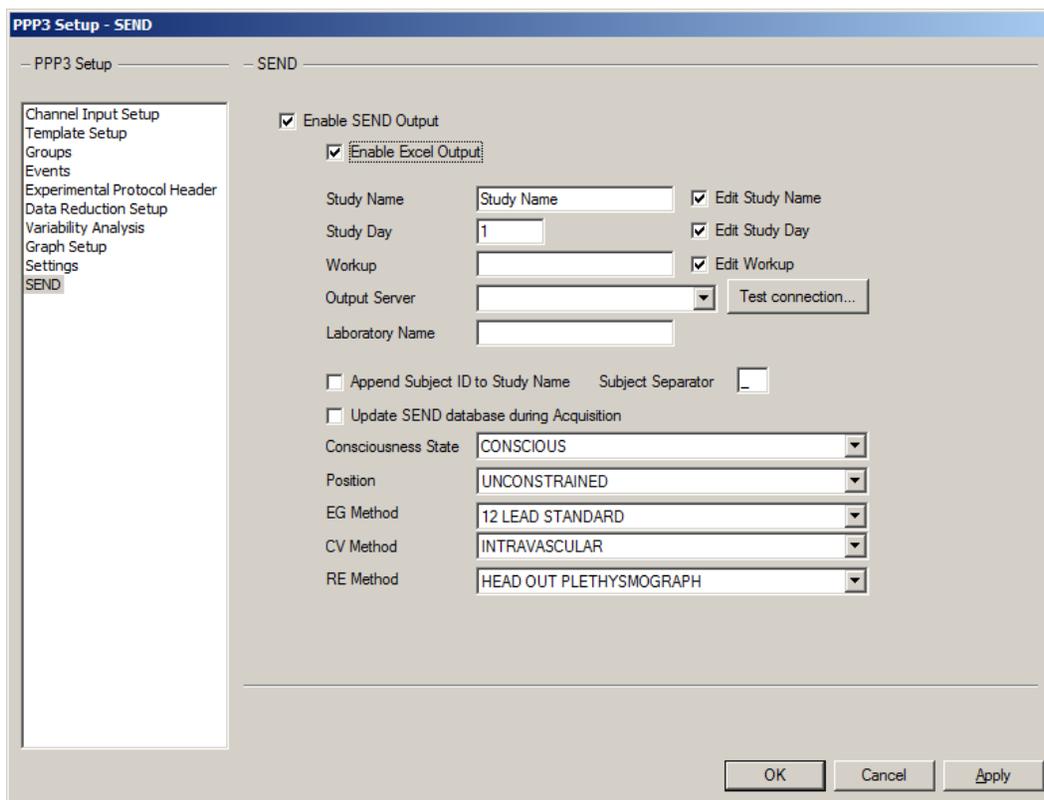


Figure 12. – Edit Study Name, Edit Study Day, and Edit Workup allows Study information to be edited

As with the SEND dialog, the Groups dialog will also automatically fill Study Subject Name and Study Dose information. Deselecting Use Study Subjects and Use Study Doses will allow modification of the Study Protocol Option generated information. To modify, simply type in the desired information for each Subject (Group).

The information populated here is based on the information contained in the Review file opened when in the Run Study dialog or outside of Study. For example, Study Dose information will update appropriately if running a Latin Square design based on the information contained in the Review file that was opened.

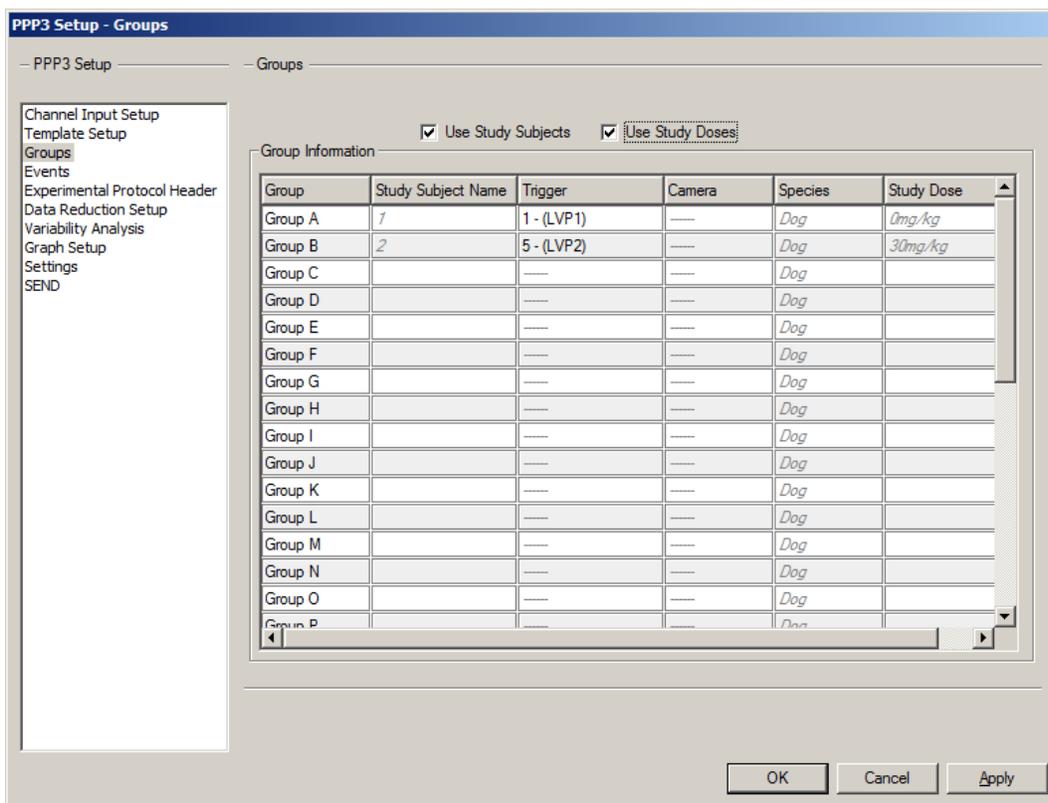


Figure 13. – Groups dialog with subject and dose information pulled from Study.

SEND MANAGER

SEND Manager is available from the SEND pull-down menu in Ponemah while in idle mode and allows you to access collections that have been previously saved. If in Review, SEND Manager will not be available (selection will be grayed out).

NOTE: It is important to note that this functionality is only available if the SEND SQL Server Option has been enabled in the license file. If only SEND Output is enabled, this feature will not be available.

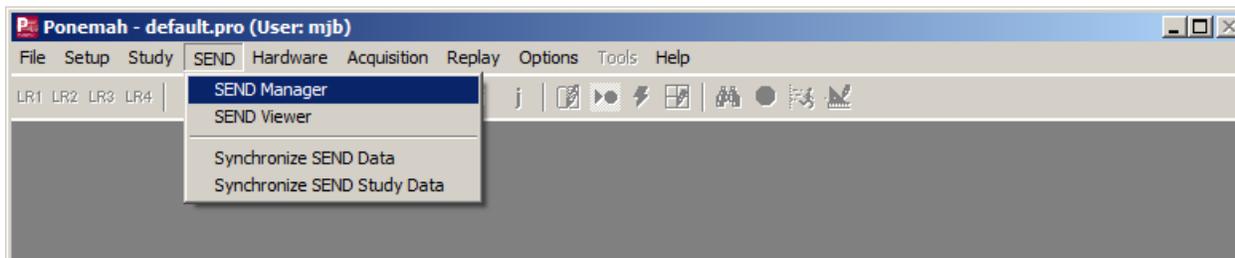


Figure 14. – SEND Manager access point from the SEND pull-down menu.

In addition to listing the collections available with SEND information, SEND Manager has backup and restore capabilities as described below. Each time a unique SEND study name is created, and data is saved, that study name will be listed in the dialog below.

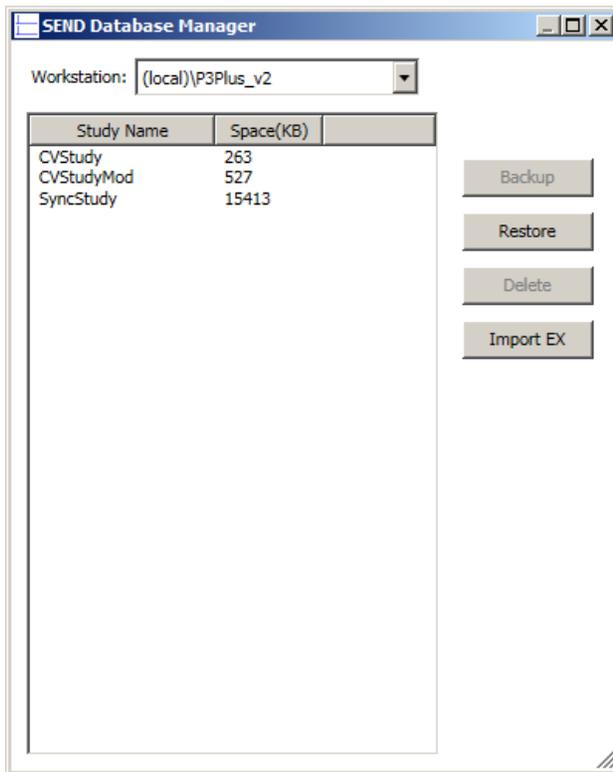


Figure 15. – SEND Manager dialog.

1 – Workstation

Lists the local Ponemah SQL Instance, P3Plus_v2, by default. If networked, users can select the P3Plus_v2 Instances on other networked workstations.

2 – Study Name and Space

Lists all available study names with previously saved SEND information and the amount of disk space that each study occupies.

3 – Backup

Allows the studies to be backed up and moved to other workstations or for archival purposes. The backup dialog is shown below. Use the typical Windows navigation to browse to the desired folder location and specify a file name for the backup file. The file type will automatically default to “.sendbkup”.

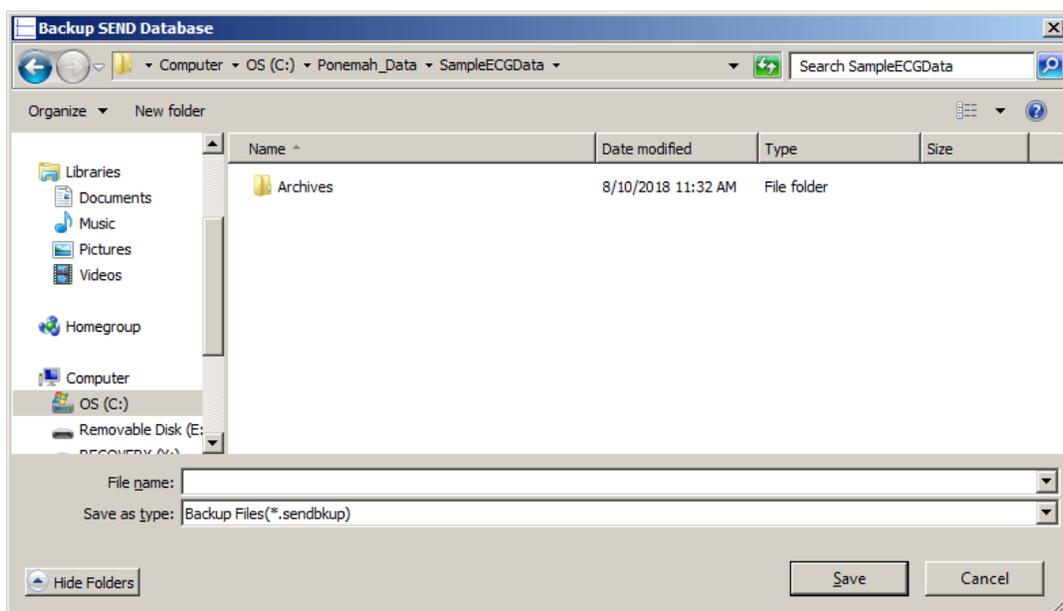


Figure 16. – Example SEND backup file being saved.

4 – Restore

Allows studies to be restored from other workstations or reopened from an archival location. This feature functions similar to that described under Backup shown above.

5 – Delete

The Delete function will remove the Study from the list in the SEND Manager dialog and will also remove any associated data with the study from the SQL database.

6 – Import EX

The ability to import information is possible by creating a CSV file and using the Import EX function in SEND Database Manager. The required columns for successful import are STUDYID, USUBJID, EXSTDTC, EXSTDY, and optionally EXENDTC may be provided. Exposure domain information is not available in Ponemah, although some Variables require exposure data. Exposure data can be brought into Ponemah by importing an EX domain that has been exported from a LIMS system or other repository.

The following variables will report NULL unless an Exposure domain containing actual dose times is imported (DY, ENDY, RFTDTC). The EX domain can be imported at any point prior to exporting the CSV.

SEND VIEWER AND OUTPUT

The following sections detail the output options available for SEND generated data and the tools which can be used to view and modify the output.

SEND EXPORT TO EXCEL

To generate SEND data to Excel, the Enable Excel Output function must be selected in the SEND setup dialog. Excel output is available for both license file options, SEND Output and SEND SQL Server Output. Additionally, users must have enabled Data Reduction and selected derived parameters in the Channel Input Setup menu that correspond to output designated by CDISC (refer to SEND Setup from PPP3 Setup section for configuration information).

SEND output will be generated and included in the Microsoft Excel workbook created upon Saving a Marks Section (Run) or Derived Data within a Review session along with other Ponemah data related worksheets. Note that not all Ponemah derived parameters are defined by CDISC. Only those CDISC defined variables will be present in the output generated.

	C	D	E	F	G	H	I	J	K	L	M	
1	USUBJID	CVSEQ	CVTESTCD	CVTEST	CVPOS	CVORRES	CVORRESU	CVSTRESSC	CVSTRESSN	CVSTRESSU	CVSTAT	CVRE
2	CV	Group A	LVSYSBP	Left Ventricular Systolic Pressure	UNCONSTRAINED	118.40	'mmHg'	118.40	118.40	'mmHg'	NULL	NULL
3	CV	Group A	LVEDP	Left Ventricular End Diastolic Pressure	UNCONSTRAINED	5.5708	'mmHg'	5.5708	5.5708	'mmHg'	NULL	NULL
4	CV	Group A	HR	Heart Rate	UNCONSTRAINED	74.319	'bpm'	74.319	74.319	'bpm'	NULL	NULL
5	CV	Group A	DPDTAVG	dP/dt Average	UNCONSTRAINED	3310.2	'mmHg/s'	3310.2	3310.2	'mmHg/s'	NULL	NULL
6	CV	Group A	SYSBP	Systolic Blood Pressure	UNCONSTRAINED	117.87	'mmHg'	117.87	117.87	'mmHg'	NULL	NULL
7	CV	Group A	DIABP	Diastolic Blood Pressure	UNCONSTRAINED	81.248	'mmHg'	81.248	81.248	'mmHg'	NULL	NULL
8	CV	Group A	MAP	Mean Arterial Pressure	UNCONSTRAINED	99.849	'mmHg'	99.849	99.849	'mmHg'	NULL	NULL
9	CV	Group A	PULSEPR	Pulse Pressure	UNCONSTRAINED	36.621	'mmHg'	36.621	36.621	'mmHg'	NULL	NULL
10	CV	Group A	ACTIVITY	Activity	UNCONSTRAINED	0.0164	"	0.0164	0.0164	"	NULL	NULL
11	CV	Group A	LVSYSBP	Left Ventricular Systolic Pressure	UNCONSTRAINED	111.20	'mmHg'	111.20	111.20	'mmHg'	NULL	NULL
12	CV	Group A	LVEDP	Left Ventricular End Diastolic Pressure	UNCONSTRAINED	4.7871	'mmHg'	4.7871	4.7871	'mmHg'	NULL	NULL
13	CV	Group A	HR	Heart Rate	UNCONSTRAINED	60.679	'bpm'	60.679	60.679	'bpm'	NULL	NULL
14	CV	Group A	DPDTAVG	dP/dt Average	UNCONSTRAINED	3023.4	'mmHg/s'	3023.4	3023.4	'mmHg/s'	NULL	NULL
15	CV	Group A	SYSBP	Systolic Blood Pressure	UNCONSTRAINED	112.41	'mmHg'	112.41	112.41	'mmHg'	NULL	NULL
16	CV	Group A	DIABP	Diastolic Blood Pressure	UNCONSTRAINED	76.446	'mmHg'	76.446	76.446	'mmHg'	NULL	NULL
17	CV	Group A	MAP	Mean Arterial Pressure	UNCONSTRAINED	94.795	'mmHg'	94.795	94.795	'mmHg'	NULL	NULL
18	CV	Group A	PULSEPR	Pulse Pressure	UNCONSTRAINED	35.967	'mmHg'	35.967	35.967	'mmHg'	NULL	NULL
19	CV	Group A	ACTIVITY	Activity	UNCONSTRAINED	NULL	NULL	NULL	NULL	NULL	'NOT DONE'	'INCC'
20	CV	Group A	LVSYSBP	Left Ventricular Systolic Pressure	UNCONSTRAINED	115.28	'mmHg'	115.28	115.28	'mmHg'	NULL	NULL
21	CV	Group A	LVEDP	Left Ventricular End Diastolic Pressure	UNCONSTRAINED	5.6705	'mmHg'	5.6705	5.6705	'mmHg'	NULL	NULL
22	CV	Group A	HR	Heart Rate	UNCONSTRAINED	64.359	'bpm'	64.359	64.359	'bpm'	NULL	NULL
23	CV	Group A	DPDTAVG	dP/dt Average	UNCONSTRAINED	3043.0	'mmHg/s'	3043.0	3043.0	'mmHg/s'	NULL	NULL
24	CV	Group A	SYSBP	Systolic Blood Pressure	UNCONSTRAINED	117.20	'mmHg'	117.20	117.20	'mmHg'	NULL	NULL
25	CV	Group A	DIABP	Diastolic Blood Pressure	UNCONSTRAINED	78.698	'mmHg'	78.698	78.698	'mmHg'	NULL	NULL
26	CV	Group A	MAP	Mean Arterial Pressure	UNCONSTRAINED	98.12	'mmHg'	98.12	98.12	'mmHg'	NULL	NULL
27	CV	Group A	PULSEPR	Pulse Pressure	UNCONSTRAINED	38.50	'mmHg'	38.50	38.50	'mmHg'	NULL	NULL

Figure 17. – Example output when enabling Excel in the SEND dialog.

If the collection contains data from more than one Domain, multiple worksheets will be created within the Excel workbook. Since the study shown in Figure 17 incorporated both CV and EG domain data, two worksheets are created in the Excel workbook with the appropriate information for each Domain.

SEND VIEWER

NOTE: It is important to note that this functionality is only available if the SEND SQL Server Option has been enabled in the license file. If only SEND Output is enabled, these features will not be available.

Selecting SEND Viewer from the SEND pulldown menu will display the following dialogs. In order to view SEND information, users must have enabled Data Reduction and selected derived parameters in the Channel Input Setup menu that correspond to output designated by CDISC (refer to SEND Setup from PPP3 Setup section for configuration information).

SEND Viewer allows selection and filtering of the study data and information for export to a CSV formatted file to be imported into other systems. Figure 18 shows the SEND Viewer dialog populated with data from the CV Domain and the associated CDISC controlled terminology displayed in the column headers. Scroll bars on the bottom and right side of the dialog allow the user to scroll and view the entire contents of the window.

The Workstation pulldown menu lists the local Ponemah SQL instance, P3Plus_v2, by default along with any designated Output Server that has been configured. If networked, users can select P3Plus_v2 Instances on other networked workstations or servers. This allows access to the local Ponemah SEND database as well as other Ponemah SEND databases on workstations that are accessible on the network. Selecting a workstation will display any previously collected information (studies) that is available for export. The pull-down menu will be populated with workstation information if studies have been created using the Study Protocol Option. All workstations that were configured for a given study will be loaded and available via the pull-down menu. The list of workstations available may change based on the study that is selected in the Manage Studies dialog.

Selecting the Options in the upper left-hand side of the dialog will expand and display filters that allow you to select the desired information for export. This includes the study, subjects, days and CDISC defined parameters.

SEND Viewer

Workstation: (local) P3Plus Export to CSV

STUDYID	DOMAIN	USUBJID	CVSEQ	CVTESTCD	CVTEST	CVPOS	CVORRES	CVORRESU
CV Test1	CV	Group A	2321	LVSYSBP	Left Ventricular Systolic Pressure	UNCONSTRAINED	118.37	mmHg
CV Test1	CV	Group A	2322	LVEDP	Left Ventricular End Diastolic Pressure	UNCONSTRAINED	5.5588	mmHg
CV Test1	CV	Group A	2323	HR	Heart Rate	UNCONSTRAINED	74.33	bpm
CV Test1	CV	Group A	2324	DPDTAVG	dP/dt Average	UNCONSTRAINED	3310.4	mmHg/s
CV Test1	CV	Group A	2325	SYSBP	Systolic Blood Pressure	UNCONSTRAINED	117.83	mmHg
CV Test1	CV	Group A	2326	DIABP	Diastolic Blood Pressure	UNCONSTRAINED	81.212	mmHg
CV Test1	CV	Group A	2327	MAP	Mean Arterial Pressure	UNCONSTRAINED	99.818	mmHg
CV Test1	CV	Group A	2328	PULSEPR	Pulse Pressure	UNCONSTRAINED	36.623	mmHg
CV Test1	CV	Group A	2329	DPDTMAX	Left Ventricular Maximum Positive dP/dt	UNCONSTRAINED	10566	mmHg/s
CV Test1	CV	Group A	2330	DPDTMIN	Left Ventricular Minimum Positive dP/dt	UNCONSTRAINED	2565.3	mmHg/s
CV Test1	CV	Group A	2331	LVSYSBP	Left Ventricular Systolic Pressure	UNCONSTRAINED	111.21	mmHg
CV Test1	CV	Group A	2332	LVEDP	Left Ventricular End Diastolic Pressure	UNCONSTRAINED	4.7925	mmHg
CV Test1	CV	Group A	2333	HR	Heart Rate	UNCONSTRAINED	60.651	bpm
CV Test1	CV	Group A	2334	DPDTAVG	dP/dt Average	UNCONSTRAINED	3023.2	mmHg/s
CV Test1	CV	Group A	2335	SYSBP	Systolic Blood Pressure	UNCONSTRAINED	112.43	mmHg
CV Test1	CV	Group A	2336	DIABP	Diastolic Blood Pressure	UNCONSTRAINED	76.463	mmHg
CV Test1	CV	Group A	2337	MAP	Mean Arterial Pressure	UNCONSTRAINED	94.816	mmHg
CV Test1	CV	Group A	2338	PULSEPR	Pulse Pressure	UNCONSTRAINED	35.969	mmHg
CV Test1	CV	Group A	2339	DPDTMAX	Left Ventricular Maximum Positive dP/dt	UNCONSTRAINED	3589.6	mmHg/s
CV Test1	CV	Group A	2340	DPDTMIN	Left Ventricular Minimum Positive dP/dt	UNCONSTRAINED	2594.9	mmHg/s
CV Test1	CV	Group A	2341	LVSYSBP	Left Ventricular Systolic Pressure	UNCONSTRAINED	115.27	mmHg
CV Test1	CV	Group A	2342	LVEDP	Left Ventricular End Diastolic Pressure	UNCONSTRAINED	5.6718	mmHg
CV Test1	CV	Group A	2343	HR	Heart Rate	UNCONSTRAINED	64.312	bpm
CV Test1	CV	Group A	2344	DPDTAVG	dP/dt Average	UNCONSTRAINED	3042.7	mmHg/s
CV Test1	CV	Group A	2345	SYSBP	Systolic Blood Pressure	UNCONSTRAINED	117.19	mmHg
CV Test1	CV	Group A	2346	DIABP	Diastolic Blood Pressure	UNCONSTRAINED	78.691	mmHg
CV Test1	CV	Group A	2347	MAP	Mean Arterial Pressure	UNCONSTRAINED	98.115	mmHg
CV Test1	CV	Group A	2348	PULSEPR	Pulse Pressure	UNCONSTRAINED	38.502	mmHg
CV Test1	CV	Group A	2349	DPDTMAX	Left Ventricular Maximum Positive dP/dt	UNCONSTRAINED	4766.2	mmHg/s

6235 records total

Figure 18. – SEND Viewer dialog.

Figure 19 shows the Options pin expanded. Four filters are available to tailor the SEND output to CSV format.

1 – Study

Lists previously collected studies that are present on the workstation selected. Selecting a study will load any SEND related information. Figure 19 shows CVTest1 selected and its associated data displayed.

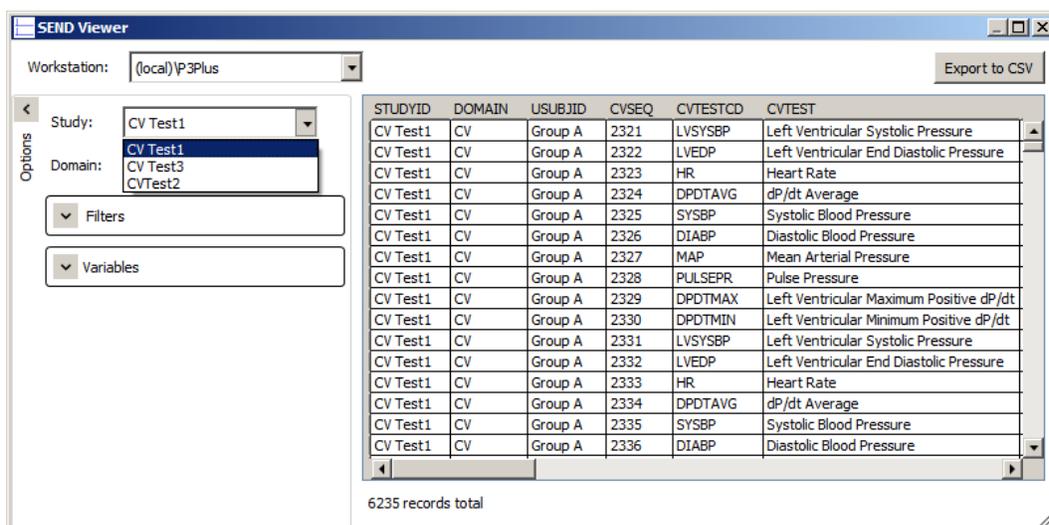


Figure 19. – SEND Viewer dialog.

2 – Domain

After the Study has been chosen, the correct SEND Domain must be selected. If an incorrect Domain is selected, the appropriate data will not populate in the window.

3 – Filters

Expanding the Filters option allows the user to define what information will be displayed and available for output to CSV. This includes USUBJID, NOMDAY, and TESTCD for the given Domain.

The “All” checkbox allows the user to toggle all variables on or off. Additionally, each variable may be selected or deselected individually to create the desired output.

It is important to note that not all Ponemah generated parameters are required by CDISC (TESTCD). Only parameters defined by CDISC will be displayed.

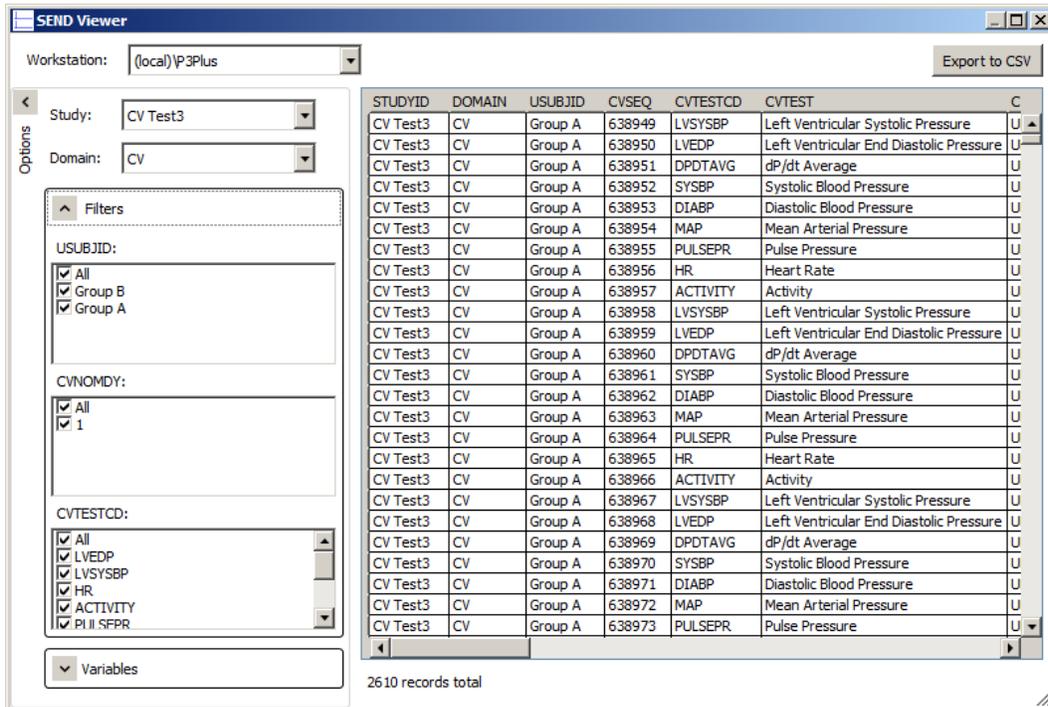


Figure 20. – SEND Viewer dialog displaying available filters.

4 – Variables

The Variables option controls which CDISC defined variables will be displayed and outputted to the CSV formatted file. Toggling specific variables will enable or disable the variables in the viewer and the associated CSV file when generated. This may be used to remove variables that are not populated with information.

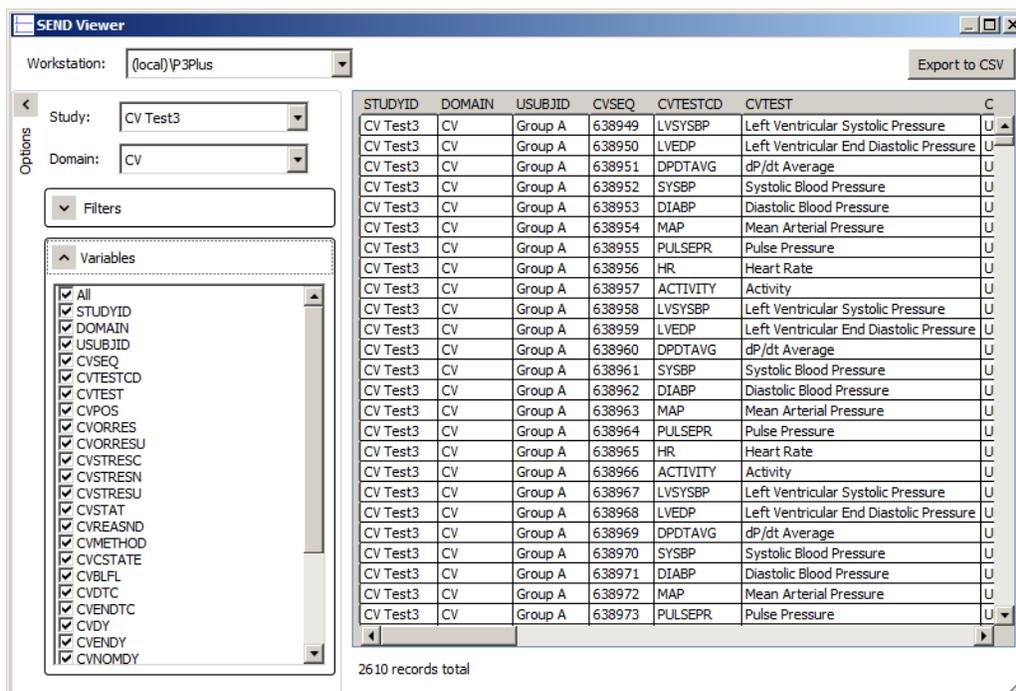


Figure 21. – SEND Viewer dialog with Variables option expanded.

In the example study above (CVTest3 from Figure 21), EG Domain data was collected along with CV Domain data. In order to get the EG Domain variables, simply select the EG Domain as shown in Figure 22 and update the remaining Options to provide the desired output.

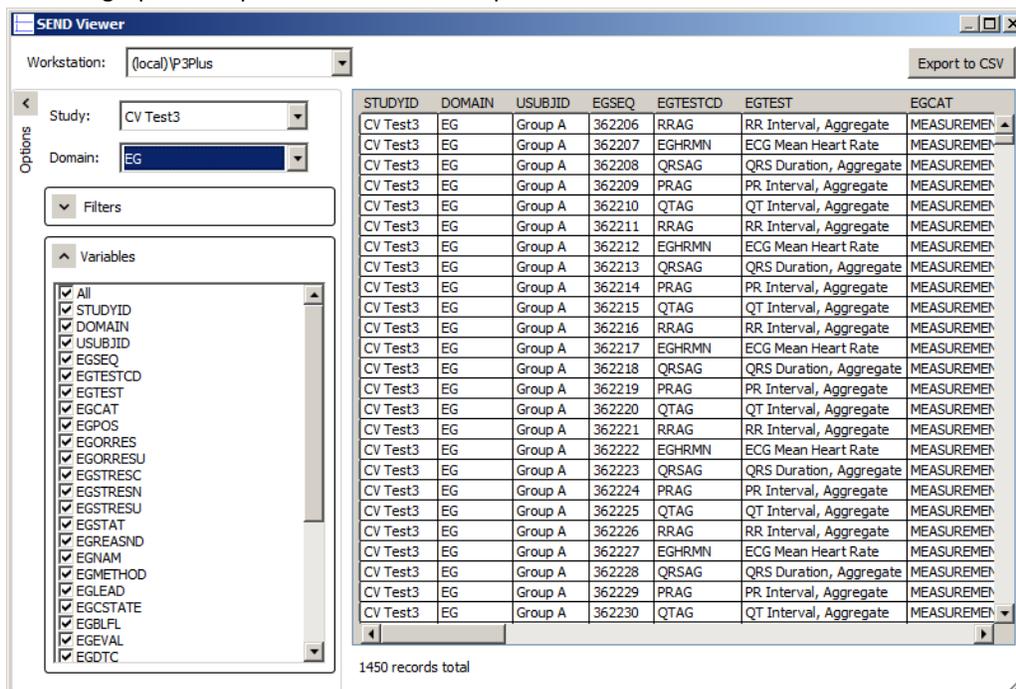


Figure 22. – SEND Viewer dialog with Variables option expanded and EG Domain selected.

SEND EXPORT TO CSV

Once the appropriate filters and variables have been defined as outlined in the SEND Viewer section, the data is ready to export to a CSV formatted file. This is accomplished by selecting the Export to CSV button in the top, right-hand portion of the dialog.

NOTE: It is important to note that this functionality is only available if the SEND SQL Server Option has been enabled in the license file. If only SEND Output is enabled, these features will not be available.

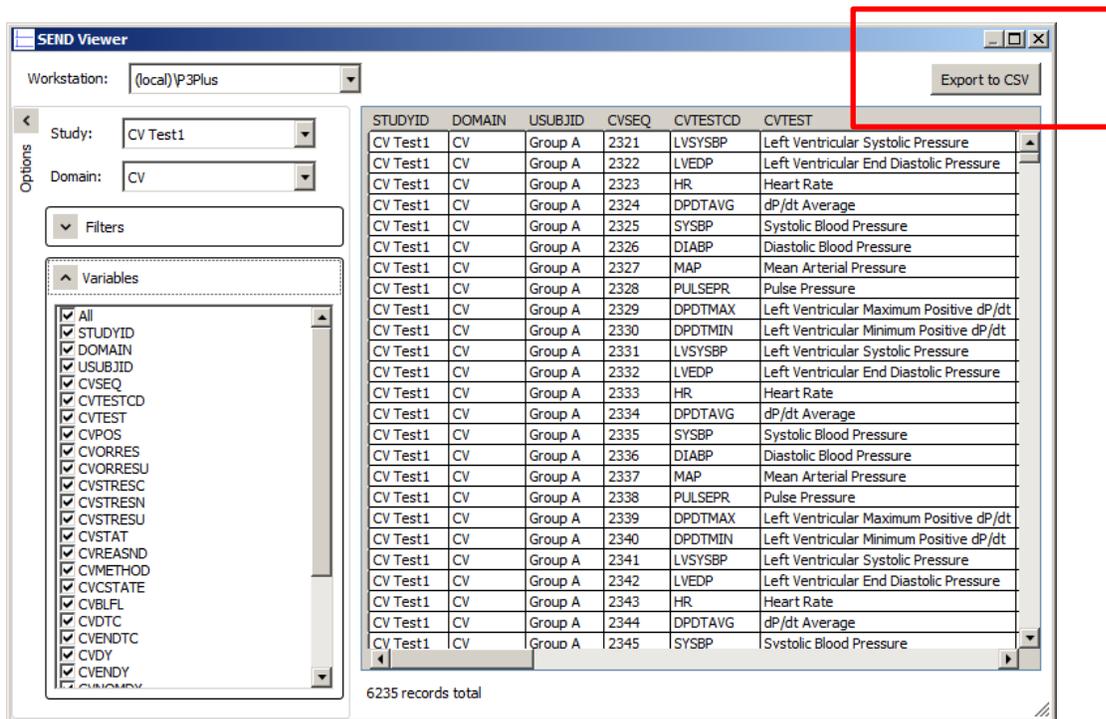


Figure 23. – SEND Viewer dialog with Export to CSV button highlighted.

A Windows dialog will be displayed with the default location set to the study folder within the Ponemah_Data directory. The file name will include the study name and date and will be configured to save in the CSV format. The name and location for saving the file may be changed by modifying the “File name” edit field and using the browse feature to specify a different location.

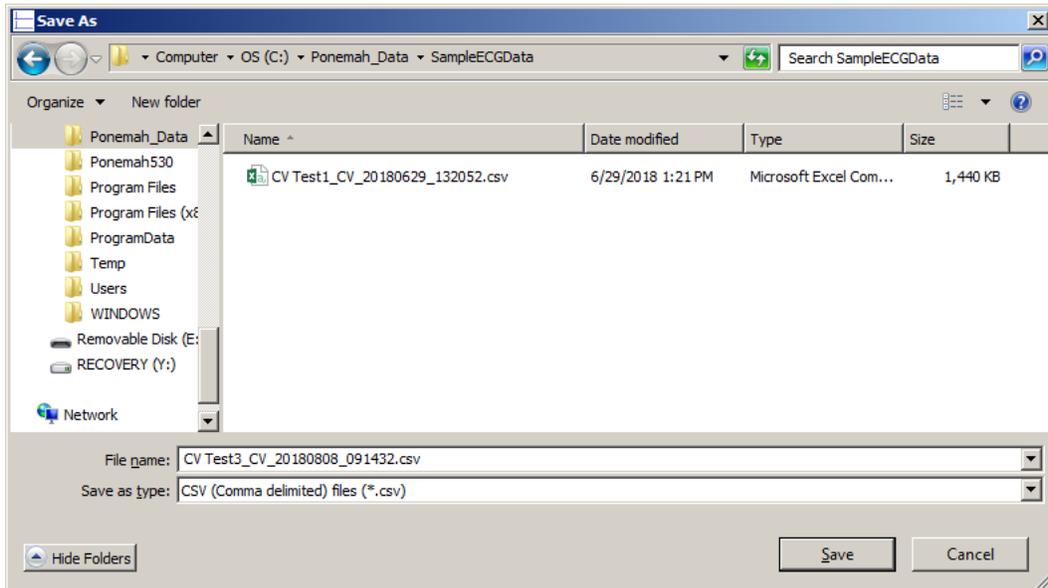


Figure 24. – Default location and file type for saving SEND output via SEND Viewer

Once the file has been saved, browsing to the location and opening the file will display the data as configured in SEND Viewer. Figure 25 shows the CSV output opened using Microsoft Excel.

STUDYID	DOMAIN	USUBJID	CVSEQ	CVTESTCD	CVTEST	CVPOS	CVORRES	CVORRES	CVSTRESC
CV Test1	CV	Group A	2321	LVSYSBP	Left Ventricular Systolic Pressure	UNCONSTRAINED	118.37	mmHg	118.37
CV Test1	CV	Group A	2322	LVEDP	Left Ventricular End Diastolic Pressure	UNCONSTRAINED	5.5588	mmHg	5.5588
CV Test1	CV	Group A	2323	HR	Heart Rate	UNCONSTRAINED	74.33	bpm	74.33
CV Test1	CV	Group A	2324	DPDTAVG	dP/dt Average	UNCONSTRAINED	3310.4	mmHg/s	3310.4
CV Test1	CV	Group A	2325	SYSBP	Systolic Blood Pressure	UNCONSTRAINED	117.83	mmHg	117.83
CV Test1	CV	Group A	2326	DIABP	Diastolic Blood Pressure	UNCONSTRAINED	81.212	mmHg	81.212
CV Test1	CV	Group A	2327	MAP	Mean Arterial Pressure	UNCONSTRAINED	99.818	mmHg	99.818
CV Test1	CV	Group A	2328	PULSEPR	Pulse Pressure	UNCONSTRAINED	36.623	mmHg	36.623
CV Test1	CV	Group A	2329	DPDTMAX	Left Ventricular Maximum Positive dP/dt	UNCONSTRAINED	10566	mmHg/s	10566
CV Test1	CV	Group A	2330	DPDTMIN	Left Ventricular Minimum Positive dP/dt	UNCONSTRAINED	2565.3	mmHg/s	2565.3
CV Test1	CV	Group A	2331	LVSYSBP	Left Ventricular Systolic Pressure	UNCONSTRAINED	111.21	mmHg	111.21
CV Test1	CV	Group A	2332	LVEDP	Left Ventricular End Diastolic Pressure	UNCONSTRAINED	4.7925	mmHg	4.7925
CV Test1	CV	Group A	2333	HR	Heart Rate	UNCONSTRAINED	60.651	bpm	60.651
CV Test1	CV	Group A	2334	DPDTAVG	dP/dt Average	UNCONSTRAINED	3023.2	mmHg/s	3023.2
CV Test1	CV	Group A	2335	CVSCP	Sustolic Blood Pressure	UNCONSTRAINED	112.43	mmHg	112.43

Figure 25. – Example of CSV output

SYNCHRONIZE SEND DATA


Send information can be synchronized to a dedicated server or workstation. There are two types of synchronization, Synchronize SEND Data and Synchronize SEND Study Data.

NOTE: It is important to note that this functionality is only available if the SEND SQL Server Option has been enabled in the license file. If only SEND Output is enabled, these features will not be available.

SYNCHRONIZE SEND DATA

The Synchronize SEND Data feature allows SEND data contained within SQL on a workstation to be synchronized with another workstation, or server, to create a single repository for SEND data. This requires configuration of the Output Server in the SEND setup dialog.

Enter the Server name or dedicated workstation name and use the Test Connection button to verify that the connection is valid. Along with the server name, the SQL Instance (P3Plus_v2) must also be defined, example “server name\P3Plus_v2”. If successful, a message will post stating that the connection to the SEND database was successful.

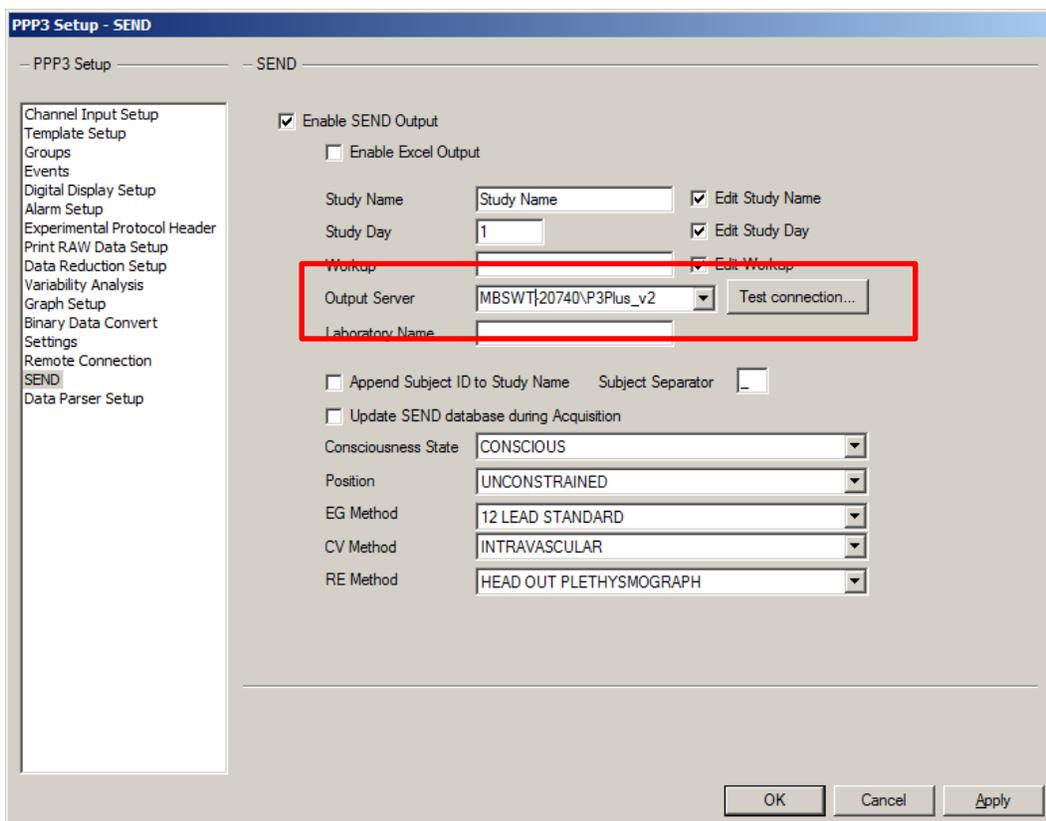

The screenshot shows the 'PPP3 Setup - SEND' dialog box. The 'SEND' tab is active. On the left is a tree view with 'SEND' selected. The main area contains the following settings:

- Enable SEND Output
- Enable Excel Output
- Study Name: Study Name Edit Study Name
- Study Day: 1 Edit Study Day
- Workshop: Edit Workshop
- Output Server: MBSWTJ20740\P3Plus_v2 - Laboratory Name:
- Append Subject ID to Study Name Subject Separator:
- Update SEND database during Acquisition
- Consciousness State: CONSCIOUS
- Position: UNCONSTRAINED
- EG Method: 12 LEAD STANDARD
- CV Method: INTRAVASCULAR
- RE Method: HEAD OUT PLETHYSMOGRAPH

Buttons at the bottom: OK, Cancel, Apply.

Figure 26. – Output Server configuration in SEND dialog

If no Output Server has been configured, attempting to synchronize study data will post a message requesting that the server be configured in P3 Setup.

When performing a synch, the Output Server will be locked out from other workstations until the current synch has been completed. If an attempt to synchronize to the Output Server is initiated by a second workstation, a message will be posted that the server (workstation) is in use and attempts to synchronize the data should be retried at a later time. In addition, error checking is performed during the synchronization. If conflicts between the data on the Output Server and the workstation performing the synch, conflict resolution dialogs will be presented to allow the appropriate information to be updated.

After configuring the Output Server, navigate to the Synchronize SEND Data as shown in Figure 27. The SEND data for the currently loaded Review file will be automatically synchronized to the SEND database on the Output Server (Synchronize SEND Study Data is not available within a Review session). Only data from the current Review session will be synchronized to the Output Server when in Review.

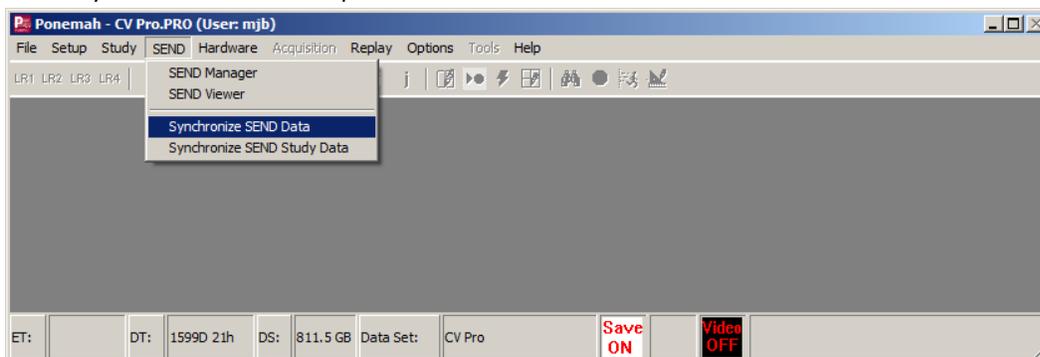


Figure 27. – Synchronize SEND Data dialog selection from SEND pull-down menu

The Synchronize Study Data function can also be used when outside of a Review session. Figure 28 shows the pop-up dialog that allows selection of the Study and Output Server. Clicking the OK button will synchronize the selected study to the designated Output Server defined. This will include all SEND data associated with that study, whereas when in a Review session, only the data for the currently loaded Review file will be synchronized.

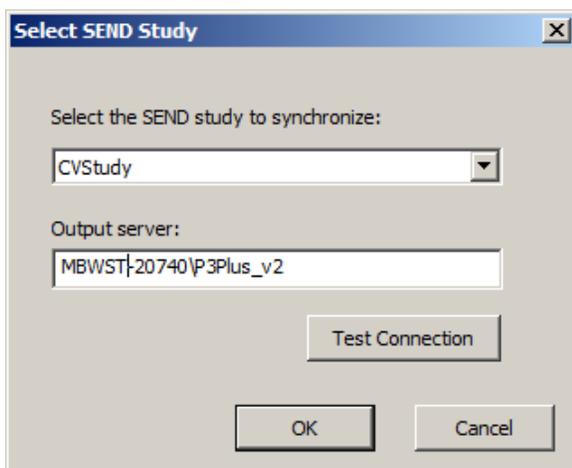


Figure 28. – Synchronize SEND Data dialog outside of Review session

SYNCHRONIZE SEND STUDY DATA

NOTE: It is important to note that this functionality is only available if the SEND SQL Server Option has been enabled in the license file. If only SEND Output is enabled, this feature will not be available.

The Synchronize SEND Study Data will synchronize from all workstations connected and visible on the network to the workstation performing the synch and to the Output Server. Both the workstation performing the synch and the Output server will then contain the same SEND information.

The Output Server defined in the SEND dialog must be part of the Study configuration, otherwise a message will post stating that they sync cannot be performed to the Output Server.

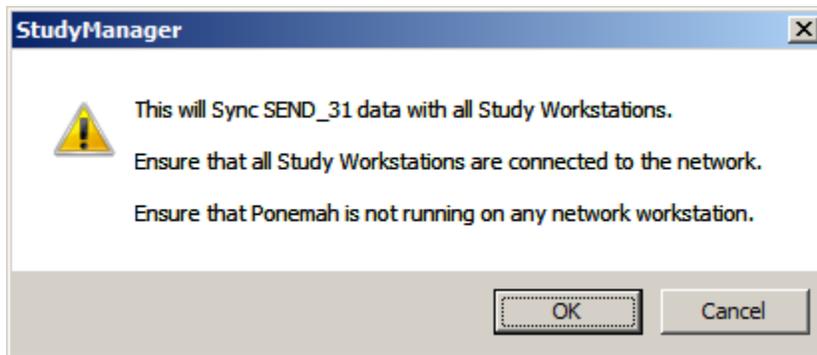


Figure 29. – Synchronize SEND Study Data notification.

The Synchronize SEND Study Data function is only available in idle mode and will not be available while in a Review session.

HIGH LEVEL PROCESS FLOW FOR GENERATING SEND OUTPUT

The previous sections have described how to configure and generate SEND output. This section will provide two scenarios and the high-level process flow needed to generate SEND output. No detailed configuration information is provided below, refer to previous sections for additional information.

While the following two examples provides guidance on creating SEND output, which includes configuration information, once a SEND configuration has been completed and saved in the protocol, only saving of the derived data is necessary to generate SEND output.

Generating SEND Output to Excel (outside of the Study Protocol Option)

- Ensure SEND Output or SEND SQL Server Output option is enabled in the license file
- Open a Review file (.RVW)
- Configure Data Reduction, if not previously configured, with CDISC defined variables
 - Ensure the event to trigger Data Reduction has been initiated

- Update Groups tab with appropriate Name and Dose information for each subject
- Enable the SEND dialog under P3 Setup and update the appropriate fields
 - Select the checkbox for Enable Excel Output
- Perform a Save Marks Section or Saved Derived Data to generate the Data Reduction output to SEND
- Navigate to the Ponemah_Data directory or the location where the Excel file was saved
 - Open Excel workbook and verify that the appropriate worksheets exist for the Domains used

Generating SEND Output to CSV using the Study Protocol Option


- Ensure SEND SQL Server Output option is enabled in the license file in order to generate CSV files
- Select the desired Study from the Manage Studies dialog
- Enter the Run Study dialog and select the desired Run from the Previous Runs list
- Data Reduction should have been previously configured and saved as part of the Study configuration (protocol file) with CDISC defined variables. Modification of Data Reduction is not permitted after data collection within the Study Protocol Option
 - Ensure the event to trigger Data Reduction has been initiated
- Groups should have been configured with the appropriate Name and Dose information for each subject as part of the Study configuration.
 - Select Use Study Subjects and Use Study Doses to automatically pull information from the Study configuration.
 - Uncheck Use Study Subjects and Use Study Doses if edits are necessary to correct entry errors from Study configuration
- Enable the SEND dialog under P3 setup and updated the appropriate fields if not previously saved in the protocol
 - Uncheck the following to allow previously defined Study information to be pulled into SEND automatically
 - Edit Study Name
 - Edit Study Day
 - Edit Study Workup
 - If desired, uncheck these edit boxes to allow for manual edits to better define the output being generated or correct for entry errors when creating the Study
- Perform a Save Run (Marks Section) to generate the Data Reduction Output to SEND
- Navigate to the Send Viewer and select the desired Study to generate a CSV file output
 - Select the appropriate Domain, Filters, and Variables
 - Use the Export to CSV button to generate the SEND output file
- Optionally, use the Synchronize SEND Data and Synchronize SEND Study Data to synchronize data across workstations or servers

TROUBLESHOOTING

This section outlines common messages encountered with incorrect settings when SEND is enabled. Each item will list the issue or message that may be posted along with a resolution.

MESSAGES POSTED DUE TO INCORRECT SEND CONFIGURATION

1 – Incorrect ECG labels

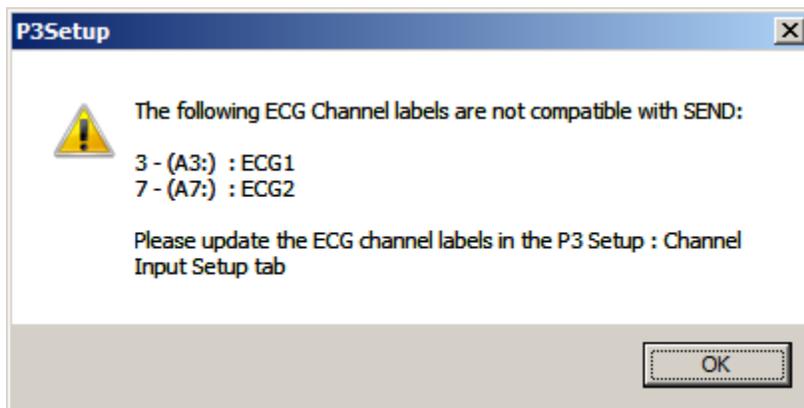

Issue: CDISC dictates specific labeling terminology for ECG lead presentations. If the label entered in the Channel Input Setup menu does not match that used by CDISC, the above message will be posted.

Resolution: Enter the Channel Input Setup menu and update the labeling for the ECG channels identified in the message posted. When you select the Label column, you will have a drop-down list that allows you to select the correct lead presentation.

2 – Duplicate TESTCDs



Issue: A CDISC TESTCD can only be reported once for a subject in a given interval in a single domain. The above error will be displayed if duplicate derived parameters exist across multiple analysis algorithms. In the example above, the HR (Heart Rate) parameter is enabled in both the Blood Pressure and Left Ventricular Pressure analysis algorithms.

Resolution: Enter the Channel Input Setup menu and update the derived parameters list for the Subject(s) listed in the warning message. Resave the file to generate the desired output. If you say OK to the message, the software automatically uses the TESTCD (in this case HR) from the first channel where the conflict was identified. In this case, the hardware configuration had LV pressure as channel 1 and BP (systemic blood pressure) defined as channel 2 for both Group A and Group B. The HR parameter was reported for the LV channel and removed from the BP channel.

3– NO SEND data in Excel or SEND Viewer

Issue: Data not present in the Excel file or in SEND Viewer after performing a save of the Marks Section/Run or Derived Data.

Resolution: Several reasons may exist which will result in no data available in the output file.


- SEND has not been enabled in P3 Setup
 - If using SEND to Excel, ensure checkbox in SEND dialog has been enabled
- Data Reduction has not been configured
- CDISC TESTCDs are not enabled in Data Reduction. Not all derived parameters (Data Reduction) are currently defined by CDISC
- Ensure the Event to trigger Data Reduction has been placed
- If SEND does not appear in your P3 Setup menu, the current Ponemah license file likely does not have the SEND Output or SEND SQL Server Output option enabled. This information can be found under the View

License File dialog from the Help pull-down menu. These features will be listed under the Options box. The SEND pull-down menu while in idle mode is only available (accessible) if the SEND SQL Server Output option is enabled. Contact Technical Support for assistance.

4 – No Output Sever Defined



Or

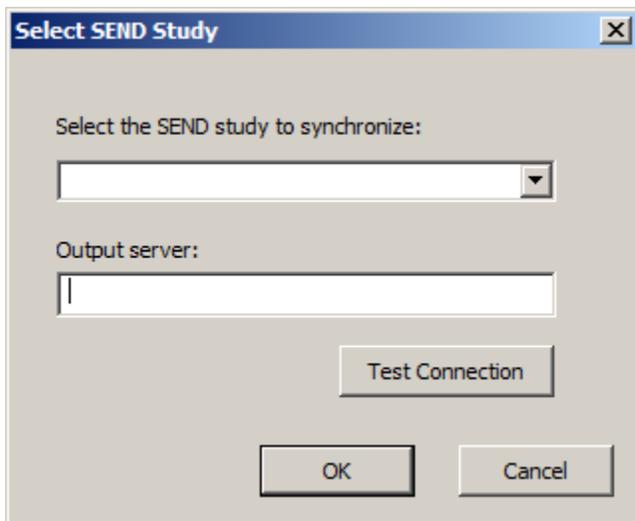

Issue: No Output Sever has been configured


Resolution: The first message will post when in an active Review session and the second message when in idle mode. Enter the appropriate server/workstation in the SEND dialog under P3 Setup. Ensure that the Ponemah Instance has been included with the server/workstation name.

5 – No subject ID for USUBJID TESTCD

Issue: No user defined subject ID present in the SEND output for USUBJID.

Resolution: When not using the Study Protocol Option and pulling in Subject information automatically, the default label (Group A, Group B, etc.) will be placed in the SEND output. From P3 Setup, enter the Groups dialog and modify the Name column. Perform a Save Derived Data or Save Marks Section to update.

6 – No Data in SEND Viewer

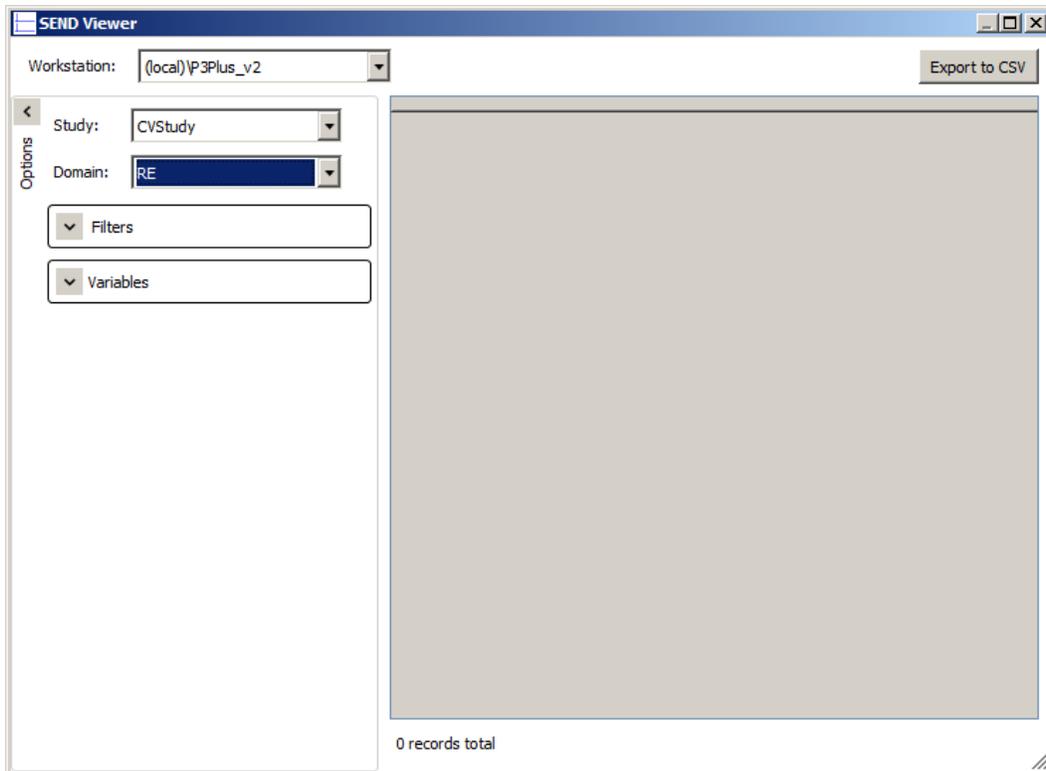

Issue: No data is present when opening the SEND Viewer dialog for a given study when all information has been configured properly (SEND, Data Reduction, etc.) and a save of the data has been performed.

Resolution: Ensure that the correct Domain is listed for the study that has been selected. Data currently exists for the CVStudy example shown above. However, no data in the RE Domain exists. Therefore, no data will populate the window. Select the correct Domain, EG and CV in this example, and the SEND information will be displayed.

TECHNICAL SUPPORT

DSI™ is available to help you with your questions and concerns. Should you hit a road block or need some additional training, please feel free to contact us. We are happy to help!

DSI TECHNICAL SUPPORT—NORTH AMERICA

Email: Support@datasci.com

Toll-free in U.S. and Canada

Phone: 1-800-262-9687

Monday through Friday: 8 AM to 5 PM CST
(except Holidays)

DSI TECHNICAL SUPPORT—EUROPE

Email: Europe-support@datasci.com

Phone: +44 1359 259400

Monday through Friday: 8 AM to 5 PM CET

DSI TECHNICAL SUPPORT—ALL OTHER COUNTRIES

Phone: +1-651-481-7400

APPENDIX 1

Not all derived parameters calculated by Ponemah are currently defined by CDISC. Below lists the currently accepted parameters by general signal type. Please note that derived parameters may exist in multiple Analysis algorithms. For example, the derived parameter “Sys” is reported by both the Blood Pressure (BP) and Left Ventricular Pressure (LVP) analysis algorithms.

In addition to enabling only the CDISC defined parameters, Data Reduction calculations selected must also be defined by CDISC or no SEND output will be generated. The accepted Data Reductions calculations are Average (AVG), Maximum (Max), and Minimum (Min).

ECG

- Mean Heart Rate
- P-H
- PWdth
- PP-I
- PR-I
- QRS
- QT-I
- QTcb
- QTcf
- QTcv
- R-H
- RR-I
- ST-I
- ST-E
- T-H
- T-A

- Tpe-I

Activity

- A_TA

Pressure

- Dia
- +dP/dt
- -dP/dt
- HR
- LVEDP
- Sys
- Mean
- PH
- Q-A
- TPR

Respiratory

- TV
- Res
- RT
- ET
- IT
- MV
- Penh
- Max (PCR analysis module)
- PIF
- Min
- Cdyn
- PEF
- BPM
- RBpm

Temperature

- Mean