



SEND 3.1: Standard for Exchange of Nonclinical Data



Data Sciences International

**119 14TH STREET NW
SUITE 100**

ST. PAUL, MN 55112 USA

(651)481-7400 • 1(800)262-9687

www.datasci.com

PN: MU00285-001 REV02

OVERVIEW

This manual highlights how DSI's SEND 3.1 can be used to produce output needed for data submission to the FDA. This document will provide an overview of the system components and basic operation.

Table of Contents

Welcome	3
Background	3
SEND Setup from PPP3 Setup	4
SEND dialog	4
Groups	7
Data Reduction	9
Using Study Protocol Option and SEND	15
SEND Configuration When Using Study	15
SEND Manager	18
SEND Viewer and Output	22
SEND Export To Excel	22
SEND Viewer	23
Send Export to CSV	28
SEND Log	29
Controlled Terminology	32
Synchronize SEND Data	35
Send Remote Server Synchronization Setup	35
Synchronize SEND Data	38
Synchronize SEND Study Data	39
High Level Process Flow for Generating SEND Output	41
Troubleshooting	43
Messages Posted Due to Incorrect SEND Configuration	43

Technical Support.....	47
Appendix 1	47
Appendix 2	49

WELCOME

The structure of this manual was designed to sequentially guide you through setup of the various SEND components and ultimately generating the standardized output needed for data submission. This manual references functionality that may not be fully detailed here. It may be necessary to reference additional Ponemah manuals for detailed feature information. This information can be found under the Help menu within Ponemah.

WHAT YOU WILL BE LEARNING

1. Determine what functions are needed in order to generate SEND information
2. Understand how to configure SEND specific features
3. Generate SEND output

BACKGROUND

SEND (Standard for Exchange of Nonclinical Data) is a standardized method to exchange data between organizations and for data submission to the FDA or other regulatory bodies for non-clinical data. Detailed information on formatting, standards, terminology and other information can be found at <https://www.cdisc.org/>.

In order to produce output using the SEND module, users must ensure the following:

- SEND Output or SEND SQL Server Output has been enabled in license file
- Derived Parameters selected
 - Not all derived parameters generated by Ponemah are currently defined by CDISC. Derived parameters will only be present in SEND output if those parameters defined by CDISC have been enabled in the Channel Input Setup menu.
 - See Appendix 1 for a list of currently accepted parameters
 - Data reduction enabled for subjects.
 - The currently defined calculations are Average (Avg), Maximum (Max), and Minimum (Min)
- The Controlled Terminology dialog has the desired mappings configured
 - The Controlled Terminology dialog provides a mapping of the terminology between the Ponemah derived nomenclature and the expected SEND controlled terminology. This dialog can be modified by the user to include additional Ponemah parameters. Additionally, as parameters are updated by CDISC, this dialog can be easily modified to include the new information.

This manual assumes users have read and understand the core components of the Ponemah software application. In some instances, SEND is built upon existing menus or functions in the software. If additional information is

needed on functionality outside of the SEND application, please review the manuals located under the Help pull-down menu in the Ponemah application.

SEND SETUP FROM PPP3 SETUP

In order to create SEND output, users must have enabled Data Reduction, selected derived parameters in the Channel Input Setup menu that correspond to output designated by CDISC, and performed a Save of the data while in Review (Save Marks Section, Save Derived Data, or Save Run).

SEND output can be generated within our outside of the Study Protocol Option (Study). This section will describe how to set up components that are used both with and without Study. Specific use with Study will be described in a later section.

SEND DIALOG

SEND can be configured in idle mode and saved as part of the protocol or modified after collection in Review. To modify SEND settings, select P3 Setup from the Setup pull-down menu and click on SEND in the tree-view menu on the left-hand side of the dialog. By default, SEND will be disabled. In order to edit, select “Enable SEND Output”. Once enabled, fields will become editable.

Definitions for the functions displayed in the SEND dialog are listed below.

PPP3 Setup - SEND

PPP3 Setup - SEND

Channel Input Setup
Template Setup
Groups
Events
Experimental Protocol Header
Data Reduction Setup
Variability Analysis
Graph Setup
Settings
SEND

Enable SEND Output

Enable Excel Output

Study Name: Study Name Edit Study Name

Study Day: 1 Edit Study Day

Workup: Edit Workup

Output Server: Test connection...

Laboratory Name:

Append Subject ID to Study Name Subject Separator:

Update SEND database during Acquisition

Consciousness State: CONSCIOUS

Position: UNCONSTRAINED

EG Method: 12 LEAD STANDARD

CV Method: INTRAVASCULAR

RE Method: HEAD OUT PLETHYSMOGRAPH

OK Cancel Apply

Figure 1. – SEND dialog from the PPP3 Setup menu.

1 – Enable SEND Output

Toggles the ability to modify the editable fields displayed in Figure 1 and controls whether or not SEND output is turned on or off.

2 – Enable Excel Output

SEND output can be ported to Microsoft Excel along with the derived, log, and subject information currently created when saving a Marks section, Derived Data, or a Run (within Study). If enabled, a new worksheet for each SEND domain will be created within the Excel workbook that contains SEND information along with the expected Derived, Data Reduction, and Experimental Log information. Refer to section SEND Export To Excel for details on the output generated when enabling this feature.

If this option is not checked, SEND output can be generated by creating a CSV file using SEND Viewer (see SEND Viewer below).

3 – Study Name

Allows the user to create a unique study name for the SEND output. If using the Study Protocol Option, this information will be automatically pulled by the software. However, the option to modify the Study name is possible by checking the box next to the edit field. This may be needed to further define a set of data or update an incorrect Study name that may have been created.

4 – Study Day


The day of collection may be entered here. Note that CDISC does not recognize Day 0. If using the first day of the study collection, a “1” should be entered. If using the Study Protocol Option, this information will be automatically pulled by the software. However, the option to modify the Study Day is possible by checking the box next to the edit field. This may be needed to further define a set of data or update an incorrect Study entry.

5 – Workups

Workups can be manually defined, or if using the Study option, this information can be automatically pulled by the software. If pulled from Study, users can modify the workup name used in Study by checking the box next to the edit field.

6 – Output Server

Permits defining a dedicated server or workstation as the primary repository for SEND data. Additionally, the SEND instance name, as it is installed with the application, must be appended to the server/workstation name. Example naming convention would be “server name\P3Plus_V2”. P3Plus_V2 is the SQL Instance used by Ponemah for SEND related data.



NOTE: It is important to note that this functionality is only available if the SEND SQL Server Option has been enabled in the license file. If only SEND Output is enabled, this feature will not be available.

7 – Laboratory Name

Allows a unique identifier to be manually entered which will populate the EGNAM variable in the SEND output for the EG Domain

8 – Append Study ID to Subject Name

Will append the Subject ID to the Study Name in the SEND output. By default, the separator is “_” but can be changed. This modifies USUBJID in the SEND output for the domain.

9 – Update SEND database during acquisition

Generation of SEND data will typically be generated during Review after post processing the data. However, there may be instances where post analysis of the data is not needed such as in short duration, Tox applications. Enabling this checkbox will allow data generated during acquisition to be used by SEND.

10 – Consciousness State

Pull-down field allows selection of CDISC controlled terminology that represents the state of the animal, such as conscious or unconscious. These definitions are managed in the Controlled Terminology dialog.

11 – Position

Presents a pull-down menu allowing the selection of CDISC controlled terminology that represents the subjects position from unconstrained to various restrained states. These definitions are managed in the Controlled Terminology dialog.

12 – EG Method

Allows the selection of the CDISC controlled terminology that represents the ECG lead presentation used for collection. These definitions are managed in the Controlled Terminology dialog.

13 – CV Method

Allows the selection of CDISC controlled terminology for definition of pressure measurements. These definitions are managed in the Controlled Terminology dialog.

14 – RE Method

Permits selection of CDISC controlled terminology such as Head Out Plethysmograph, Whole Body Plethysmograph, or Invasive to describe the respiratory system being used. These definitions are managed in the Controlled Terminology dialog.

GROUPS

The Groups dialog allows the user to control the Group name and dose which is associated with a given Subject. In order to edit the Study Subject Name or Study Dose, click in the edit field under the appropriate column. Figure 2 shows subjects one and two (Group A and Group B) with the names and doses modified.

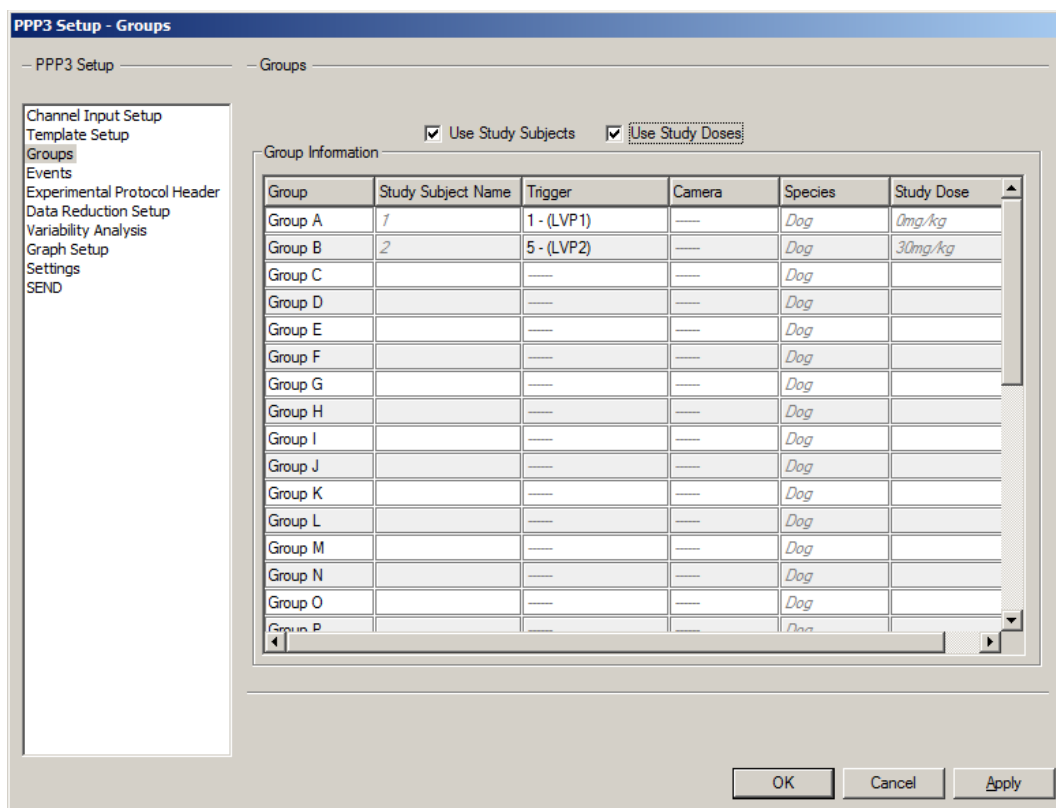


Figure 2. – Groups dialog from the PPP3 Setup menu.

Providing a unique name in the Groups dialog will propagate the subject information to SEND. Below shows example SEND output where Group A Study Subject Name was changed in the Groups tab (Figure 2). This identifier, ID 123, is shown under USUBJID in Figure 3.

Two check boxes are available at the top of the Groups dialog, Use Study Subjects and Use Study Doses. These two features are related to data generated using the Study Protocol Option. If enabled, Subject and Dose information entered in Study will be automatically pulled into the Groups tab. When these features are enabled, manually editing Study Subject Name and Study Dose will not be permitted. However, it may be necessary to update or correct information that was entered incorrectly into the Study Protocol Option. If needed, uncheck these boxes and enter the appropriate information manually. When these boxes are unchecked, the column headers revert back to the default values of Name and Dose. If not using the Study Protocol Option, these features do not need to be enabled.

1 – Use Study Subjects

Subject information defined within the Study Protocol Option can be automatically pulled into SEND output by selecting this checkbox

2 – Use Study Doses

Dose information entered within the Study Protocol Option will be pulled into SEND by selecting this checkbox

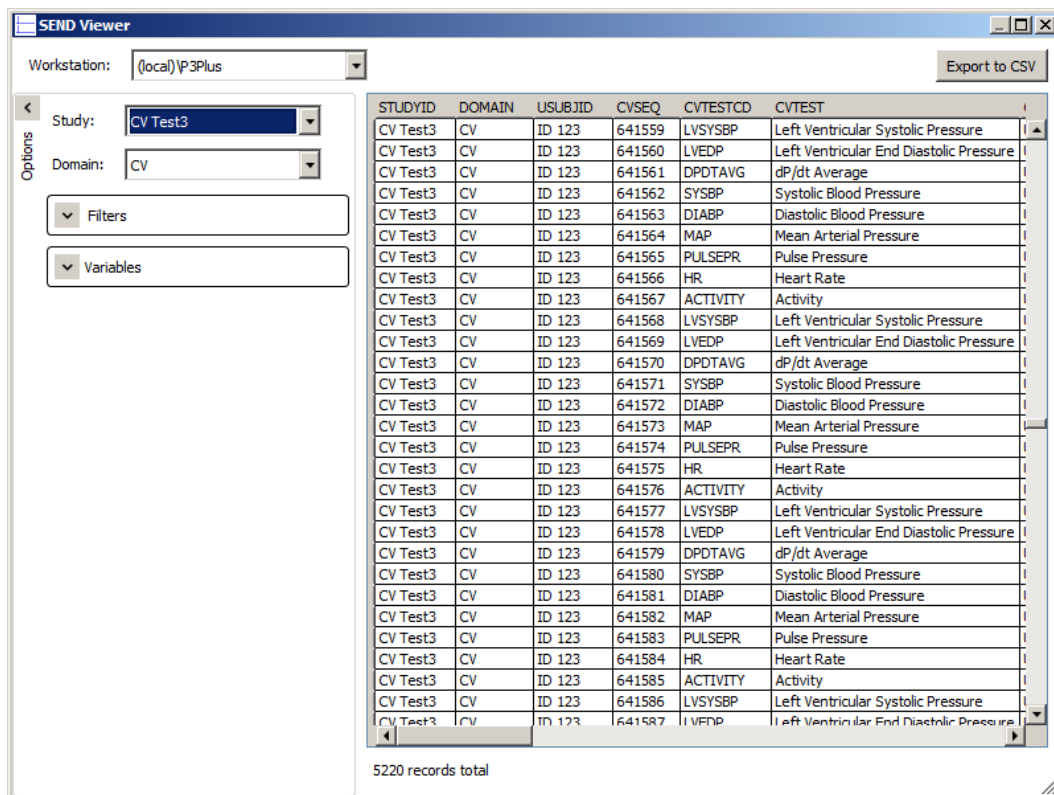


Figure 3. – Example SEND Viewer dialog showing Group information pulled into SEND (USUBJID).

DATA REDUCTION

Data Reduction must be enabled for all desired subjects in order to generate SEND output. Only parameters defined by CDISC will be available in the SEND output.

With SEND enabled, users can no longer create user specified labels as that will be handled automatically to conform to SEND requirements. When SEND has been enabled, Data Reduction will display “SEND compatibility enabled” at the top of the dialog to notify users that they will be complying with SEND requirements and some functions will not be able to be modified.

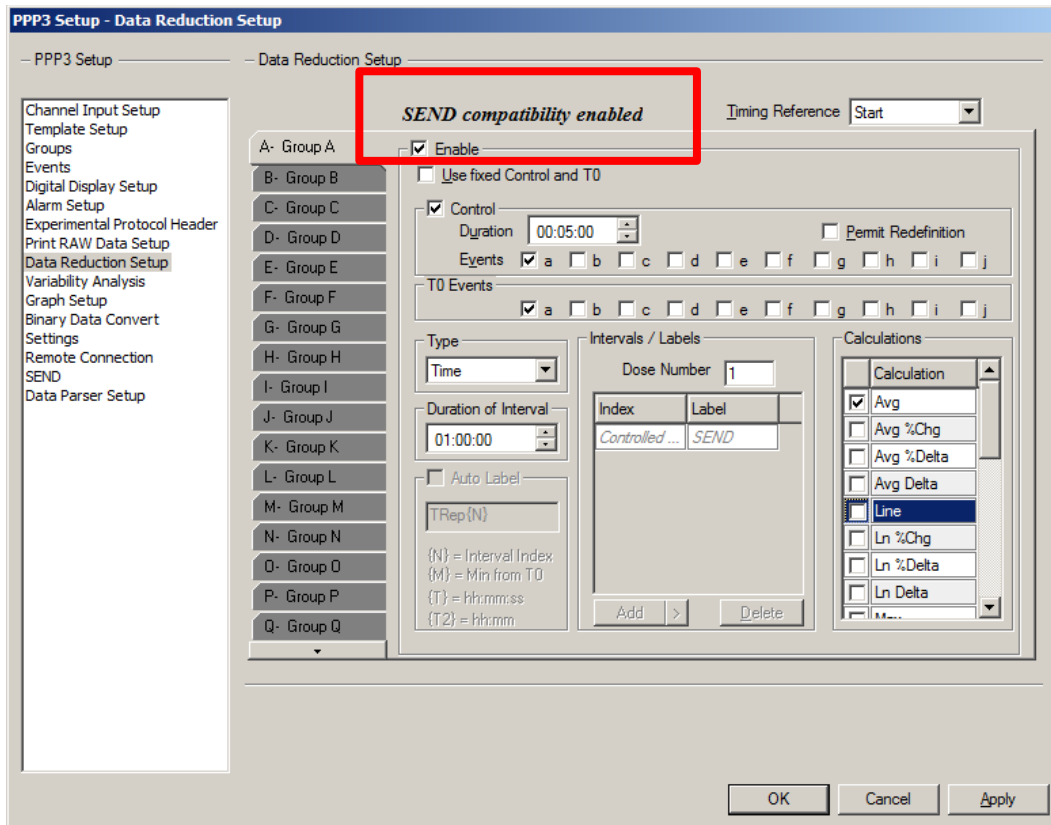


Figure 4. – Data Reduction dialog showing SEND enabled.

Several features have been added to the Data Reduction dialog to provide details on the data generated for SEND output.

1 – Timing Reference

The Timing Reference selection determines the point on each data reduction interval that will be used to report timing information for the interval (Figure 5). This is used to define Data Reduction labels along with TPT and ELTM.

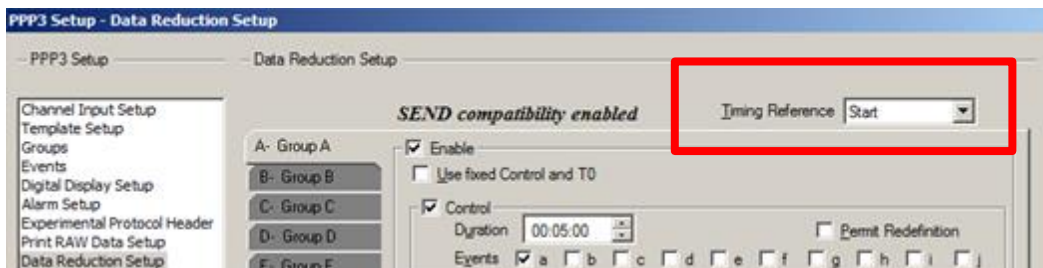


Figure 5. – Data Reduction showing Timing Reference selection.

Three selections are available which include Start, Middle, and End. The selection chosen will update the Data Reduction label to display the time information at the start of the logging interval, the middle of the interval, or the end of the interval as shown in Figures 6 and 7.

Figure 6 shows labels starting at 0s and incrementing by 10 minutes (10m, 20m, 30m) as dictated by the “Duration of Interval” chosen by the user.

Label	Start Time	End Time	Duration	1: Sys	1: LVEDP	1: +dP/dt	2: Sys	2: Dia	2: Mean	2: PH	2: HR
T0	00:00:00										
0s Postdose	00:00:00	00:10:00	0000:10:00	118.40	5.5708	3310.2	117.87	81.248	99.849	36.621	74.423
Avg											
10m Postdose	00:10:00	00:20:00	0000:10:00	111.20	4.7871	3023.4	112.41	76.446	94.795	35.967	60.68
Avg											
20m Postdose	00:20:00	00:30:00	0000:10:00	115.28	5.6705	3043.0	117.20	78.698	98.12	38.50	64.35
Avg											
30m Postdose	00:30:00	00:40:00	0000:10:00	117.80	6.0854	3062.5	119.69	81.78	100.54	37.906	69.436
Avg											
40m Postdose	00:40:00	00:50:00	0000:10:00	127.26	7.4468	3525.0	128.53	89.39	109.08	39.142	86.194
Avg											
50m Postdose	00:50:00	01:00:00	0000:10:00	121.04	6.2823	3302.2	121.49	82.506	102.14	38.985	74.726
Avg											
1h Postdose	01:00:00	01:10:00	0000:10:00	143.49	12.489	4083.5	142.69	98.744	120.67	43.951	106.27
Avg											
1h 10m Postd...	01:10:00	01:20:00	0000:10:00	119.46	7.1452	3384.1	120.36	83.135	102.49	37.225	76.565
Avg											
1h 20m Postd...	01:20:00	01:30:00	0000:10:00								

Figure 6. – Data Reduction label showing “Start” selected for the Timing Reference.

Figure 7 shows the same Start Time, End Time, Duration and averages for the data, but the label has been modified to display the “Middle” selection. Since the duration of the averaging interval is 10 minutes, selecting Middle will update the label information to specify the middle of that logging interval. Figure 7 shows those updated labels starting at 5m and incrementing every 10 minutes (5m, 15m, 25m, etc.).

Ponemah - SampleECGData.RVW (User: mjb) - [Data Reduction: a]

Functions Setup SEND Data Parser Options Window

Label	Start Time	End Time	Duration	1: Sys	1: LVEDP	1:+dP/dt	2: Sys	2: Dia	2: Mean	2: PH	2: HR
T0	00:00:00	00:00:00	0000:10:00								
5m Postdose Avg	00:00:00	00:10:00	0000:10:00	118.40	5.5708	3310.2	117.87	81.248	99.849	36.621	74.423
15m Postdose Avg	00:10:00	00:20:00	0000:10:00	111.20	4.7871	3023.4	112.41	76.446	94.795	35.967	60.68
25m Postdose Avg	00:20:00	00:30:00	0000:10:00	115.28	5.6705	3043.0	117.20	78.698	98.12	38.50	64.35
35m Postdose Avg	00:30:00	00:40:00	0000:10:00	117.80	6.0854	3062.5	119.69	81.78	100.54	37.906	69.436
45m Postdose Avg	00:40:00	00:50:00	0000:10:00	127.26	7.4468	3525.0	128.53	89.39	109.08	39.142	86.194
55m Postdose Avg	00:50:00	01:00:00	0000:10:00	121.04	6.2823	3302.2	121.49	82.506	102.14	38.985	74.726
1h 5m Postdose Avg	01:00:00	01:10:00	0000:10:00	143.49	12.489	4083.5	142.69	98.744	120.67	43.951	106.27
1h 15m Postd... Avg	01:10:00	01:20:00	0000:10:00	119.46	7.1452	3384.1	120.36	83.135	102.49	37.225	76.565
1h 25m Postd... Avg	01:20:00	01:30:00	0000:10:00								

Current Group Name Change Reference Group Name Change Raw File C:\Ponemah_Data\SampleECGData\SampleECGDat

Figure 7. – Data Reduction label showing “Middle” selected for the Timing Reference.

2 – Dose Number

Dose Number is manually edited and refers to the dose that will be referenced in the SEND output. The dose number listed here will be placed in the SEND output under TPT. By default, the value is “1” but can be modified to represent the desired dose if multiple dosing events occur in a single day. If SEND is not enabled in the SEND dialog, this field will not be editable.

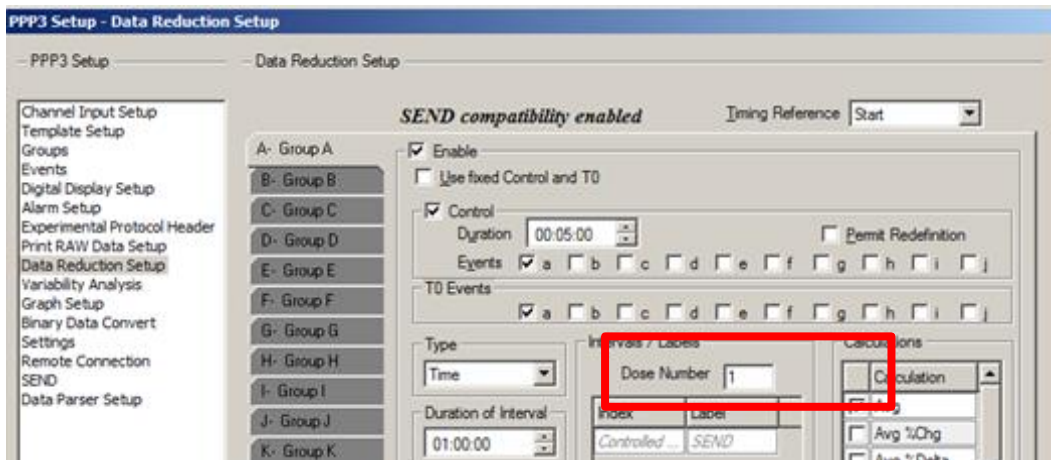


Figure 8. – Data Reduction dialog showing Dose Number edit field.

3 – Use Fixed Control and T0

The Use Fixed Control and T0 function allows the designation of the start of the Data Reduction without the use of an event. Entering in date and time into the T0 Time edit fields will trigger the start of Data Reduction. When using Time as the Data Reduction “Type”, all labels will be controlled by Ponemah for SEND compliance.

A time of 9:00AM was entered into the T0 Time field. In the list view dialog, the “0s Postdose” label is entered and controlled by Ponemah and the first logged line of data is at 9:00AM. All subsequent intervals are based on the Duration of Interval configured.

The image shows two screenshots. The top screenshot is the 'PPP3 Setup - Data Reduction Setup' dialog box. It features a left-hand menu with options like 'Channel Input Setup', 'Template Setup', 'Groups', 'Events', 'Experimental Protocol Header', 'Data Reduction Setup', 'Variability Analysis', 'Graph Setup', 'Settings', and 'SEND'. The main area is titled 'SEND compatibility enabled' and 'Timing Reference Start'. It includes a list of groups (A-Q) on the left. The 'Enable' checkbox is checked, along with 'Use fixed Control and T0'. The 'Control' section has a 'Duration' of 00:05:00 and a 'Control Time' of 5/17/2007 9:00:00 AM. The 'T0' section has a 'T0 Time' of 5/17/2007 9:00:00 AM. The 'Type' is set to 'Time' and the 'Duration of Interval' is 00:10:00. The 'Intervals / Labels' table shows 'Controlled ...' and 'SEND'. The 'Calculations' list includes 'Avg', 'Avg %Chg', 'Avg %Delta', 'Avg Delta', 'Line', 'Ln %Chg', 'Ln %Delta', and 'Ln Delta'. The bottom screenshot shows the 'Ponemah - CV Pro.RVW' software interface with a 'Data Reduction: a' table. The table has columns for Label, Start Time, End Time, Duration, and various data points (1: Sys, 1: LVEDP, 1: HR, 1: +dP/dt, 2: Sys, 2: Dia, 2: Mean). The data starts at 09:00:00 on 05/17/2007.

Label	Start Time	End Time	Duration	1: Sys	1: LVEDP	1: HR	1: +dP/dt	2: Sys	2: Dia	2: Mean
T0	0001:01:12 09:00:00 05/17/2007									
0s Postdose	0001:01:12 09:00:00 05/17/2007	0001:11:12 09:10:00 05/17/2007	0000:10:00							
Avg				143.64	12.826	106.97	4080.8	142.74	99.035	120.88
10m Post...	0001:11:12 09:10:00 05/17/2007	0001:21:12 09:20:00 05/17/2007	0000:10:00	119.18	6.5803	77.269	3435.4	119.90	82.338	101.84
Avg				116.34	3.4051	74.239	3672.5	115.41	75.577	95.801
20m Post...	0001:21:12 09:20:00 05/17/2007	0001:31:12 09:30:00 05/17/2007	0000:10:00							
Avg				119.50	3.1541	79.70	3804.6	117.94	75.991	96.84
30m Post...	0001:31:12 09:30:00 05/17/2007	0001:41:12 09:40:00 05/17/2007	0000:10:00							
Avg				118.41	6.4415	66.17	3441.6	118.18	79.706	98.018
40m Post...	0001:41:12 09:40:00 05/17/2007	0001:51:12 09:50:00 05/17/2007	0000:10:00							
Avg				121.06	5.8841	73.593	3605.2	119.84	79.766	99.597
50m Post...	0001:51:12 09:50:00 05/17/2007	0002:01:12 10:00:00 05/17/2007	0000:10:00							
Avg				137.34	5.8307	96.607	4829.7	136.31	89.754	112.78
1h Postdose	0002:01:12 10:00:00 05/17/2007	0002:11:12 10:10:00 05/17/2007	0000:10:00							
Avg				137.34	5.8307	96.607	4829.7	136.31	89.754	112.78
1h 10m Po...	0002:11:12 10:10:00 05/17/2007	0002:21:12 10:20:00 05/17/2007	0000:10:00							
Avg				137.34	5.8307	96.607	4829.7	136.31	89.754	112.78

Figure 9. – Data Reduction with Use fixed Control and T0 enabled.

Parser Segments can also be configured and used for Data Reduction. When using Parser Segments, the ability to add user defined labels is permitted. Figure 10 shows the Add button expanded under the Intervals/Labels field. All labels manually entered here will be fixed in the Data Reduction list view and any additional reduction line labels will be handled by Ponemah.

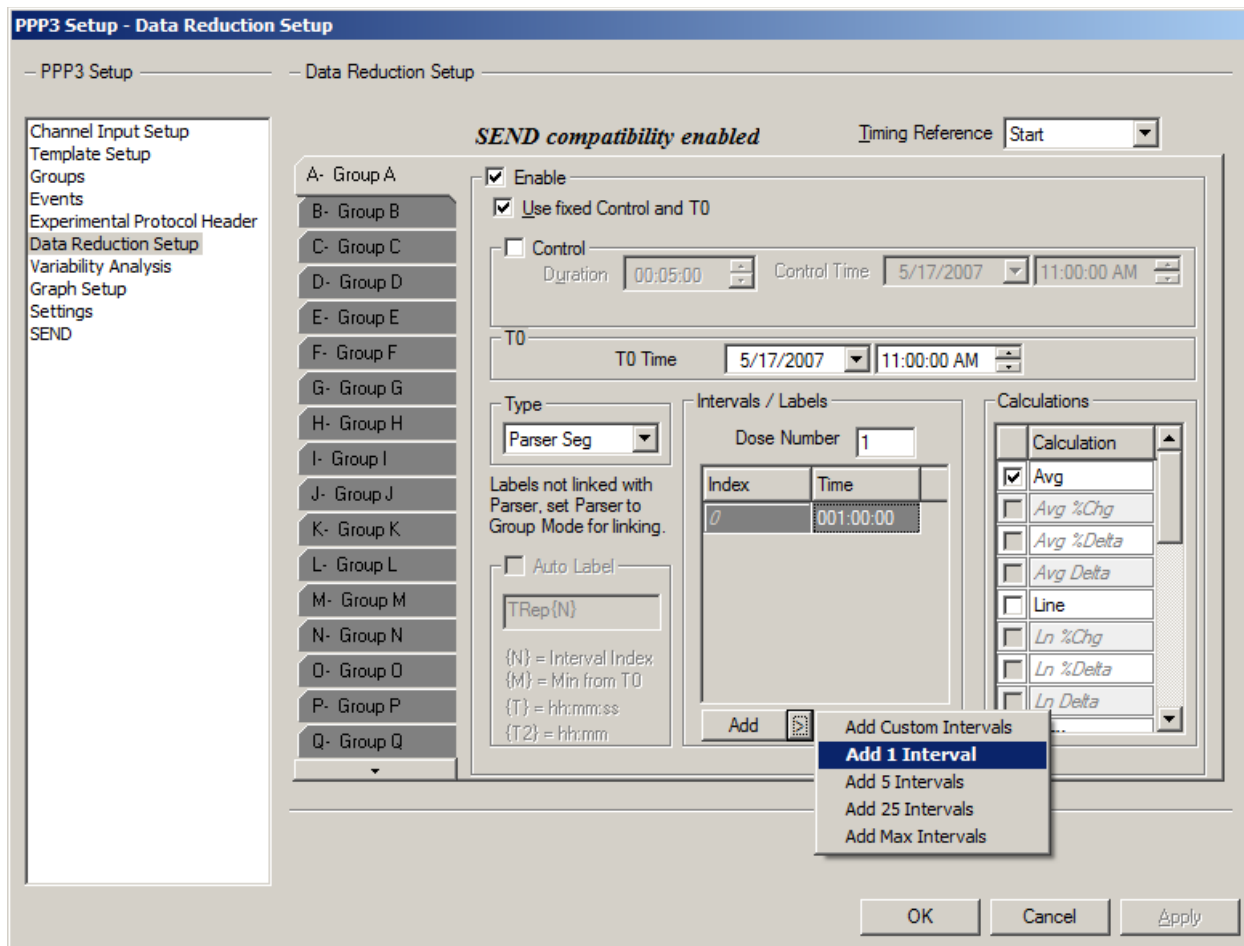


Figure 10. – Use of Parser Segments with Fixed Control enabled.

If Use fixed Control and T0 has been enabled, any parser segment prior to your T0 Time will be listed as a “Predose” time point. If Control has been enabled, the parser segment prior to your T0 Time will be used as your control value. If parser segments exist before the Control segment, they will be labeled as Predose.

Label	Start Time	End Time	Duration	1: Sys	1: Dia	1: Mean	1: PH
2h 38m 14s Predose	0000:00:0...	0000:01:0...	0000:01:00				
Avg				126.23	73.776	93.613	52.458
1h 38m 14s Predose	0001:00:0...	0001:01:0...	0000:01:00				
Avg				149.14	95.849	116.43	53.288
38m 14s Predose	0002:00:0...	0002:01:0...	0000:01:00				
Avg				163.28	104.51	126.84	58.773
T0	0002:39:1...						
21m 46s Postdose	0003:00:0...	0003:01:0...	0000:01:00				
Avg				134.39	81.159	101.41	53.232
1h 21m 46s Postdose	0004:00:0...	0004:01:0...	0000:01:00				
Avg				136.86	84.099	103.77	52.76
2h 21m 46s Postdose	0005:00:0...	0005:01:0...	0000:01:00				
Avg				138.31	81.749	102.57	56.566
3h 21m 46s Postdose	0006:00:0...	0006:01:0...	0000:01:00				

Figure 11. – Parser Segments prior to T0 Time shown as Predose, Control not enabled.

USING STUDY PROTOCOL OPTION AND SEND

The following outlines the differences in SEND options when using the Study Protocol Option. Refer to previous sections outlining the setup required for the SEND, Groups and Data Reduction dialogs to ensure requirements for generating SEND output have been configured.

SEND CONFIGURATION WHEN USING STUDY

If using the Study Protocol Option, the study configuration has been previously configured and saved as part of the Study prior to any collections. This includes Data Reduction, Subject IDs and other information. When using the Study Protocol Option, information can be automatically pulled into the SEND dialog and other menus. However, it may be necessary to modify information prior to generating SEND output to correct entry errors or further define a set of SEND generated data.

Functions specific to the Study Protocol Option in the SEND dialog include the ability to automatically load the Study Name, Study Day, and Workup. In order for the Study information to be displayed in the SEND dialog, you must have accessed a Review file within the Run Study dialog. With the Review file open, navigating to the SEND dialog will display the information for the Study that is currently accessed. Additionally, a Review file created within Study can be opened outside of Study with access to this information.

When the checkboxes have been disabled as in Figure 12, the information created using the Study option will be pulled into the SEND dialog. However, no modification of these fields will be allowed. Figure 12 shows CVStudy and Study Day 1 pulled from the Study configuration. Workup has no information since Workups were not configured as part of this Study.

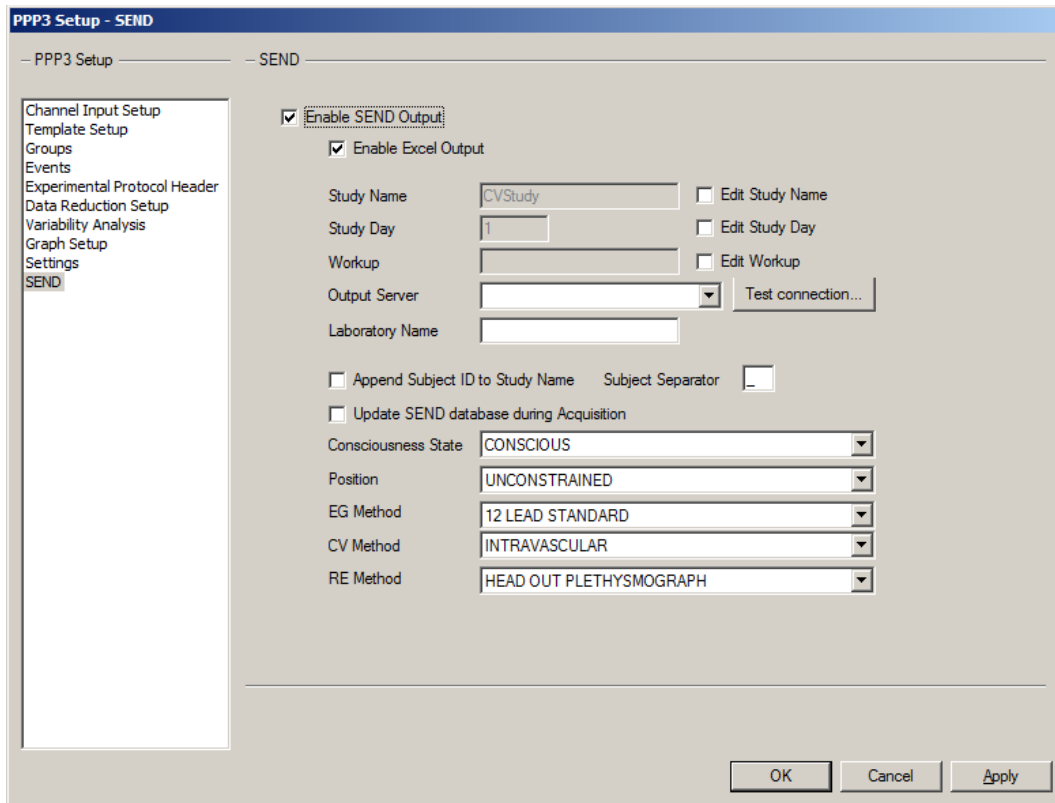


Figure 12. – Study Name, Study Day, and Workup pulled from the Study Protocol Option.

To modify Study information, place a check in the Edit Study Name, Edit Study Day, or Edit Workup checkbox. If the check marks beside these options are enabled as shown in Figure 13, these fields may be manually modified from their original values from when the Study was created.

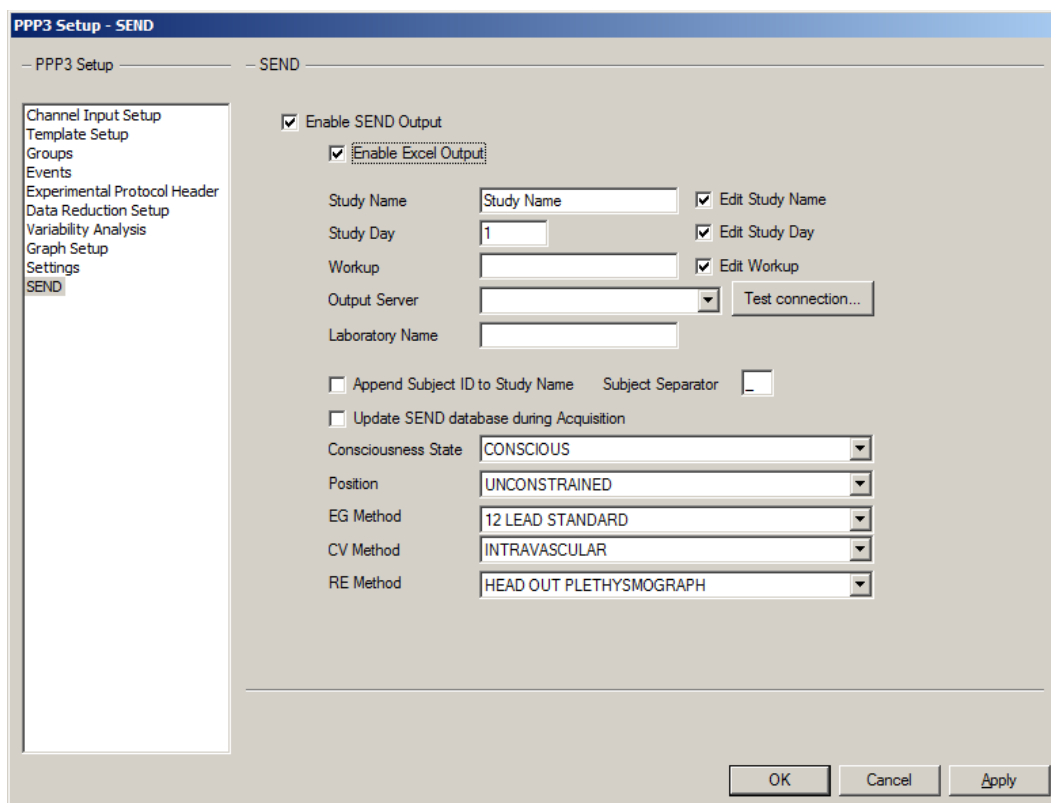


Figure 13. – Edit Study Name, Edit Study Day, and Edit Workup allows Study information to be edited.

As with the SEND dialog, the Groups dialog will also automatically fill Study Subject Name and Study Dose information. Deselecting Use Study Subjects and Use Study Doses will allow modification of the Study Protocol Option generated information. To modify, simply type in the desired information for each Subject (Group).

The information populated here is based on the information contained in the Review file opened when in the Run Study dialog or outside of Study. For example, Study Dose information will update appropriately if running a Latin Square design based on the information contained in the Review file that was opened.

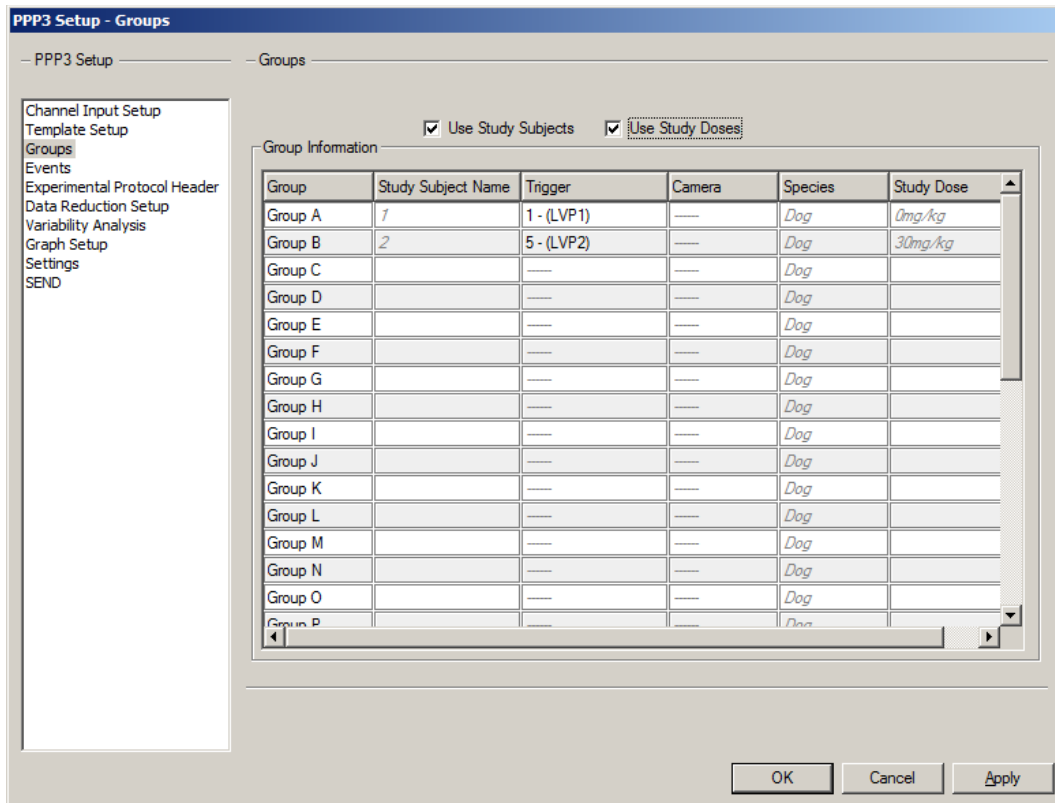


Figure 14. – Groups dialog with subject and dose information pulled from Study.

SEND MANAGER

SEND Manager is available from the SEND pull-down menu in Ponemah while in idle mode and allows you to access collections that have been previously saved. If in Review, SEND Manager will not be available (selection will be grayed out).

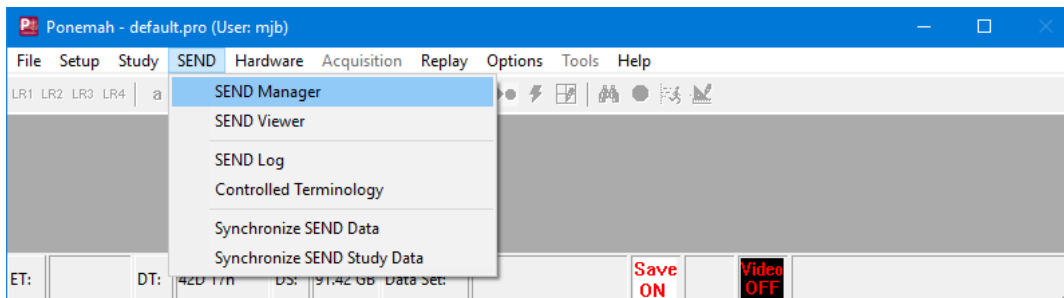


Figure 15. – SEND Manager access point from the SEND pull-down menu.

In addition to listing the collections available with SEND information, SEND Manager has backup and restore capabilities as described below. Each time a unique SEND study name is created, and data is saved, that study name will be listed in the dialog below.

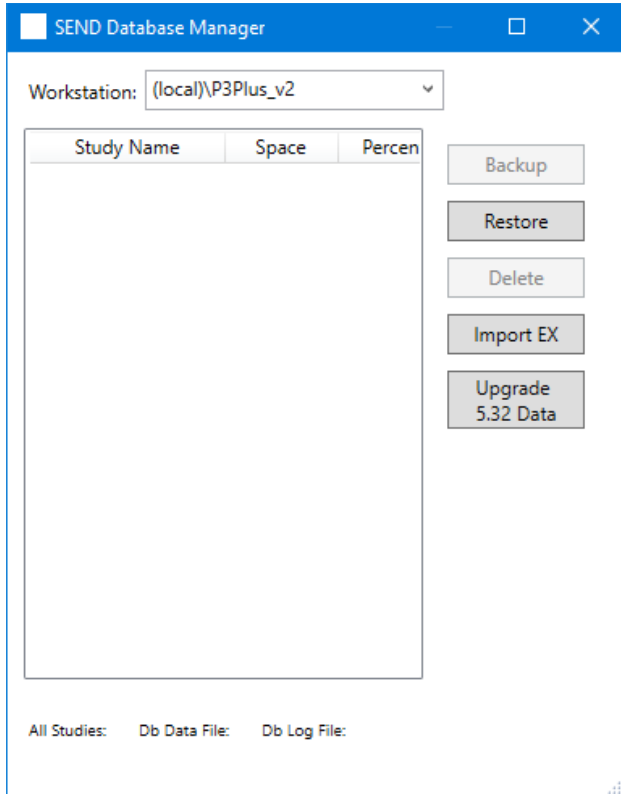


Figure 16. – SEND Manager dialog.

1 – Workstation

Lists the local Ponemah SQL Instance, P3Plus_v2, by default. If networked, users can select the P3Plus_v2 Instances on other networked workstations.

2 – Study Name and Space

Lists all available study names with previously saved SEND information and the amount of disk space that each study occupies.

3 – Backup

Allows the studies to be backed up and moved to other workstations or for archival purposes. The backup dialog is shown below. Use the typical Windows navigation to browse to the desired folder location and specify a file name for the backup file. The file type will automatically default to “.sendbkup”.

NOTE: If migrating the SEND database from the version found in 5.32 to 5.40, you may first want to backup all SEND databases. See Upgrade 5.32 Data below.

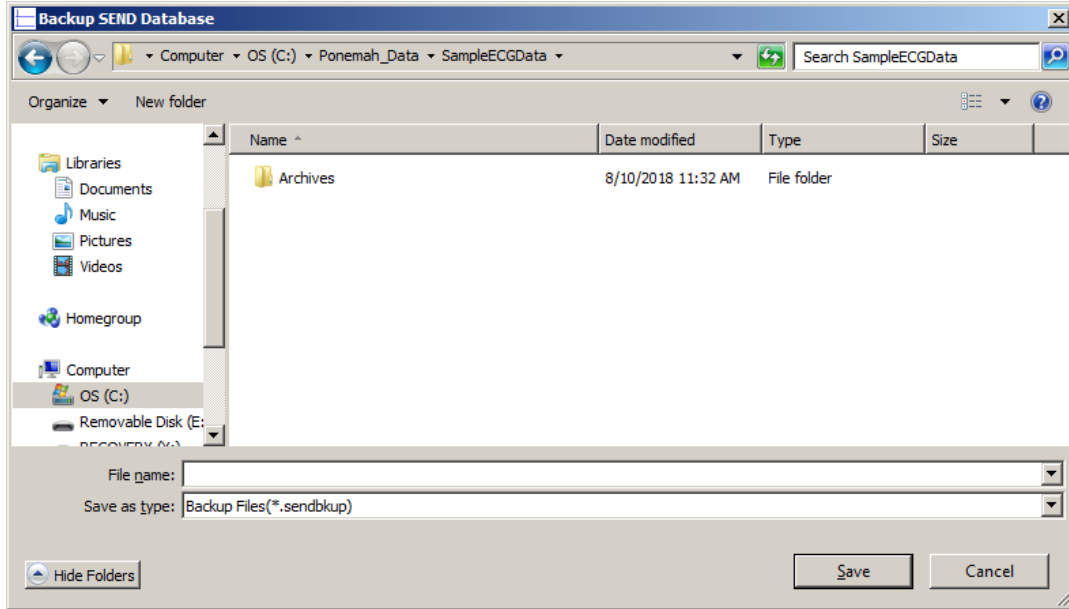


Figure 17. – Example SEND backup file being saved.

4 – Restore

Allows studies to be restored from other workstations or reopened from an archival location. This feature functions similar to that described under Backup shown above.

5 – Delete

The Delete function will remove the Study from the list in the SEND Manager dialog and will also remove any SEND associated data (SEND database) with the study from the SQL database. Size information (All Studies, Db Data File, and Db Log File) is provided at the bottom of the window. This change will be reflected under All Studies as the data size will decrease. Even though the SEND data was deleted, the space allocated for that data will remain and may appear as if the size of the database has not changed for the Db Data and Log Files. When new SEND data is generated, that space will be utilized by the new data.

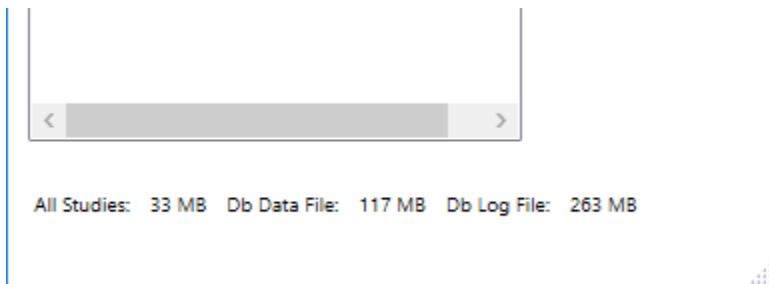


Figure 18. – SEND Manager Data size information.

6 – Import EX

The ability to import information is possible by creating a CSV file and using the Import EX function in SEND Database Manager. The required columns for successful import are STUDYID, USUBJID, EXSTDTC, EXSTDY, and optionally EXENDTC may be provided. Exposure domain information is not available in Ponemah, although some Variables require exposure data. Exposure data can be brought into Ponemah by importing an EX domain that has been exported from a LIMS system or other repository.

The following variables will report NULL unless an Exposure domain containing actual dose times is imported (DY, ENDY, RFTDTC). The EX domain can be imported at any point prior to exporting the CSV.

7 – Upgrade 5.32 Data

If a 5.32 SEND database exists, the user will be prompted to upgrade the SEND database in order to be compatible with Ponemah version 5.40. This is also true if attempting to access a remote database. Launching Ponemah 5.40 with a previous version of the SEND database being utilized will display the following message and allow the conversion of the 5.32 SEND database to be completed. Selecting Cancel will not allow the upgrade and automatically close Ponemah. The Upgrade 5.32 Data button, in SEND Manager, should be used to update the SEND database.

NOTE: It is highly recommended to back up the SEND database prior to upgrading to version 5.40

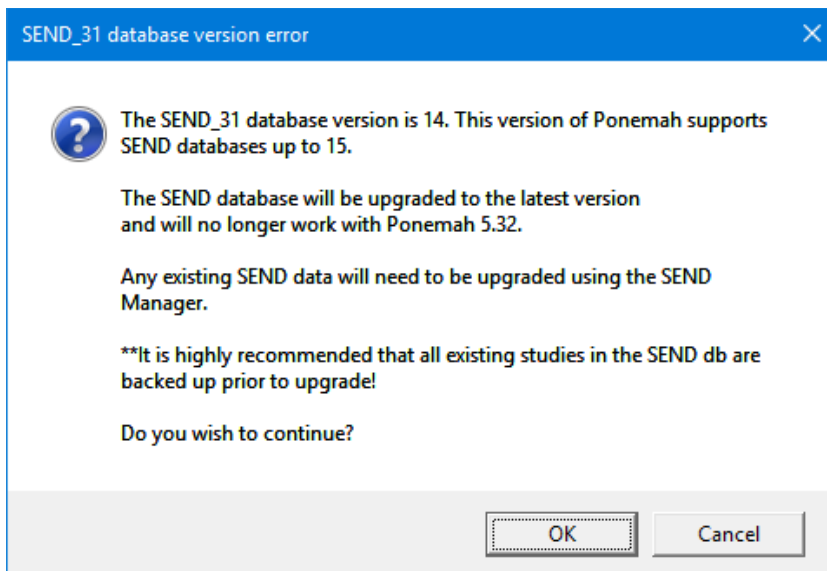



Figure 19. – Message confirming that the SEND database will be upgraded for version 5.40.

A dialog will post listing the data being converted along with a progress dialog.



Once the SEND database has been converted, the Upgrade 5.32 Data button will no longer be displayed in the SEND Manager dialog.

SEND VIEWER AND OUTPUT

The following sections detail the output options available for SEND generated data and the tools which can be used to view and modify the output.

SEND EXPORT TO EXCEL

To generate SEND data to Excel, the Enable Excel Output function must be selected in the SEND setup dialog. Excel output is available for both license file options, SEND Output and SEND SQL Server Output. Additionally, users must have enabled Data Reduction and selected derived parameters in the Channel Input Setup menu that correspond to output designated by CDISC (refer to SEND Setup from PPP3 Setup section for configuration information).

SEND output will be generated and included in the Microsoft Excel workbook created upon Saving a Marks Section (Run) or Derived Data within a Review session along with other Ponemah data related worksheets. Note that not all Ponemah derived parameters are defined by CDISC. Only those CDISC defined variables will be present in the output generated.

	C	D	E	F	G	H	I	J	K	L	M	
1	USUBJID	CVSEQ	CVTESTCD	CVTEST	CVPOS	CVORRES	CVORRESU	CVSTRESSC	CVSTRESN	CVSTRESU	CVSTAT	CVRE
2	CV	Group A	LVSYSBP	Left Ventricular Systolic Pressure	UNCONSTRAINED	118.40	'mmHg'	118.40	118.40	'mmHg'	NULL	NULL
3	CV	Group A	LVEDP	Left Ventricular End Diastolic Pressure	UNCONSTRAINED	5.5708	'mmHg'	5.5708	5.5708	'mmHg'	NULL	NULL
4	CV	Group A	HR	Heart Rate	UNCONSTRAINED	74.319	'bpm'	74.319	74.319	'bpm'	NULL	NULL
5	CV	Group A	DPDTAVG	dP/dt Average	UNCONSTRAINED	3310.2	'mmHg/s'	3310.2	3310.2	'mmHg/s'	NULL	NULL
6	CV	Group A	SYSBP	Systolic Blood Pressure	UNCONSTRAINED	117.87	'mmHg'	117.87	117.87	'mmHg'	NULL	NULL
7	CV	Group A	DIABP	Diastolic Blood Pressure	UNCONSTRAINED	81.248	'mmHg'	81.248	81.248	'mmHg'	NULL	NULL
8	CV	Group A	MAP	Mean Arterial Pressure	UNCONSTRAINED	99.849	'mmHg'	99.849	99.849	'mmHg'	NULL	NULL
9	CV	Group A	PULSEPR	Pulse Pressure	UNCONSTRAINED	36.621	'mmHg'	36.621	36.621	'mmHg'	NULL	NULL
10	CV	Group A	ACTIVITY	Activity	UNCONSTRAINED	0.0164	"	0.0164	0.0164	"	NULL	NULL
11	CV	Group A	LVSYSBP	Left Ventricular Systolic Pressure	UNCONSTRAINED	111.20	'mmHg'	111.20	111.20	'mmHg'	NULL	NULL
12	CV	Group A	LVEDP	Left Ventricular End Diastolic Pressure	UNCONSTRAINED	4.7871	'mmHg'	4.7871	4.7871	'mmHg'	NULL	NULL
13	CV	Group A	HR	Heart Rate	UNCONSTRAINED	60.679	'bpm'	60.679	60.679	'bpm'	NULL	NULL
14	CV	Group A	DPDTAVG	dP/dt Average	UNCONSTRAINED	3023.4	'mmHg/s'	3023.4	3023.4	'mmHg/s'	NULL	NULL
15	CV	Group A	SYSBP	Systolic Blood Pressure	UNCONSTRAINED	112.41	'mmHg'	112.41	112.41	'mmHg'	NULL	NULL
16	CV	Group A	DIABP	Diastolic Blood Pressure	UNCONSTRAINED	76.446	'mmHg'	76.446	76.446	'mmHg'	NULL	NULL
17	CV	Group A	MAP	Mean Arterial Pressure	UNCONSTRAINED	94.795	'mmHg'	94.795	94.795	'mmHg'	NULL	NULL
18	CV	Group A	PULSEPR	Pulse Pressure	UNCONSTRAINED	35.967	'mmHg'	35.967	35.967	'mmHg'	NULL	NULL
19	CV	Group A	ACTIVITY	Activity	UNCONSTRAINED	NULL	NULL	NULL	NULL	NULL	'NOT DONE'	'INCC'
20	CV	Group A	LVSYSBP	Left Ventricular Systolic Pressure	UNCONSTRAINED	115.28	'mmHg'	115.28	115.28	'mmHg'	NULL	NULL
21	CV	Group A	LVEDP	Left Ventricular End Diastolic Pressure	UNCONSTRAINED	5.6705	'mmHg'	5.6705	5.6705	'mmHg'	NULL	NULL
22	CV	Group A	HR	Heart Rate	UNCONSTRAINED	64.359	'bpm'	64.359	64.359	'bpm'	NULL	NULL
23	CV	Group A	DPDTAVG	dP/dt Average	UNCONSTRAINED	3043.0	'mmHg/s'	3043.0	3043.0	'mmHg/s'	NULL	NULL
24	CV	Group A	SYSBP	Systolic Blood Pressure	UNCONSTRAINED	117.20	'mmHg'	117.20	117.20	'mmHg'	NULL	NULL
25	CV	Group A	DIABP	Diastolic Blood Pressure	UNCONSTRAINED	78.698	'mmHg'	78.698	78.698	'mmHg'	NULL	NULL
26	CV	Group A	MAP	Mean Arterial Pressure	UNCONSTRAINED	98.12	'mmHg'	98.12	98.12	'mmHg'	NULL	NULL
27	CV	Group A	PULSEPR	Pulse Pressure	UNCONSTRAINED	38.50	'mmHg'	38.50	38.50	'mmHg'	NULL	NULL

Figure 20. – Example output when enabling Excel in the SEND dialog.

If the collection contains data from more than one Domain, multiple worksheets will be created within the Excel workbook. Since the study shown in Figure 20 incorporated both CV and EG domain data, two worksheets are created in the Excel workbook with the appropriate information for each Domain.

SEND VIEWER

Selecting SEND Viewer from the SEND pulldown menu will display the following dialogs. In order to view SEND information, users must have enabled Data Reduction and selected derived parameters in the Channel Input Setup menu that correspond to output designated by CDISC (refer to SEND Setup from PPP3 Setup section for configuration information).

SEND Viewer allows selection and filtering of the study data and information for export to a CSV formatted file to be imported into other systems. Figure 21 shows the SEND Viewer dialog populated with data from the CV Domain and the associated CDISC controlled terminology displayed in the column headers. Scroll bars on the bottom and right side of the dialog allow the user to scroll and view the entire contents of the window.

The Workstation pulldown menu lists the local Ponemah SQL instance, P3Plus_v2, by default along with any designated Output Server that has been configured. If networked, users can select P3Plus_v2 Instances on other

networked workstations or servers. This allows access to the local Ponemah SEND database as well as other Ponemah SEND databases on workstations that are accessible on the network. Selecting a workstation will display any previously collected information (studies) that is available for export. The pull-down menu will be populated with workstation information if studies have been created using the Study Protocol Option. All workstations that were configured for a given study will be loaded and available via the pull-down menu. The list of workstations available may change based on the study that is selected in the Manage Studies dialog.

Selecting the Options in the upper left-hand side of the dialog will expand and display filters that allow you to select the desired information for export. This includes the study, subjects, days and CDISC defined parameters.

STUDYID	DOMAIN	USUBJID	CVSEQ	CVTESTCD	CVTEST	CVPOS	CVORRES	CVORRESU
CV Test1	CV	Group A	2321	LVSYSBP	Left Ventricular Systolic Pressure	UNCONSTRAINED	118.37	mmHg
CV Test1	CV	Group A	2322	LVEDP	Left Ventricular End Diastolic Pressure	UNCONSTRAINED	5.588	mmHg
CV Test1	CV	Group A	2323	HR	Heart Rate	UNCONSTRAINED	74.33	bpm
CV Test1	CV	Group A	2324	DPDTAVG	dP/dt Average	UNCONSTRAINED	3310.4	mmHg/s
CV Test1	CV	Group A	2325	SYSBP	Systolic Blood Pressure	UNCONSTRAINED	117.83	mmHg
CV Test1	CV	Group A	2326	DIABP	Diastolic Blood Pressure	UNCONSTRAINED	81.212	mmHg
CV Test1	CV	Group A	2327	MAP	Mean Arterial Pressure	UNCONSTRAINED	99.818	mmHg
CV Test1	CV	Group A	2328	PULSEPR	Pulse Pressure	UNCONSTRAINED	36.623	mmHg
CV Test1	CV	Group A	2329	DPDTMAX	Left Ventricular Maximum Positive dP/dt	UNCONSTRAINED	10566	mmHg/s
CV Test1	CV	Group A	2330	DPDTMIN	Left Ventricular Minimum Positive dP/dt	UNCONSTRAINED	2565.3	mmHg/s
CV Test1	CV	Group A	2331	LVSYSBP	Left Ventricular Systolic Pressure	UNCONSTRAINED	111.21	mmHg
CV Test1	CV	Group A	2332	LVEDP	Left Ventricular End Diastolic Pressure	UNCONSTRAINED	4.7925	mmHg
CV Test1	CV	Group A	2333	HR	Heart Rate	UNCONSTRAINED	60.651	bpm
CV Test1	CV	Group A	2334	DPDTAVG	dP/dt Average	UNCONSTRAINED	3023.2	mmHg/s
CV Test1	CV	Group A	2335	SYSBP	Systolic Blood Pressure	UNCONSTRAINED	112.43	mmHg
CV Test1	CV	Group A	2336	DIABP	Diastolic Blood Pressure	UNCONSTRAINED	76.463	mmHg
CV Test1	CV	Group A	2337	MAP	Mean Arterial Pressure	UNCONSTRAINED	94.816	mmHg
CV Test1	CV	Group A	2338	PULSEPR	Pulse Pressure	UNCONSTRAINED	35.969	mmHg
CV Test1	CV	Group A	2339	DPDTMAX	Left Ventricular Maximum Positive dP/dt	UNCONSTRAINED	3589.6	mmHg/s
CV Test1	CV	Group A	2340	DPDTMIN	Left Ventricular Minimum Positive dP/dt	UNCONSTRAINED	2594.9	mmHg/s
CV Test1	CV	Group A	2341	LVSYSBP	Left Ventricular Systolic Pressure	UNCONSTRAINED	115.27	mmHg
CV Test1	CV	Group A	2342	LVEDP	Left Ventricular End Diastolic Pressure	UNCONSTRAINED	5.6718	mmHg
CV Test1	CV	Group A	2343	HR	Heart Rate	UNCONSTRAINED	64.312	bpm
CV Test1	CV	Group A	2344	DPDTAVG	dP/dt Average	UNCONSTRAINED	3042.7	mmHg/s
CV Test1	CV	Group A	2345	SYSBP	Systolic Blood Pressure	UNCONSTRAINED	117.19	mmHg
CV Test1	CV	Group A	2346	DIABP	Diastolic Blood Pressure	UNCONSTRAINED	78.691	mmHg
CV Test1	CV	Group A	2347	MAP	Mean Arterial Pressure	UNCONSTRAINED	98.115	mmHg
CV Test1	CV	Group A	2348	PULSEPR	Pulse Pressure	UNCONSTRAINED	38.502	mmHg
CV Test1	CV	Group A	2349	DPDTMAX	Left Ventricular Maximum Positive dP/dt	UNCONSTRAINED	4766.7	mmHg/s

6235 records total

Figure 21. – SEND Viewer dialog.

Figure 22 shows the Options pin expanded. Four filters are available to tailor the SEND output to CSV format.

1 – Study

Lists previously collected studies that are present on the workstation selected. Selecting a study will load any SEND related information. Figure 22 shows CVTest1 selected and its associated data displayed.

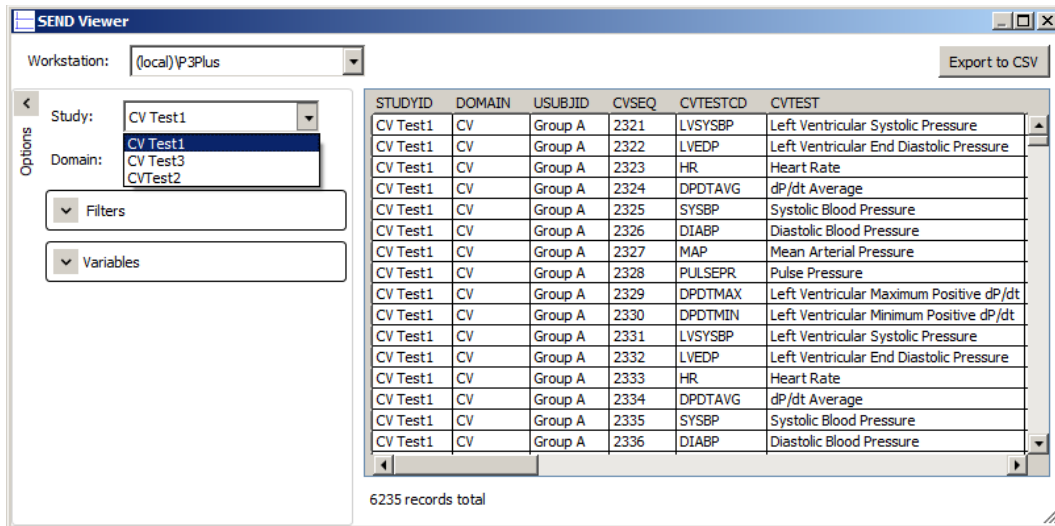


Figure 22. – SEND Viewer dialog.

2 – Domain

After the Study has been chosen, the correct SEND Domain must be selected. If an incorrect Domain is selected, the appropriate data will not populate in the window.

3 – Filters

Expanding the Filters option allows the user to define what information will be displayed and available for output to CSV. This includes USUBJID, NOMDAY, and TESTCD for the given Domain.

The “All” checkbox allows the user to toggle all variables on or off. Additionally, each variable may be selected or deselected individually to create the desired output.

It is important to note that not all Ponemah generated parameters are required by CDISC (TESTCD). Only parameters defined by CDISC will be displayed.

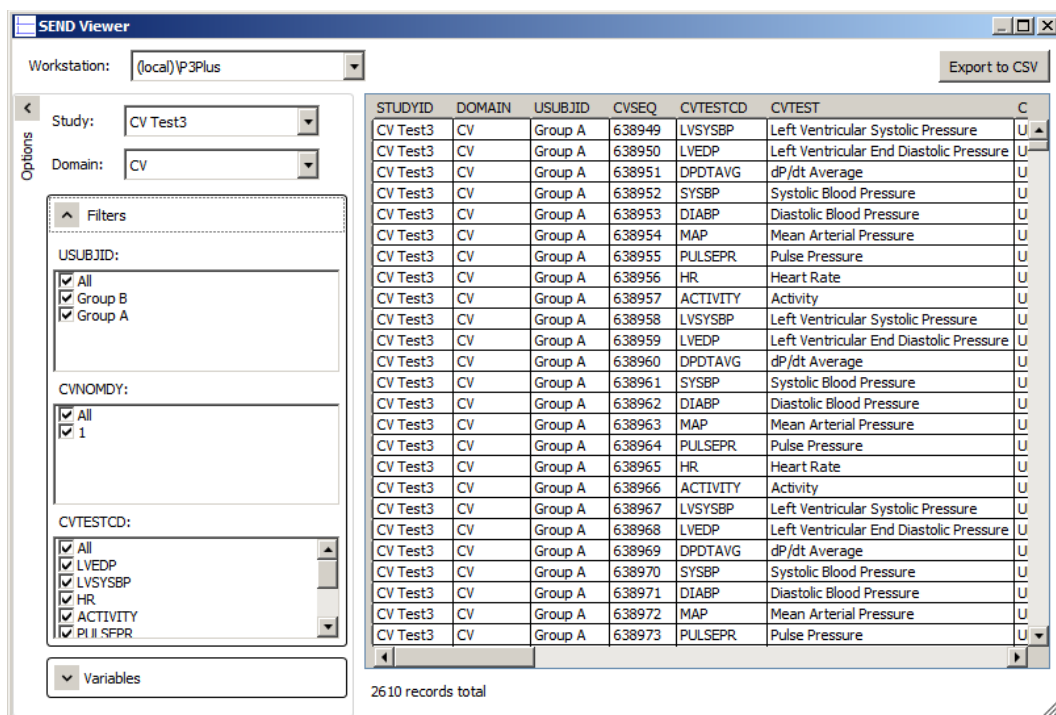


Figure 23. – SEND Viewer dialog displaying available filters.

4 – Variables

The Variables option controls which CDISC defined variables will be displayed and outputted to the CSV formatted file. Toggling specific variables will enable or disable the variables in the viewer and the associated CSV file when generated. This may be used to remove variables that are not populated with information.

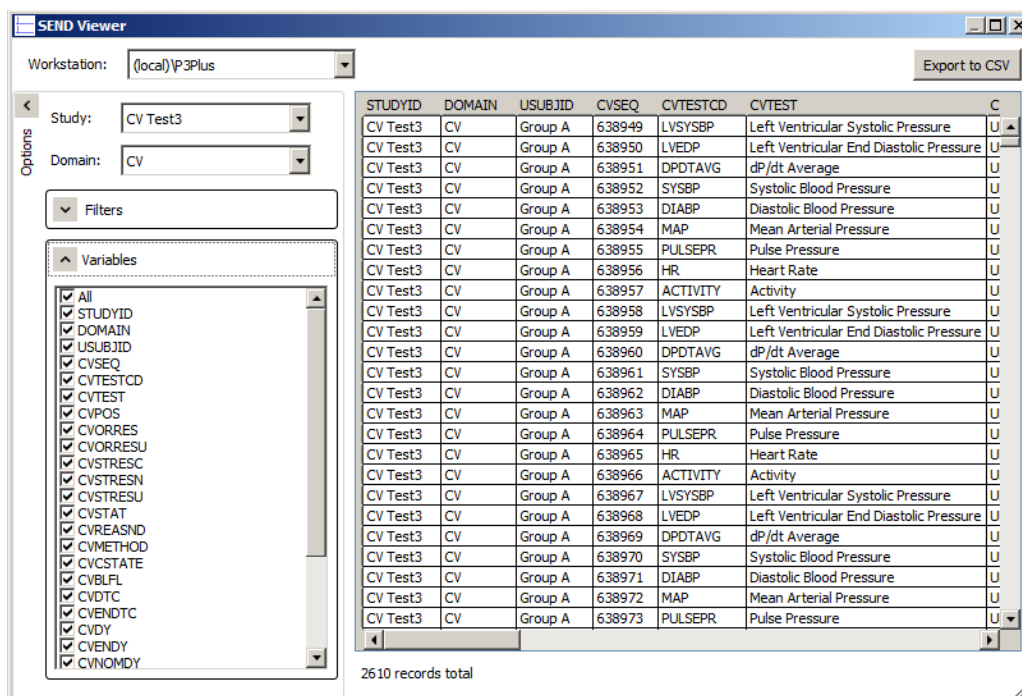


Figure 24. – SEND Viewer dialog with Variables option expanded.

In the example study above (CVTest3 from Figure 24), EG Domain data was collected along with CV Domain data. In order to get the EG Domain variables, simply select the EG Domain as shown in Figure 25 and update the remaining Options to provide the desired output.

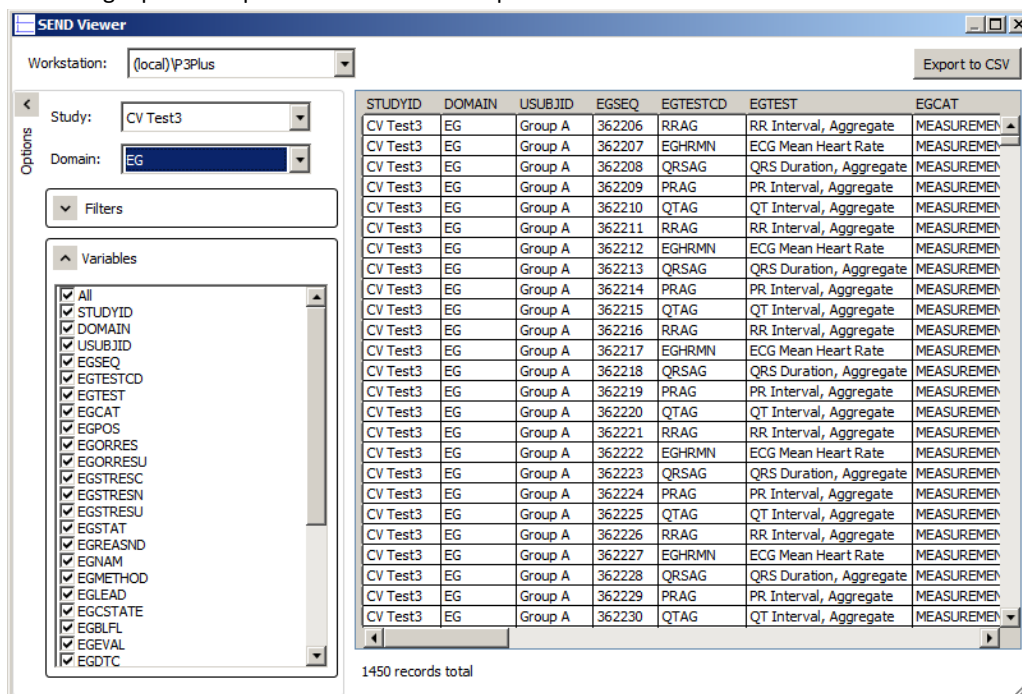


Figure 25. – SEND Viewer dialog with Variables option expanded and EG Domain selected.

SEND EXPORT TO CSV

Once the appropriate filters and variables have been defined as outlined in the SEND Viewer section, the data is ready to export to a CSV formatted file. This is accomplished by selecting the Export to CSV button in the top, right-hand portion of the dialog.

NOTE: It is important to note that this functionality is only available if the SEND SQL Server Option has been enabled in the license file. If only SEND Output is enabled, these features will not be available.

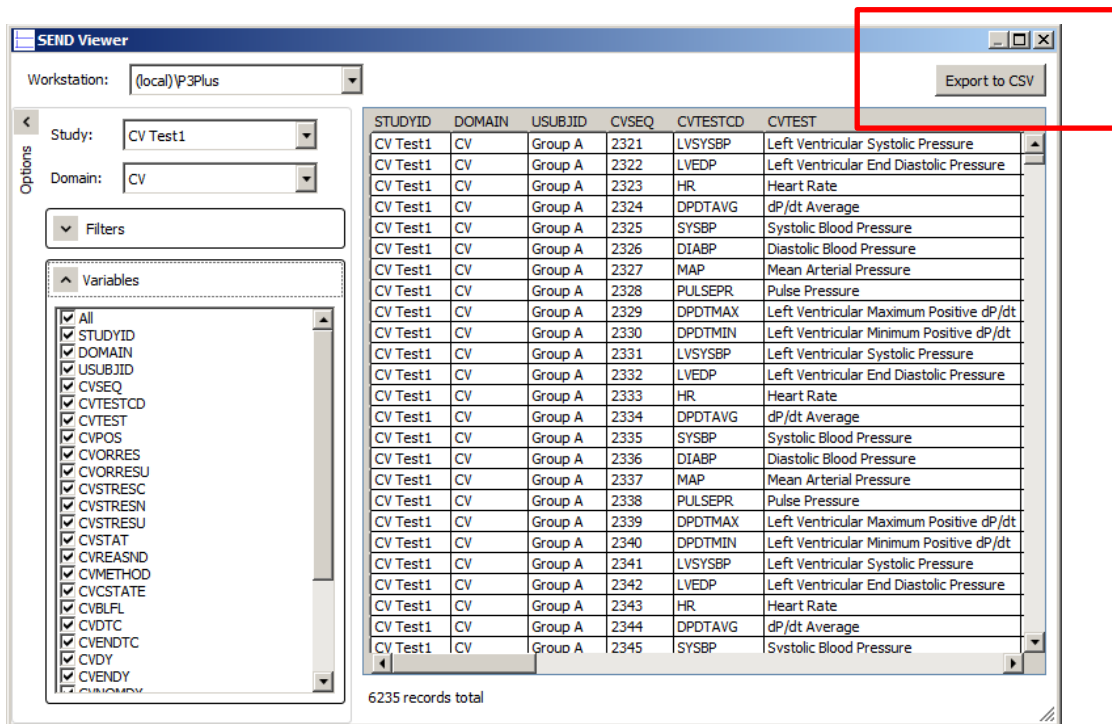


Figure 26. – SEND Viewer dialog with Export to CSV button highlighted.

A Windows dialog will be displayed with the default location set to the study folder within the Ponemah_Data directory. The file name will include the study name and date and will be configured to save in the CSV format. The name and location for saving the file may be changed by modifying the “File name” edit field and using the browse feature to specify a different location.

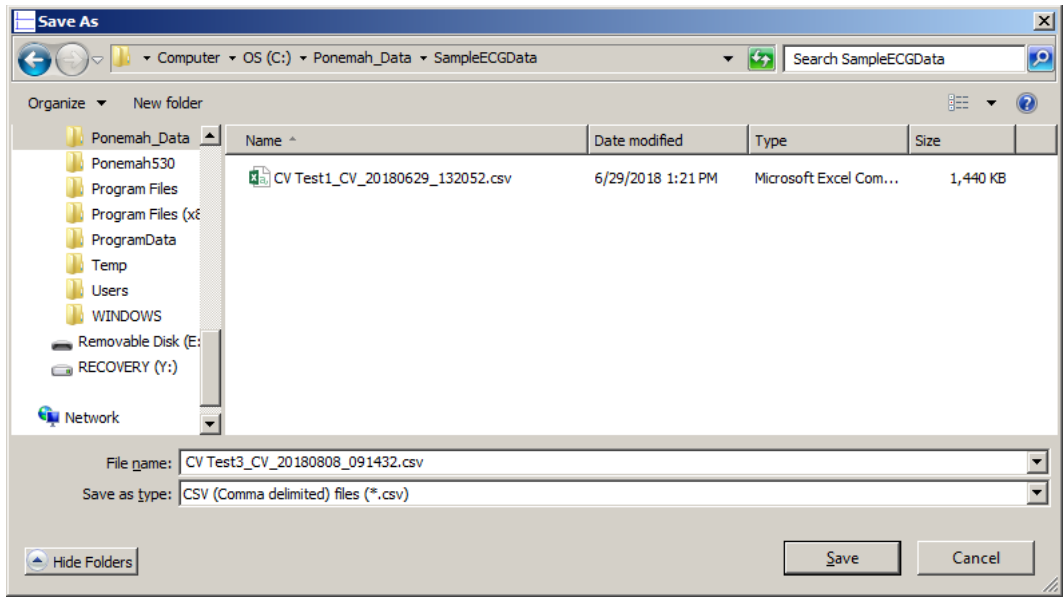


Figure 27. – Default location and file type for saving SEND output via SEND Viewer.

Once the file has been saved, browsing to the location and opening the file will display the data as configured in SEND Viewer. Figure 28 shows the CSV output opened using Microsoft Excel.

	A	B	C	D	E	F	G	H	I	J
1	STUDYID	DOMAIN	USUBJID	CVSEQ	CVTESTCD	CVTEST	CVPOS	CVORRES	CVORRESI	CVSTRESC
2	CV Test1	CV	Group A	2321	LVSYSBP	Left Ventricular Systolic Pressure	UNCONSTRAINED	118.37 mmHg		118.37
3	CV Test1	CV	Group A	2322	LVEDP	Left Ventricular End Diastolic Pressure	UNCONSTRAINED	5.5588 mmHg		5.5588
4	CV Test1	CV	Group A	2323	HR	Heart Rate	UNCONSTRAINED	74.33 bpm		74.33
5	CV Test1	CV	Group A	2324	DPDTAVG	dP/dt Average	UNCONSTRAINED	3310.4 mmHg/s		3310.4
6	CV Test1	CV	Group A	2325	SYSBP	Systolic Blood Pressure	UNCONSTRAINED	117.83 mmHg		117.83
7	CV Test1	CV	Group A	2326	DIABP	Diastolic Blood Pressure	UNCONSTRAINED	81.212 mmHg		81.212
8	CV Test1	CV	Group A	2327	MAP	Mean Arterial Pressure	UNCONSTRAINED	99.818 mmHg		99.818
9	CV Test1	CV	Group A	2328	PULSEPR	Pulse Pressure	UNCONSTRAINED	36.623 mmHg		36.623
10	CV Test1	CV	Group A	2329	DPDTMAX	Left Ventricular Maximum Positive dP/dt	UNCONSTRAINED	10566 mmHg/s		10566
11	CV Test1	CV	Group A	2330	DPDTMIN	Left Ventricular Minimum Positive dP/dt	UNCONSTRAINED	2565.3 mmHg/s		2565.3
12	CV Test1	CV	Group A	2331	LVSYSBP	Left Ventricular Systolic Pressure	UNCONSTRAINED	111.21 mmHg		111.21
13	CV Test1	CV	Group A	2332	LVEDP	Left Ventricular End Diastolic Pressure	UNCONSTRAINED	4.7925 mmHg		4.7925
14	CV Test1	CV	Group A	2333	HR	Heart Rate	UNCONSTRAINED	60.651 bpm		60.651
15	CV Test1	CV	Group A	2334	DPDTAVG	dP/dt Average	UNCONSTRAINED	3023.2 mmHg/s		3023.2
16	CV Test1	CV	Group A	2335	CVSBP	Systolic Blood Pressure	UNCONSTRAINED	112.43 mmHg		112.43

Figure 28. – Example of CSV output.

SEND LOG

The SEND Log provides information regarding changes to the data stored in the SEND database on the local system or the server.

NOTE: Migrating from version 5.32 to 5.40 will result in the SEND database from version 5.32 being converted to conform to the current schema used in Ponemah version 5.40. After conversion, the database will no longer be functional for version 5.32. Therefore, consideration must be taken to determine if separate storage (servers or workstations) and Ponemah 5.32 acquisition and Post systems should be maintained separately from 5.4 workstations and servers. Or, if updating the 5.32 SEND database is applicable to your given situation. Ponemah versions 5.32 and 5.4 cannot be used simultaneously for generating SEND output. If a migration has been performed, this will be listed in the SEND log as “Upgrade” under the Action column. See also Upgrade 5.32 Data. Only log information pertaining to the most recently saved data for a given collection is preserved in the SEND Log. This is due to the fact that the SEND database only contains the most recently saved information for a given subject or collection of subjects from a collection. Data is not appended in the SEND database. Only data that matches the Study and subject for the associated acquisition (day and time) is replaced. For example, if Day 7 data is analyzed and SEND data is generated, only Day 7 data would be affected, not previously saved SEND data for prior collection days.

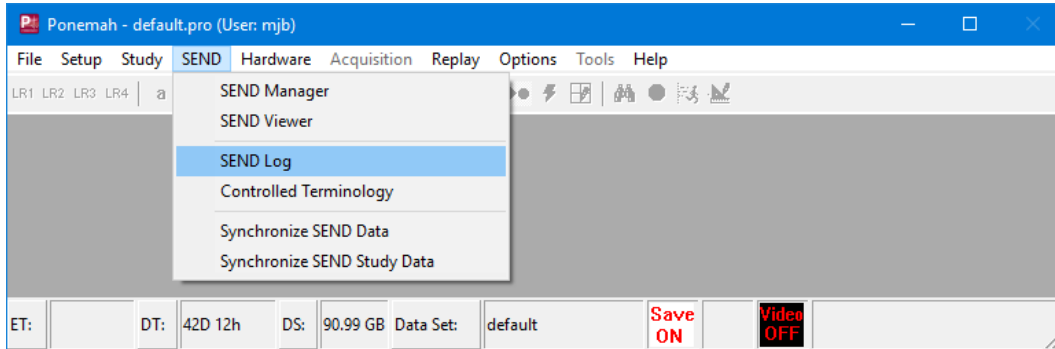


Figure 29. – SEND Log access point from the SEND pull-down menu.

The SEND Log displayed below will provide information about changes to the SEND database. Additional details may be found in the Application Log depending on the function performed.

The screenshot shows the SEND Log window with a table of records. The workstation is set to '(local)\P3Plus_v2'. The table has columns for Domain, NOMDY, USUBJID, Review File Name, RAW File Start, RAW File End, User, Wc, Action Date, and Action. There are 4 records total.

Domain	NOMDY	USUBJID	Review File Name	RAW File Start	RAW File End	User	Wc	Action Date	Action
CV	1	Group A	CVStudy1 (Day 1) (Group 1) (80830009).RVW	2019-09-19T16:15:02	2019-09-19T16:18:24	mjb	MB	2019-09-23T09:26:05.570	Save - Review
CV	1	Group B	CVStudy1 (Day 1) (Group 1) (80830009).RVW	2019-09-19T16:15:02	2019-09-19T16:18:24	mjb	MB	2019-09-23T09:26:05.840	Save - Review
CV	1	Group A	CVStudy1 (Day 1) (Group 1) (80830009).RVW	2019-09-19T16:15:02	2019-09-19T16:18:24	mjb	MB	2019-09-23T10:32:00.863	Delete
CV	1	Group A	CVStudy1 (Day 1) (Group 1) (80830009).RVW	2019-09-19T16:15:02	2019-09-19T16:18:24	mjb	MB	2019-09-23T11:33:17.213	Delete

Figure 30. – SEND Log.

1 – Workstation

Defaults to the local instance. However, if workstations are networked, users can access the SEND Log on other workstations.

2 – Options

The Options flyout is available in the left hand edge of the menu. Users can filter on Domain, Subjects, Review File Name, and all available columns in the log itself as shown below.

The screenshot shows the SEND Log application window. On the left, there is an 'Options' flyout menu with sections for 'Filters' and 'Columns'. The 'Filters' section includes checkboxes for 'Domain' (All, CV), 'USUBID' (All, Group A), 'NOMDY' (All, 1), 'Review File Name' (All, CVStudy1 (Day 1) (Group 1) (80830015).RVW), and 'User' (All, mjib). The 'Columns' section includes checkboxes for All, Action, Domain, NOMDY, RAW File End, RAW File Start, Review File Name, User, USUBID, and Workstation Name. The main area displays a table with 2 records. The table has columns: Domain, NOMDY, USUBID, Review File Name, RAW File Start, RAW File End, User, Wo, Action Date, and Action. The data rows are: CV | 1 | Group A | CVStudy1 (Day 1) (Group 1) (80830015).RVW | 2019-09-19T16:15:02 | 2019-09-19T16:18:24 | mjib | MB | 2019-09-30T15:03:13.790 | Save - Review; and CV | 1 | Group B | CVStudy1 (Day 1) (Group 1) (80830015).RVW | 2019-09-19T16:15:02 | 2019-09-19T16:18:24 | mjib | MB | 2019-09-30T15:03:14.800 | Save - Review. An 'Export to CSV' button is located in the top right corner. At the bottom left of the table area, it says '2 records total'.

Domain	NOMDY	USUBID	Review File Name	RAW File Start	RAW File End	User	Wo	Action Date	Action
CV	1	Group A	CVStudy1 (Day 1) (Group 1) (80830015).RVW	2019-09-19T16:15:02	2019-09-19T16:18:24	mjib	MB	2019-09-30T15:03:13.790	Save - Review
CV	1	Group B	CVStudy1 (Day 1) (Group 1) (80830015).RVW	2019-09-19T16:15:02	2019-09-19T16:18:24	mjib	MB	2019-09-30T15:03:14.800	Save - Review

Figure 31. – SEND Log showing available Options for filtering information presented in the Log.

3 – Export To CSV

Allows the log information to be exported to a .CSV file. The default location is Ponemah_Data but the file can be saved to any location using Windows browse function.

4 – Log Information

Displays information on the Domain, Day of collection, file information, User and the Action initiated which resulted in the log being created.

CONTROLLED TERMINOLOGY

The Controlled Terminology menu allows the user to view or modify the terminology used for SEND submission. Select Controlled Terminology from the SEND pull-down menu.

NOTE: If installing Ponemah version 5.40 over an existing 5.32 installation, Ponemah will automatically pull in the existing information contained in the SEND.xml file from version 5.32. Therefore, any modifications to the Controlled terminology, EG Methods, etc. from 5.32 will be preserved and loaded into 5.40. If 5.32 has been uninstalled, the SEND.xml simply needs to be present in the directory that 5.40 will be installed.

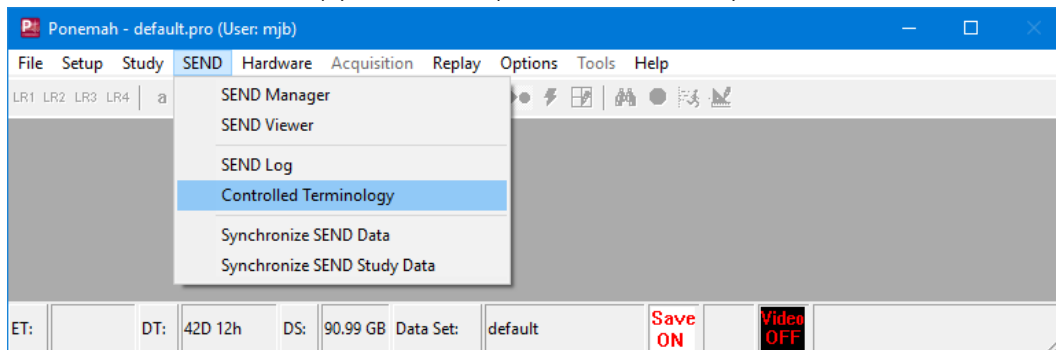


Figure 32. – Access point for Controlled Terminology.

The currently approved CDISC terminology has been prepopulated in the Test Codes menu shown below. Please note that this menu does not determine SEND output when data is saved. This is a table of the parameters and terminology that will be available in the SEND Viewer menu.

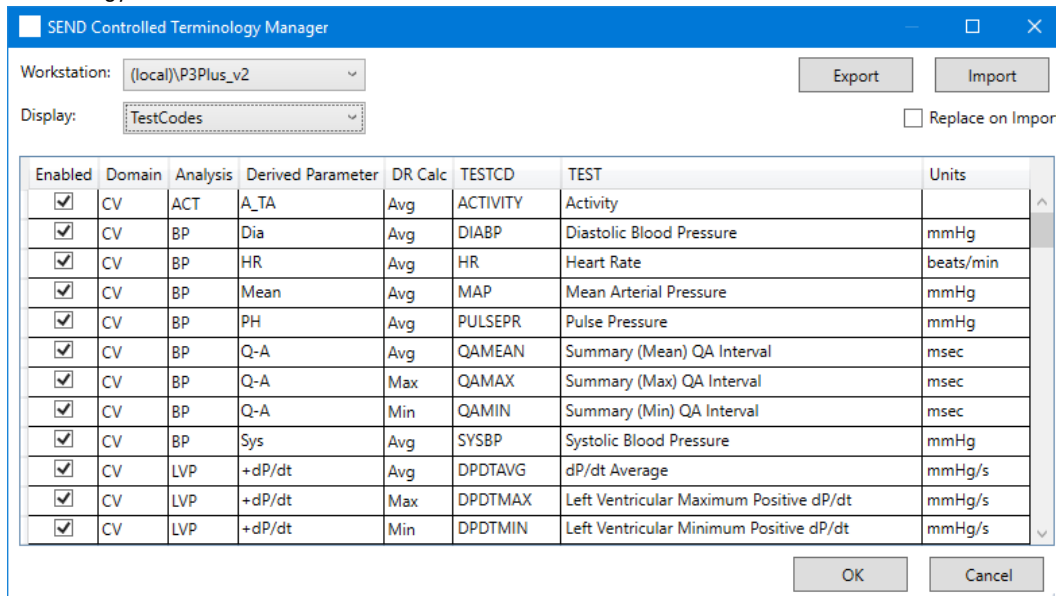


Figure 33. – Test Codes listing within Controlled Terminology.

1 – Workstation

Defaults to the local instance. However, if workstations are networked, users can access Controlled Terminology menus on other workstations.

2 – Display

Allows the selection of the following menu which can be modified by the user. The information entered here will be the available options displayed in SEND Viewer for generating SEND output.

- TestCodes
- EG Methods
- EG Leads
- CV Methods
- RE Methods
- CStates
- Position
- ECG Results

3 – Export

Allows the currently saved settings to be exported to a .xml file for importing into other workstations.

4 – Import

Allows previously exported .xml files to be imported. This will automatically update the current settings to minimize manual changes on each workstation. If information differs between the .xml file currently stored and the file being imported, the differences will be appended. Information is not overwritten using this function.

Each of the options available under the Display pull-down menu can be modified as mentioned. Figure 34 shows the addition of the TTI derived parameter. To add a parameter that is available in the Ponemah derived parameter list (but not defined by CDISC), scroll to the bottom of the list and click in the available cells under the various column headers. A drop down list will be displayed to provide available selections. In this example, CV was selected for the Domain, LVP was selected for the analysis module, and TTI was the derived parameter of interest. The DR Calc column shows the cell being modified and the three available options.

When adding parameters not defined by default, TESTCD, TEST, and Units will be manually entered.

The option to disable and delete entries is also available. In the case of the TestCode menu, deselecting the check mark under the Enable column for a Test Code will preserve the entry but will remove it from the SEND Viewer menu and not allow reporting of that parameter through SEND. When a TestCode has been disabled, it will drop down to the bottom of the list for that Domain (list is sorted by Domain).

To delete an entry, right click on the desired item and the option to delete will be available. Multiple entries can be selected and deleted at once using the shift key. Note that entries will be removed from the list and will need to be manually entered if needed in the future.

It is recommended to Export the list prior to making any changes and save the .xml file for future use. When importing a file, the Controlled Terminology dialog must be closed and reopened to view changes. Any new entries will be displayed in the list but will be disabled by default.

The Replace on Import checkbox provides a method to remove all information currently available in the Controlled Terminology dialog. Check this box prior to importing an xml file. Next, use the Import button to browse to and load the desired SEND xml file. Only the contents of the xml file being loaded will be available within the Controlled Terminology dialog as the previous information will be deleted. This is useful when updating multiple workstations to ensure all workstations have the same information.

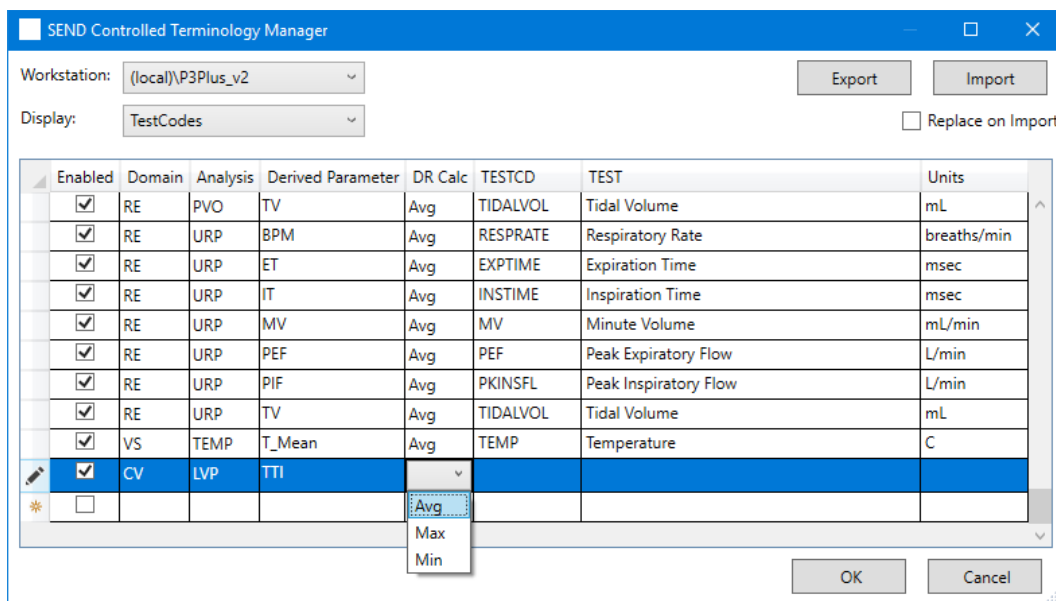


Figure 34. – Modification of Test Code calculation.

Figure 35 shows the EG Leads currently configured. All available options under the Display pull-down menu allow modification as described above.

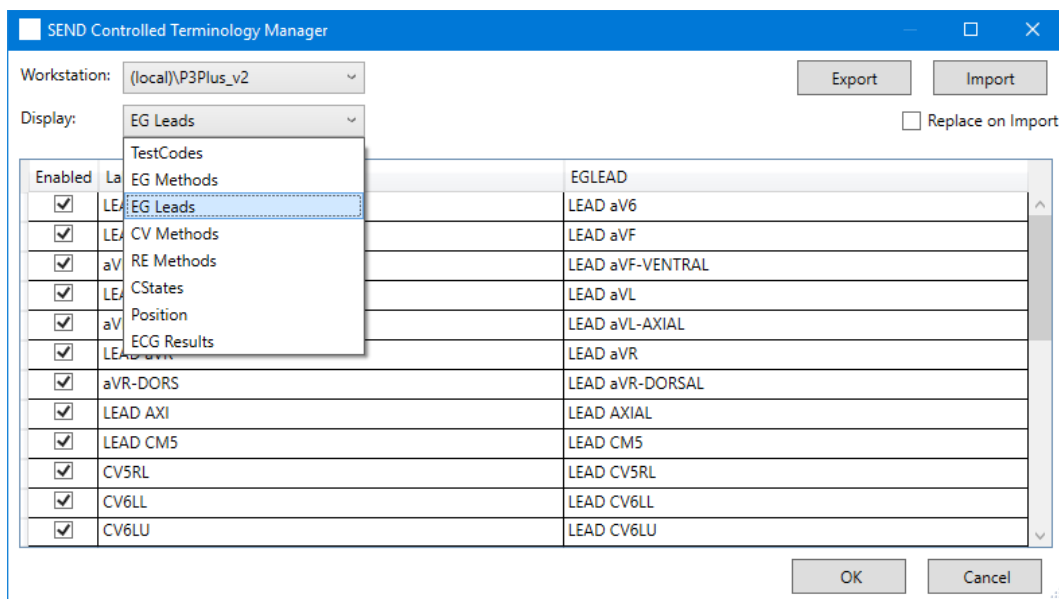


Figure 35. – EG Leads list shown.

SYNCHRONIZE SEND DATA

Send information can be synchronized to a dedicated server or workstation. There are two types of synchronization, Synchronize SEND Data and Synchronize SEND Study Data.

NOTE: It is important to note that this functionality is only available if the SEND SQL Server Option has been enabled in the license file. If only SEND Output is enabled, these features will not be available.

SEND REMOTE SERVER SYNCHRONIZATION SETUP

The following instructions require the Ponemah SEND v3.1 module and apply to Ponemah v5.31 and greater.


1 – Create a SQL Server Instance

Consult your local IT group for assistance configuring the instance required for use with SEND. On the server, configure an instance name P3Plus_v2. Add a temporary SQL user with the following rights. Do not use P3StudyUsers as the name of the temporary SQL user.

- dbcreator
- securityadmin
- Set password to never expire
- Do not require a password change on the first logon

2 – Central Server Utility

Please note that the example below uses Ponemah v5.31 but this utility can be used with Ponemah versions 5.31 and greater. Substitute v531 with the Ponemah version being used, example 532, 540, etc.

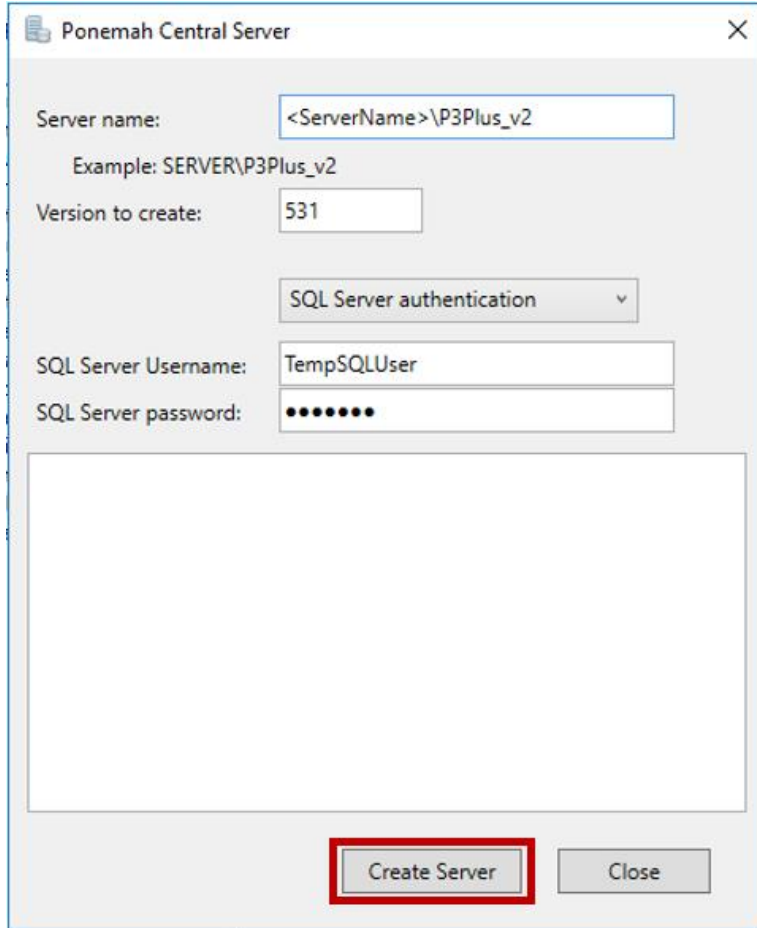

The screenshot shows a window titled "Ponemah Central Server" with the following fields and controls:

- Server name: (Example: SERVER\P3Plus_v2)
- Version to create:
- SQL Server authentication:
- SQL Server Username:
- SQL Server password:
- Buttons: "Create Server" (highlighted with a red box) and "Close"

Figure 36. – Output Server configuration using the Central Server Utility.

3– Run Ponemah.CentralServer.exe

The Ponemah Central Server utility can be found at C:\Ponemah\Utilities\Ponemah.CentralServer\. Enter the following information:

- Sever name:
 - <ServerName>\P3Plus_v2
 - <ServerName>\P3Plus_v2<port> (if a specific port is being used)
- Version to create:
 - Use the version of the Ponemah software installed, example, 531, 532, 540, etc.
 - Select SQL Server authentication
 - SQL Server Username: Name of the temporary SQL account
 - SQL Server password: Password of the temporary SQL account

- Select Create Server
- Note any messages/errors that are listed in the application
- The following actions will be taken
 - A P3StudyUsers account will be setup by the application with dbcreator rights
 - P3Plus_531 database (or appropriate version) will be created.
 - P3Studies database and tables will be created

4 – Create SEND_31 Database

Launch Ponemah and select P3 Setup from the Setup pull-down menu. In the tree view on the left hand side, select SEND and enter the server name in the Output Server edit field and select Test Connection

Figure 37. – SEND configuration tab.

Enter the Server name or dedicated workstation name and use the Test Connection button to verify that the connection is valid. Along with the server name, the SQL Instance (P3Plus_v2) must also be defined, example “server name\P3Plus_v2”. If successful, a message will post stating that the connection to the SEND database was successful.

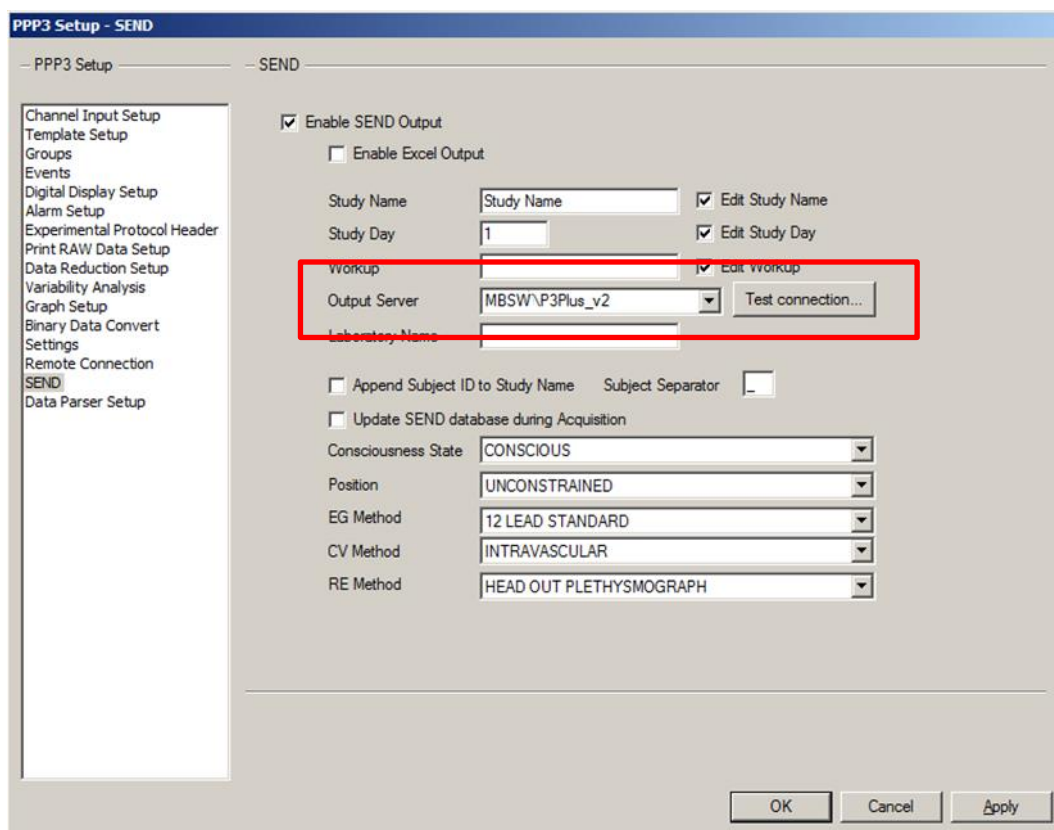


Figure 38. – Output Server configuration in SEND dialog.

If no Output Server has been configured, attempting to synchronize study data will post a message requesting that the server be configured in P3 Setup.

SYNCHRONIZE SEND DATA

The Synchronize SEND Data feature allows SEND data contained within SQL on a workstation to be synchronized with another workstation, or server, to create a single repository for SEND data. This requires configuration of the Output Server in the SEND setup dialog.

NOTE: It is important to note that when performing a sync within Review with a subset of subjects loaded, all subjects associated with the Review file will be synced. In addition, if only a single channel from a subject is loaded in review, saved, and a sync performed, all existing SEND data will be replaced for that subject in the current review file.

After configuring the Output Server, navigate to the Synchronize SEND Data as shown in Figure 39. The SEND data for the currently loaded Review file will be automatically synchronized to the SEND database on the Output Server (Synchronize SEND Study Data is not available within a Review session). Only data from the current Review session will be synchronized to the Output Server when in Review.

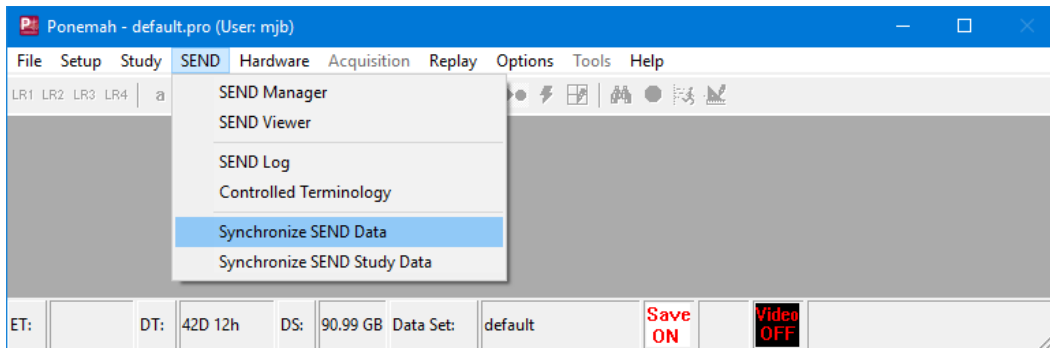


Figure 39. – Synchronize SEND Data dialog selection from SEND pull-down menu.

The Synchronize Study Data function can also be used when outside of a Review session. Figure 40 shows the pop-up dialog that allows selection of the Study and Output Server. Clicking the OK button will synchronize the selected study to the designated Output Server defined. This will include all SEND data associated with that study, whereas when in a Review session, only the data for the currently loaded Review file will be synchronized.

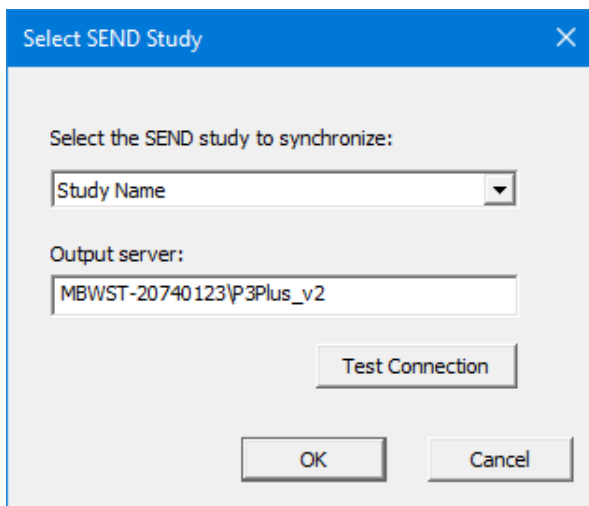


Figure 40. – Synchronize SEND Data dialog outside of Review session.

SYNCHRONIZE SEND STUDY DATA

The Synchronize SEND Study Data will synchronize from all workstations connected and visible on the network to the workstation performing the sync and to the Output Server. Both the workstation performing the sync and the Output server will then contain the same SEND information.

The Output Server defined in the SEND dialog must be part of the Study configuration, otherwise a message will post stating that they sync cannot be performed to the Output Server.

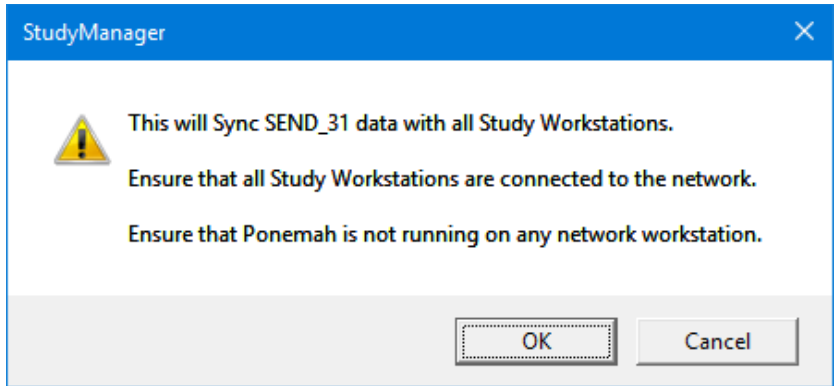


Figure 41. – Synchronize SEND Study Data notification.

The Synchronize SEND Study Data function is only available in idle mode and will not be available while in a Review session.

When performing a sync, the Output Server will be locked out from other workstations until the current sync has been completed. If an attempt to synchronize to the Output Sever is initiated by a second workstation, a message will be posted that the server (workstation) is in use and attempts to synchronize the data should be retried at a later time. In addition, error checking is performed during the synchronization. If conflicts between the data on the Output Server and the workstation performing the synch, Conflict Resolution dialog will be presented to allow the appropriate information to be updated.

Below, figure 42, is an example of the conflict resolution dialog. The Options fly out menu in the upper left hand portion of the dialog will allow you to expand or reduce the information shown. The dialog provides two columns, one column showing data from the workstation performing the sync and a second column showing a server (or second workstation) where data is being synched. In the example, both the workstation and server contain SEND data for these two subjects.

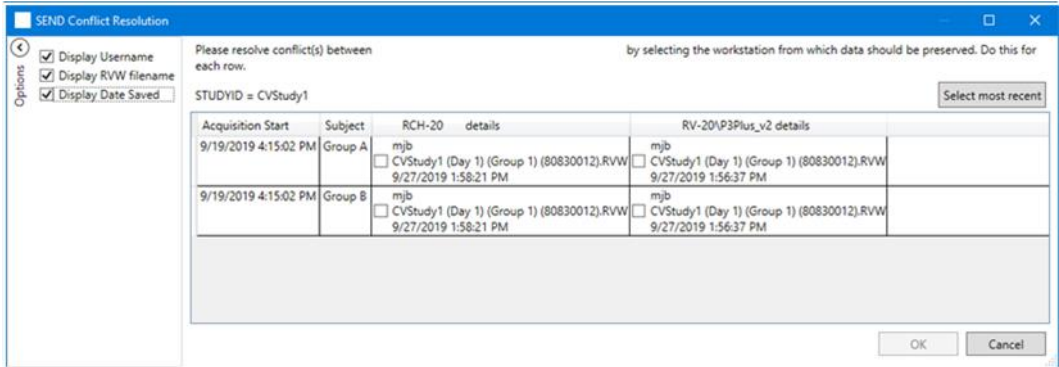


Figure 42. – Conflict Resolution dialog.

Since the information that exists in both SEND databases does not match, the Conflict Resolution dialog is posted allowing the user to select the most appropriate data for each animal. In this example, data was reanalyzed and new values for the derived data were created. Figure 43 shows the Run information along with the time that data was saved. Time 1:58:21pm is the newly analyzed data. The option to select the most recent data, or previous data, is available per subject. Once selected, the cell is highlighted in green to easily see which runs will be synchronized to the SEND databases of each.

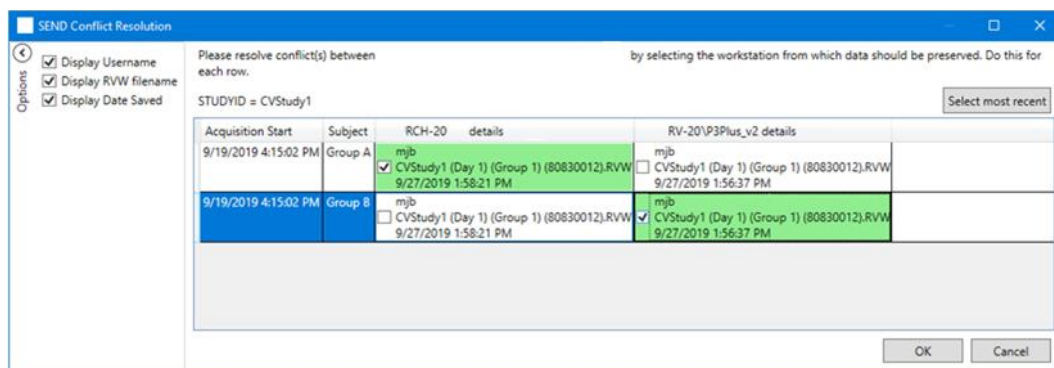


Figure 43. – Conflict Resolution dialog.

The Select most recent button (Figure 44) instructs the software to select the most recent information for each subject. Again, all highlighted subjects will be show in green.

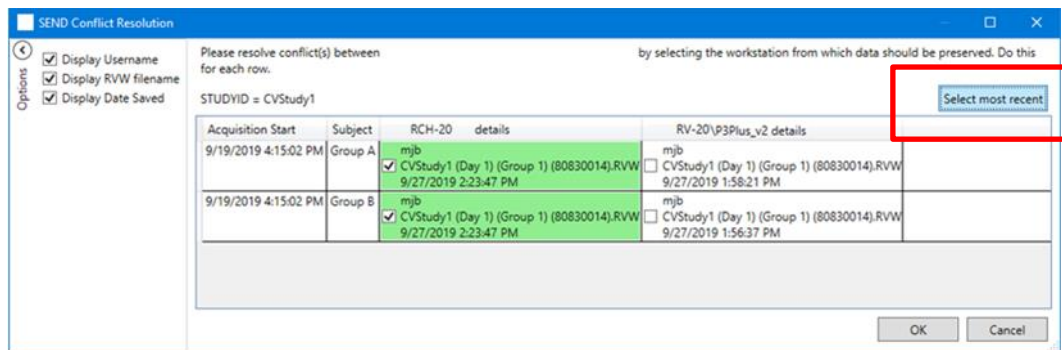


Figure 44. – Conflict Resolution dialog, most recent information highlighted.

HIGH LEVEL PROCESS FLOW FOR GENERATING SEND OUTPUT

The previous sections have described how to configure and generate SEND output. This section will provide two scenarios and the high-level process flow needed to generate SEND output. No detailed configuration information is provided below, refer to previous sections for additional information.

While the following two examples provides guidance on creating SEND output, which includes configuration information, once a SEND configuration has been completed and saved in the protocol, only saving of the derived data is necessary to generate SEND output.

Generating SEND Output to Excel (outside of the Study Protocol Option)

- Ensure SEND Output or SEND SQL Server Output option is enabled in the license file
- Open a Review file (.RVW)
- Configure Data Reduction, if not previously configured, with CDISC defined variables
 - Ensure the event to trigger Data Reduction has been initiated
- Update Groups tab with appropriate Name and Dose information for each subject
- Enable the SEND dialog under P3 Setup and update the appropriate fields
 - Select the checkbox for Enable Excel Output
- Perform a Save Marks Section or Saved Derived Data to generate the Data Reduction output to SEND
- Navigate to the Ponemah_Data directory or the location where the Excel file was saved
 - Open Excel workbook and verify that the appropriate worksheets exist for the Domains used

Generating SEND Output to CSV using the Study Protocol Option

- Ensure SEND SQL Server Output option is enabled in the license file in order to generate CSV files
- Select the desired Study from the Manage Studies dialog
- Enter the Run Study dialog and select the desired Run from the Previous Runs list
- Data Reduction should have been previously configured and saved as part of the Study configuration (protocol file) with CDISC defined variables. Modification of Data Reduction is not permitted after data collection within the Study Protocol Option
 - Ensure the event to trigger Data Reduction has been initiated
- Groups should have been configured with the appropriate Name and Dose information for each subject as part of the Study configuration.
 - Select Use Study Subjects and Use Study Doses to automatically pull information from the Study configuration.
 - Uncheck Use Study Subjects and Use Study Doses if edits are necessary to correct entry errors from Study configuration
- Enable the SEND dialog under P3 setup and updated the appropriate fields if not previously saved in the protocol
 - Uncheck the following to allow previously defined Study information to be pulled into SEND automatically
 - Edit Study Name
 - Edit Study Day
 - Edit Study Workup
 - If desired, uncheck these edit boxes to allow for manual edits to better define the output being generated or correct for entry errors when creating the Study
- Perform a Save Run (Marks Section) to generate the Data Reduction Output to SEND
- Navigate to the Send Viewer and select the desired Study to generate a CSV file output
 - Select the appropriate Domain, Filters, and Variables
 - Use the Export to CSV button to generate the SEND output file
- Optionally, use the Synchronize SEND Data and Synchronize SEND Study Data to synchronize data across workstations or servers

TROUBLESHOOTING

This section outlines common messages encountered with incorrect settings when SEND is enabled. Each item will list the issue or message that may be posted along with a resolution.

MESSAGES POSTED DUE TO INCORRECT SEND CONFIGURATION

1 – Incorrect ECG labels

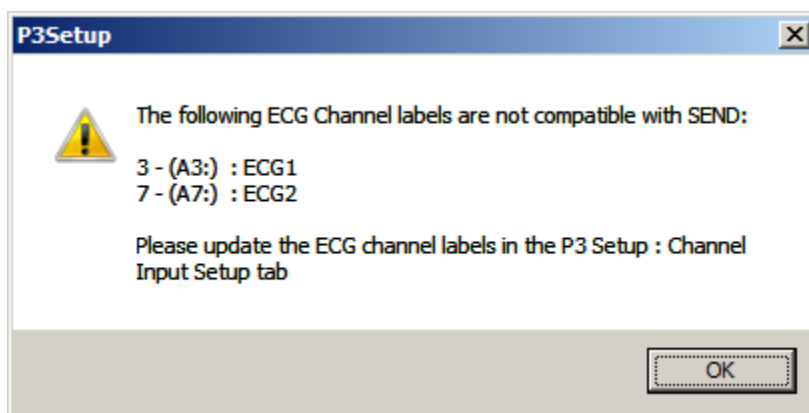


Figure 45. – Incorrect ECG labels in Channel Input Setup.

Issue: CDISC dictates specific labeling terminology for ECG lead presentations. If the label entered in the Channel Input Setup menu does not match that used by CDISC, the above message will be posted.

Resolution: Enter the Channel Input Setup menu and update the labeling for the ECG channels identified in the message posted. When you select the Label column, you will have a drop-down list that allows you to select the correct lead presentation.

2 – Duplicate TESTCDs

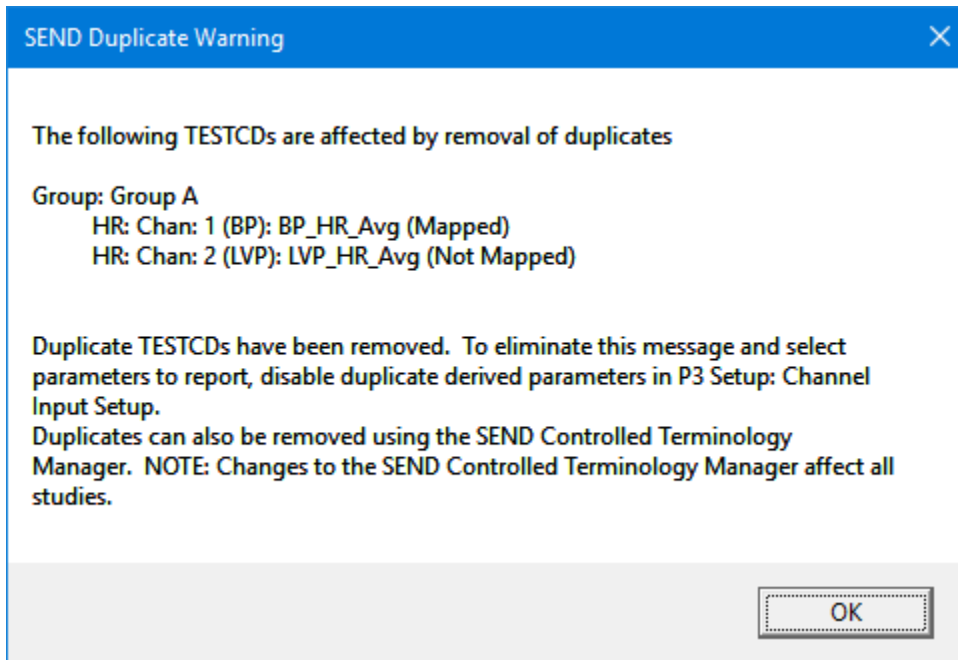


Figure 46. – Duplicate TESTCDs message.

Issue: A CDISC TESTCD can only be reported once for a subject in a given interval in a single domain. The above error will be displayed if duplicate derived parameters exist across multiple analysis algorithms. In the example above, the HR (Heart Rate) parameter is enabled in both the Blood Pressure and Left Ventricular Pressure analysis algorithms.

Resolution: Enter the Channel Input Setup menu and update the derived parameters list for the Subject(s) listed in the warning message. Resave the file to generate the desired output. If you say OK to the message, the software automatically uses the TESTCD (in this case HR) from the first channel where the conflict was identified. In this case, the hardware configuration had BP pressure as channel 1 and LV defined as channel 2. The HR parameter was reported for the BP channel and removed from the LV channel.

3– NO SEND data in Excel or SEND Viewer


Issue: Data not present in the Excel file or in SEND Viewer after performing a save of the Marks Section/Run or Derived Data.

Resolution: Several reasons may exist which will result in no data available in the output file .

- SEND has not been enabled in P3 Setup
 - If using SEND to Excel, ensure checkbox in SEND dialog has been enabled
- Data Reduction has not been configured

- CDISC TESTCDs are not enabled in Data Reduction. Not all derived parameters (Data Reduction) are currently defined by CDISC
- Ensure the Event to trigger Data Reduction has been placed
- If SEND does not appear in your P3 Setup menu, the current Ponemah license file likely does not have the SEND option enabled. This information can be found under the View License File dialog from the Help pull-down menu. These features will be listed under the Options box. The SEND pull-down menu while in idle mode is only available (accessible) if the SEND option is enabled. Contact Technical Support for assistance.

4 – No Output Sever Defined

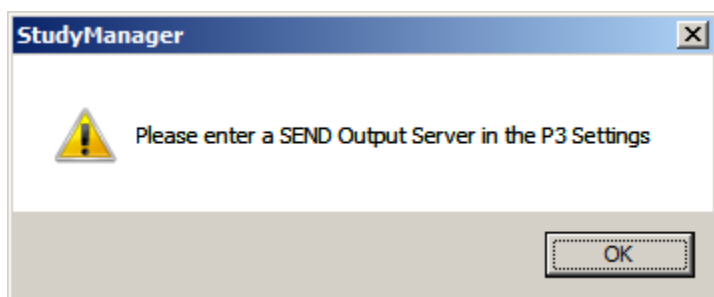

Or

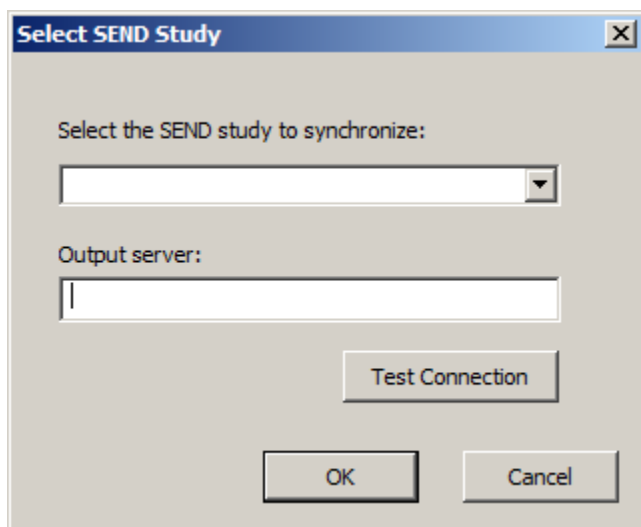


Figure 47. – Incorrect Output Server configuration.

Issue: No Output Sever has been configured

Resolution: The first message will post when in an active Review session and the second message when in idle mode. Enter the appropriate server/workstation in the SEND dialog under P3 Setup. Ensure that the Ponemah Instance has been included with the server/workstation name.

5 – No subject ID for USUBJID TESTCD

Issue: No user defined subject ID present in the SEND output for USUBJID.

Resolution: When not using the Study Protocol Option and pulling in Subject information automatically, the default label (Group A, Group B, etc.) will be placed in the SEND output. From P3 Setup, enter the Groups dialog and modify the Name column. Perform a Save Derived Data or Save Marks Section to update.

6 – No Data in SEND Viewer

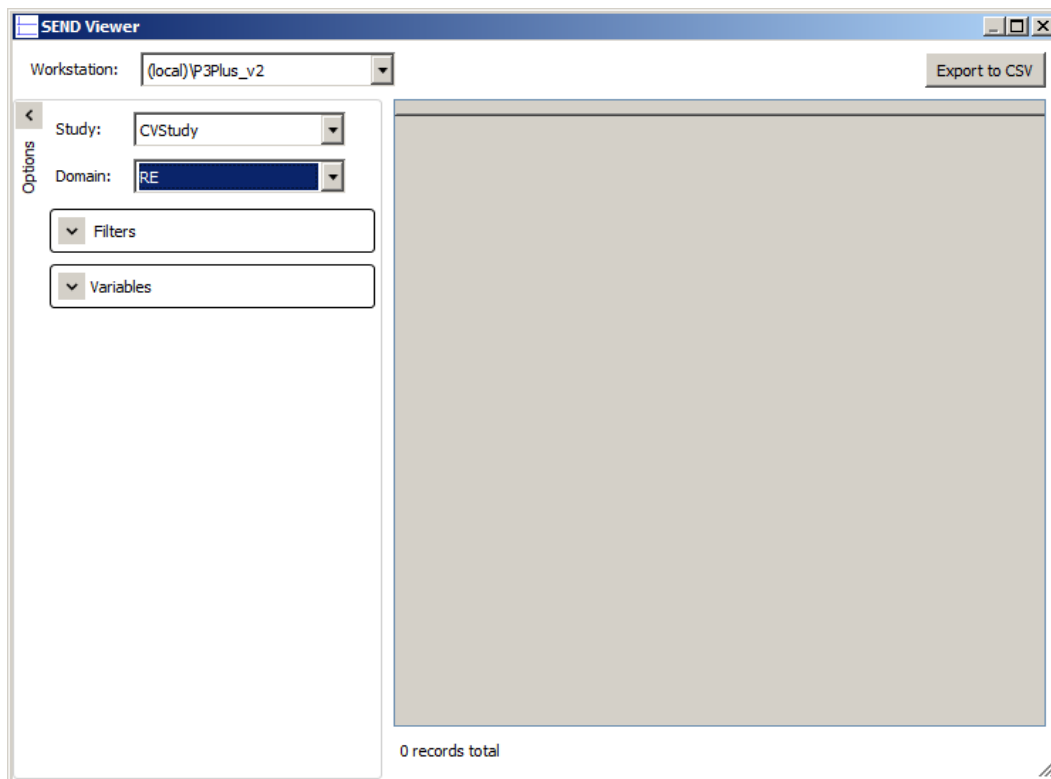


Figure 48. – No SEND data currently visible in SEND Viewer.

Issue: No data is present when opening the SEND Viewer dialog for a given study when all information has been configured properly (SEND, Data Reduction, etc.) and a save of the data has been performed.

Resolution: Ensure that the correct Domain is listed for the study that has been selected. Data currently exists for the CVStudy example shown above. However, no data in the RE Domain exists. Therefore, no data will populate the window. Select the correct Domain, EG and CV in this example, and the SEND information will be displayed.

TECHNICAL SUPPORT

DSI™ is available to help you with your questions and concerns. Should you hit a road block or need some additional training, please feel free to contact us. We are happy to help!

DSI TECHNICAL SUPPORT— NORTH AMERICA

Email: Support@datasci.com

Toll-free in U.S. and Canada

Phone: 1-800-262-9687

Monday through Friday: 8 AM to 5 PM CST
(except Holidays)

DSI TECHNICAL SUPPORT— EUROPE

Email: Europe-support@datasci.com

Phone: +44 1359 259400

Monday through Friday: 8 AM to 5 PM CET

DSI TECHNICAL SUPPORT— ALL OTHER COUNTRIES

Phone: +1-651-481-7400

APPENDIX 1

Not all derived parameters calculated by Ponemah are currently defined by CDISC. Below lists the currently accepted parameters by general signal type. Please note that derived parameters may exist in multiple Analysis algorithms. For example, the derived parameter “Sys” is reported by both the Blood Pressure (BP) and Left Ventricular Pressure (LVP) analysis algorithms.

In addition to enabling only the CDISC defined parameters, Data Reduction calculations selected must also be defined by CDISC or no SEND output will be generated. The accepted Data Reductions calculations are Average (AVG), Maximum (Max), and Minimum (Min).

ECG

- Heart Rate
- P-H
- PWdth
- PP-I
- PR-I
- QRS
- QT-I

- QTcb
- QTcf
- QTcv
- QTca
- QTcl
- QTcu
- R-H
- RR-I
- ST-I
- ST-E
- T-H
- T-A
- Tpe-I

Activity

- A_TA

Pressure (includes analysis modules BP and LVP)

- Dia
- +dP/dt
- -dP/dt
- HR
- LVEDP
- Sys
- Mean
- PH
- Q-A
- TPR

Respiratory (includes analysis modules dEMG, PAF, BPR, PAP, PCR/PCRP, PVO and URP)

- TV
- Res
- RT
- ET
- IT
- MV
- Penh
- Max (PCRP analysis module)
- Min (PCRP analysis module)
- PIF
-
- Cdyn
- PEF
- BPM
- RBpm

- DIA (PAP analysis module)
- Mean (PAP analysis module)
- SYS (PAP analysis module)


Temperature

- Mean

APPENDIX 2

The following is an example of a Multi-lead ECG configuration created using Study with Data Reduction output defined by Parser Segments. The protocol, subjects, and doses have all been created. In this example, Workups will be defined for both pre-dose and post-dose collection times.


When Workups are configured in the appropriate format, the workup label will replace the Data Reduction label and be included in the SEND output. Also, the assumption when configuring a Tox Study design is that there will be only one interval reported for an acquisition, unlike a telemetry collection where multiple Data Reduction intervals will be reported over an extended collection period.

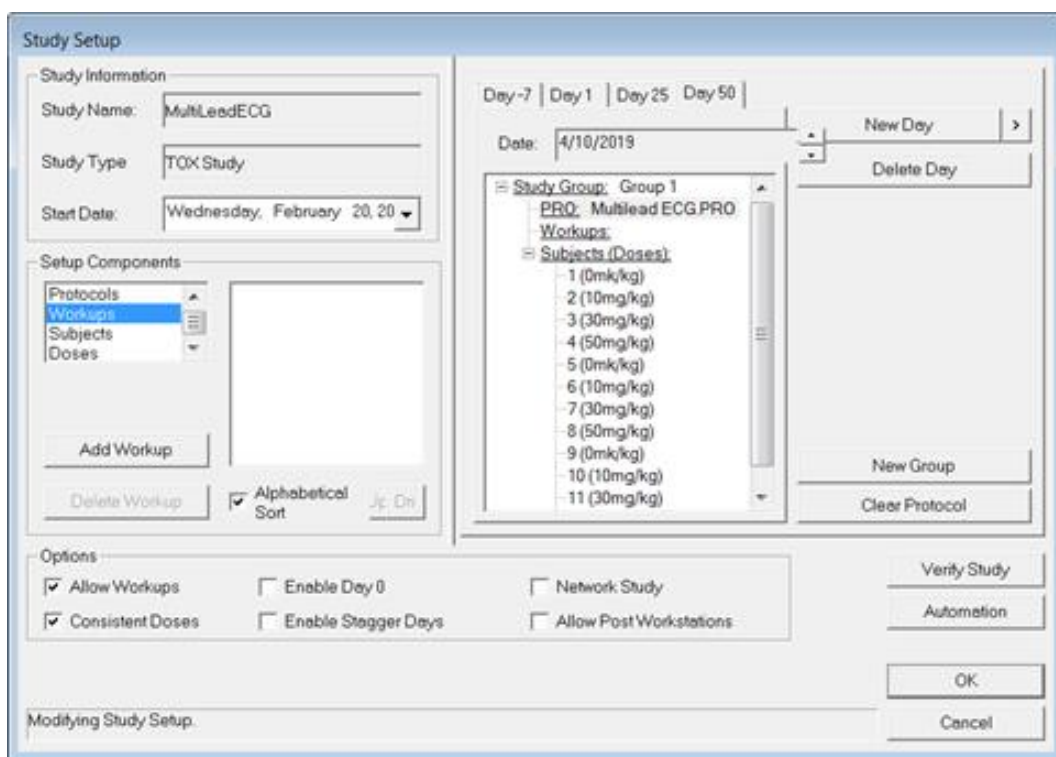


Study allows any text to be entered and used within this study context. However, only SEND compliant terminology (ISO 8601 interval format) will be able to be pulled into the SEND output. If the workups defined here do not meet the accepted format, acquisition and review sessions will still function as expected, however, the workup label entered above will not appear in the SEND tables as entered.

To define a Workup, click on the Add Workup button. The following dialog will be displayed. To define a two hour post-dose time, enter the time in hh:mm:ss format. Note that zeros used for padding can be dropped as they are not needed, example 04H would simply be 4H.

To remain compliant, a two-hour post-dose collection would include "PT" in the string. P represents the start of the duration and T represents time. The designators must precede the actual time. Therefore, the SEND compliant time point would be entered as "PT2H".

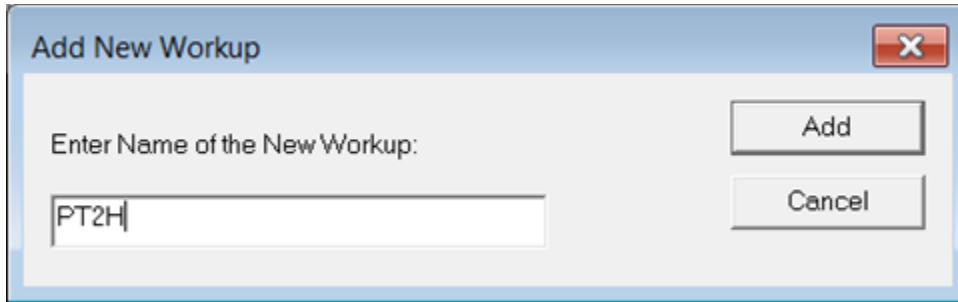


Additional timepoints have been entered to represent a two-hour pre-dose timepoint and a post-dose timepoint of 4:30:15 seconds.

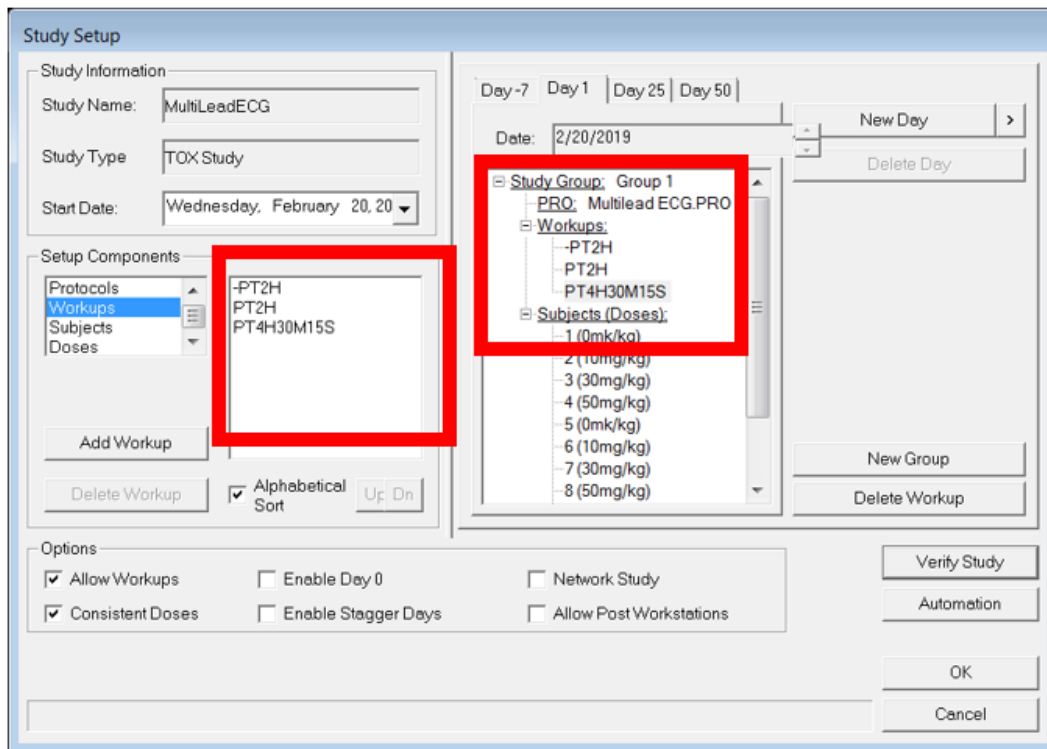


Data Review and SEND Configuration

Below shows the two-hour pre-dose Review session with a single parser segment and the AVG (average) calculated for the enabled derived parameters. It is also worth noting that derived parameters will only be present in SEND output if those parameters have been mapped to a TESTCD in the Controlled Terminology menu and enabled in the Channel Input Setup menu. By default the Controlled Terminology menu only contains mappings for results that are defined by CDISC. Ponemah provides a number of calculations that are not currently defined by CDISC.

	Time	Date	Time	Time	Time	8: RR-I	8: HR	8: QRS	8: PR-I	8: QT-I	8: QTcb
T0	0000:00:00	14:46:16	02/06/2019								
2h Predose Avg	0000:00:02	14:46:18	02/06/2019	0000:00:05	0000:00:05	999.80	60.012	74.667	171.50	349.83	349.87

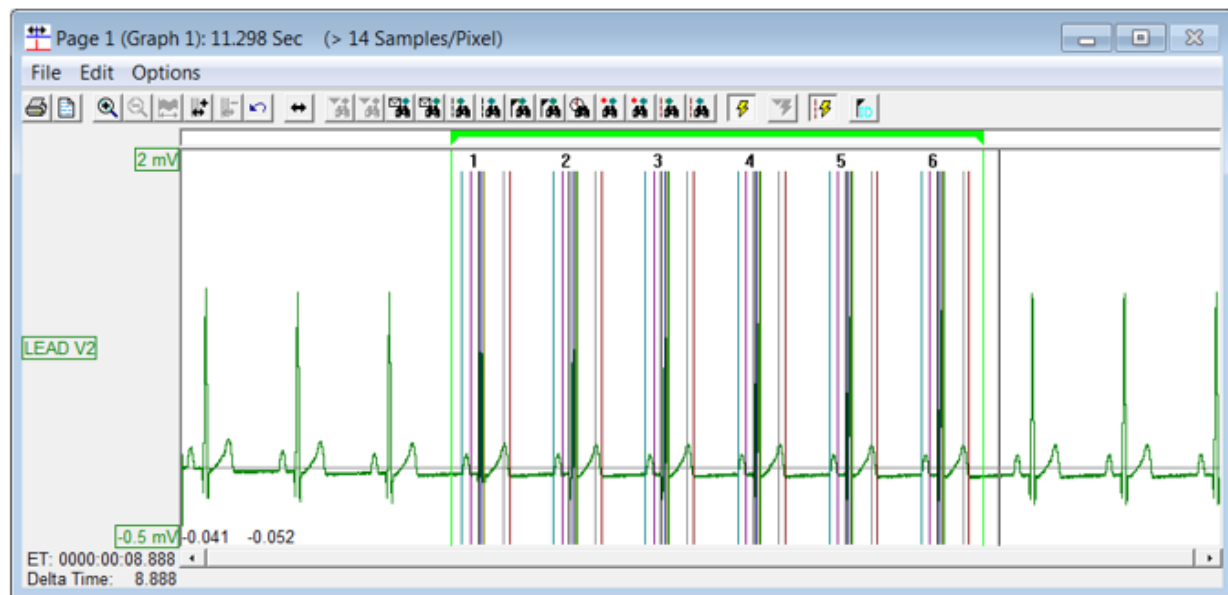

SEND information can be configured in idle mode along with other study and setup information. Below is an example of the SEND dialog after configuration. To utilize the automated features with Study, uncheck Edit Study Name, Edit Study Day, and Edit Workup. This will pull in the Study information based on the file that is currently loaded. In this example, it is the 4:30:15 timepoint.

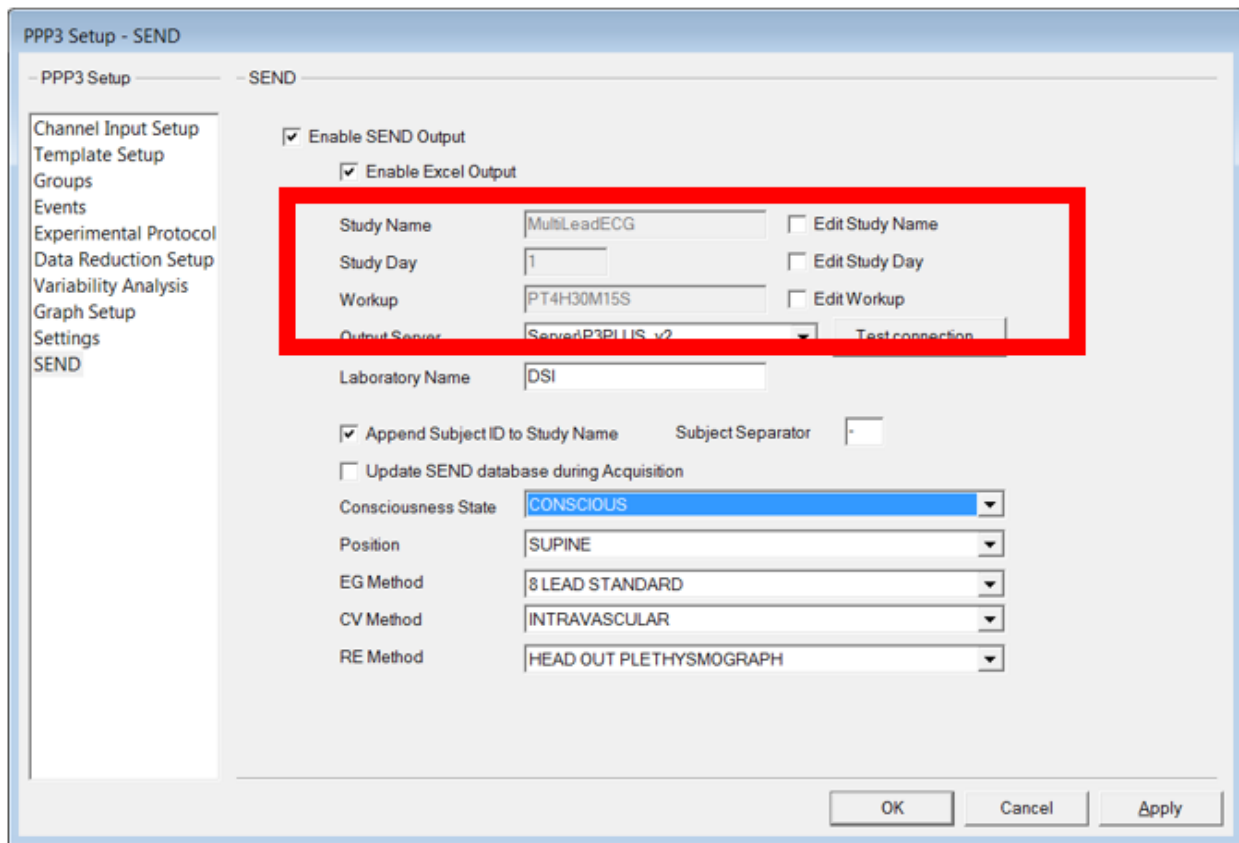

SEND Output

From the SEND pull-down menu on the main Ponemah window, select SEND Viewer. On the left side of the dialog, select the filters specific to the study and data of interest.

Workup information entered into Study is included in the SEND output under EGTPPT and EGELTM.

SEND Viewer

Workstation: (local)\P3Plus_v2 Export to CSV

Options

Study: MultiLeadECG

Domain: EG

Filters

Variables

iENDTC	EGDY	EGE	EGDY	EGNOMDY	EGTPT	EGTPTNUM	EGELTM	EGTPTREF	EG	TT
19-02-06T14:46:24				1	2h Predose	1	-PT2H	Day 1 Dose 1:0mk/kg		
19-02-06T14:46:24				1	2h Predose	1	-PT2H	Day 1 Dose 1:0mk/kg		
19-02-06T14:46:24				1	2h Predose	1	-PT2H	Day 1 Dose 1:0mk/kg		
19-02-06T14:46:24				1	2h Predose	1	-PT2H	Day 1 Dose 1:0mk/kg		
19-02-06T14:46:24				1	2h Predose	1	-PT2H	Day 1 Dose 1:0mk/kg		
19-02-06T14:46:56				1	2h Postdose	2	PT2H	Day 1 Dose 1:0mk/kg		
19-02-06T14:46:56				1	2h Postdose	2	PT2H	Day 1 Dose 1:0mk/kg		
19-02-06T14:46:56				1	2h Postdose	2	PT2H	Day 1 Dose 1:0mk/kg		
19-02-06T14:46:56				1	2h Postdose	2	PT2H	Day 1 Dose 1:0mk/kg		
19-02-06T14:46:56				1	2h Postdose	2	PT2H	Day 1 Dose 1:0mk/kg		
19-02-06T14:47:11				1	4h 30m 15s Postdose	3	PT4H30M15S	Day 1 Dose 1:0mk/kg		
19-02-06T14:47:11				1	4h 30m 15s Postdose	3	PT4H30M15S	Day 1 Dose 1:0mk/kg		
19-02-06T14:47:11				1	4h 30m 15s Postdose	3	PT4H30M15S	Day 1 Dose 1:0mk/kg		
19-02-06T14:47:11				1	4h 30m 15s Postdose	3	PT4H30M15S	Day 1 Dose 1:0mk/kg		
19-02-06T14:47:11				1	4h 30m 15s Postdose	3	PT4H30M15S	Day 1 Dose 1:0mk/kg		
19-02-06T14:47:11				1	4h 30m 15s Postdose	3	PT4H30M15S	Day 1 Dose 1:0mk/kg		

18 records total