


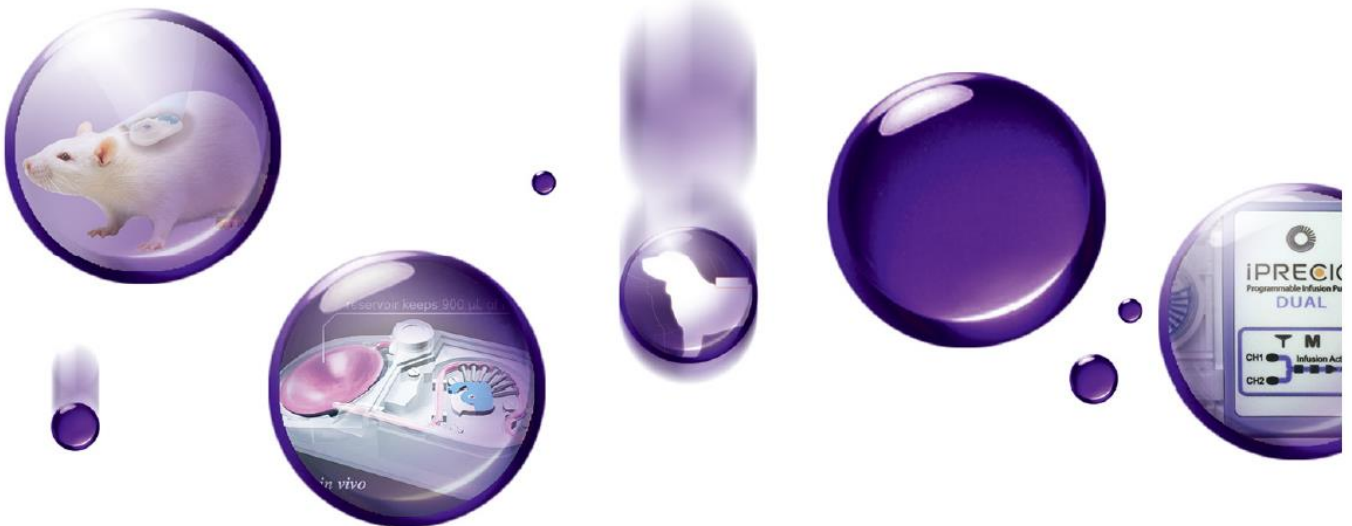

iPRECIO[®]
Micro Infusion Pump


iPRECIO Management Software Release Note

Version 1.3 Rev.532renew Release Notes
August 31, 2020



Version 1.3 Rev.532 renew

- 1) Application software updated to be able to read SMP-200 pumps manufactured in 2020.
 - Pump SN from 2010001 can only be read correctly using Version 1.3 Rev.532 renew
 - Installation Instructions for Version 1.3 Rev.532 renew (includes information on driver install)
- 2) Fixed: Fixed a bug that in the Infusion Protocol Setup window, enter the date of the time beyond 2020/12/31 23.59 in Set Date / Time and click [Next>>] to set Instant Infusion Mode / Post Recovery Mode. an error message shown below appears, and the protocol cannot be set up.

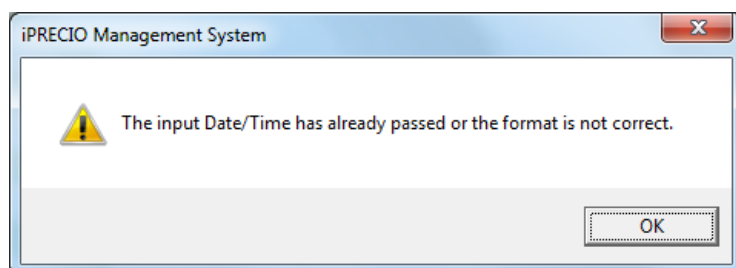

※ After this Rev.532 renew, bug fixes and improvements will be made only for Windows 10 compatible versions.

Device Driver Information

- 1) Update of USB ID on UCD-200
 - From Serial Number (SN) 02421 (June 18 2020)
 - Signed drivers will automatically install on Windows 10 PCs with internet connection
 - Drivers are also on CD for manual installation for PCs without internet connection (Folder: Drivers 2020)

For SN before 02421, USB ID on UCD-200 will required workaround to install drivers on Windows 10 PCs

- Drivers are installed during Test Signing Mode
- Detailed instructions are provided – Driver Install on Windows 10 PC (2019-12-23)
- Drivers are also on CD (Folder: Drivers)

Instructions – Administrator rights required to install and uninstall Application Software.

- 1. Make copies of important ipf and csv files from ongoing and past studies for safety before installing and uninstalling IMS-200 Application.**
2. Uninstall current version of IMS-200 Application Using Windows 10 <Apps and Features> before installing new version
3. Download full folder and copy to CD or hard disk of PC
4. Follow installation instructions

iPRECIO install CD ver13rev532 revised (August 2020) – Zipped

<https://drive.google.com/drive/folders/1Y69qouiMLQtBi4FJf9Eld8QcG19aEgMy?usp=sharing>

iPRECIO install CD ver13rev532 revised (August 2020)

https://drive.google.com/drive/folders/15naU5hf-p_pOTFyFrnBaGS2dbSNFxFrjX?usp=sharing

Past version history

Version1.3 Rev.532 (only released in Japan)

Fixed: The bug that flushing was not executed properly and subsequent administration was not performed when Total Dead Volume was set to 59.650 µl or more using "Recovery Mode" was fixed.

Version1.3 Rev.531 (Only released Internationally)

Fixed: Fixed the bug that when using up to 10 steps in "Variable Flow Mode", the error message that appears when the dosing schedule exceeds the battery level is displayed on the monitoring screen even after the software automatically corrects the schedule.

Fixed a bug that concentration units of chemicals were not displayed correctly in the exported data file.

Fixed a bug that the refill date and time notation (yy/mm/dd) is displayed incorrectly as (dd/mm/yy) in the exported eta file.

Other changes:

The folder structure in the installation disk has been changed.

The installer was fixed while unifying the folders that were independent for each OS.

Primetech Corporation, Japan. September 10th. 2020.