




## iPRECIO IMS-200 Application Software Version 1.3 Revision 532 Renew

- DISK1
- DISK2
- DISK3
- DISK4
- Documents for Returns
- Driver
- iPRECIO Workflow Examples
- Surgical Videos (links to download videos)
- Technical Notes
- User Manual - FAQ
- AUTORUN
- FirstStep
- Important Notice-Total Dead Volume
- Installation Instructions for iPRECIO install CD 13rev532 (Feb 2020)
- Readme (for Serial Number information)


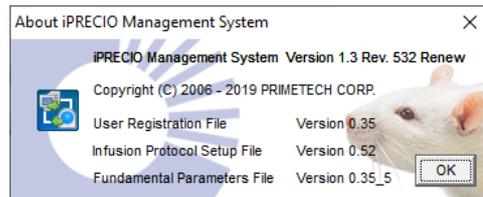


### From Install CD

1. Run <FirstStep> as Administrator (Administrator Rights Requested) for installation. Screen below <iPRECIO Management System Setup> will appear.
2. Plug UCD-200 device to PC – Required drivers will be installed (internet access required)
  - a. Driver Install Option shown below is no longer required as Windows 10 will automatically install the required drivers when UCD-200 device is plugged to PC USB port.
  - b. After completed, Run <Install> to start the installation process
3. Then follow instructions on the PC/Screen

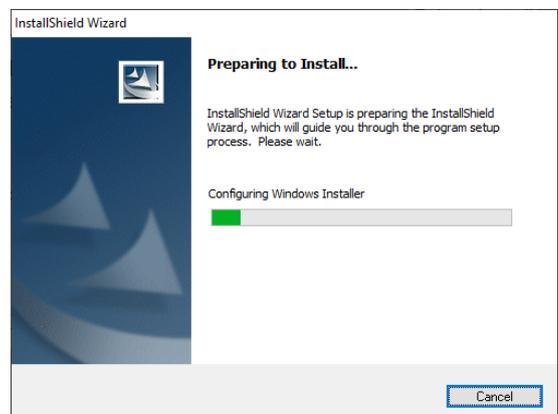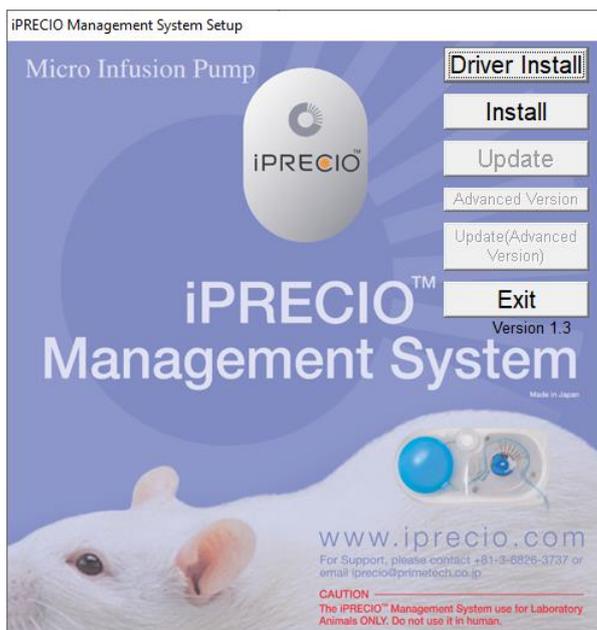


Install Shield Wizard appears, Click <Next> and then accept License Agreement and click <Next>

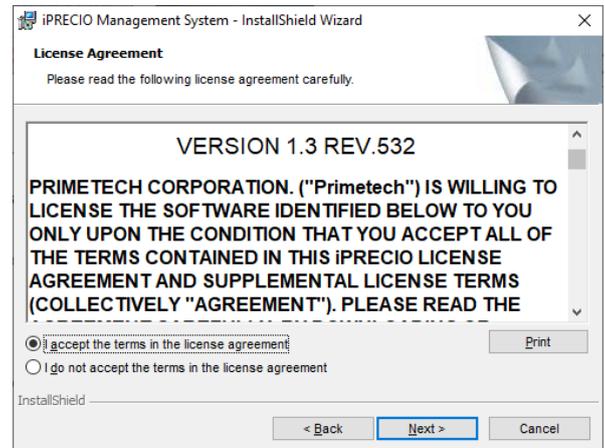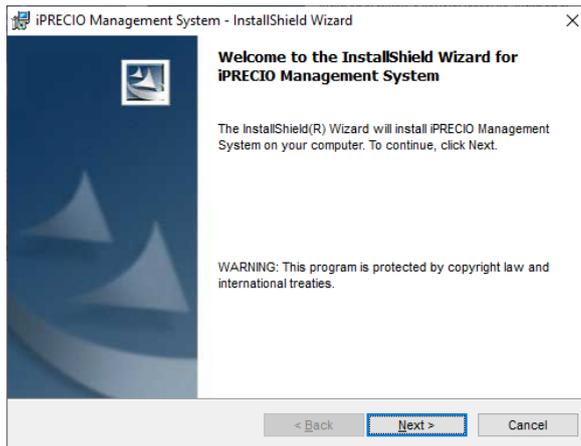


You will need to use the software key from the CD included in the package.

- The iPRECIO Software CD is labeled with serial number (S/N),
- IMSV13R532XXX where XXX is a unique number. Screenshot does not show XXX and only 532
- Refer to iPRECIO User Manual for full installation instructions. (User Manual Version is 1.3rev531)
- See page 14 for S/N details and page 13 for full installation guide. (User Manual Version is 1.3rev531)
- **Latest version is 1.3rev532**



Fill as required and click on <Next>

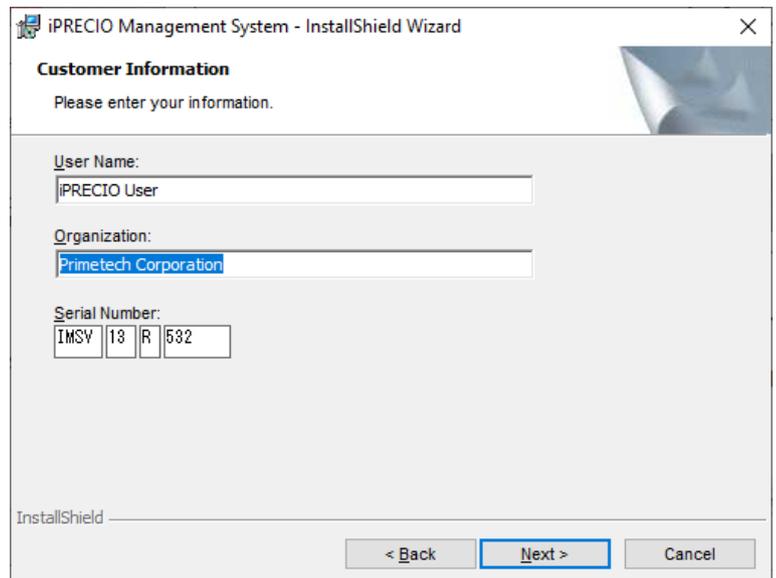

Confirm or change folder location and click on <Next> to proceed.

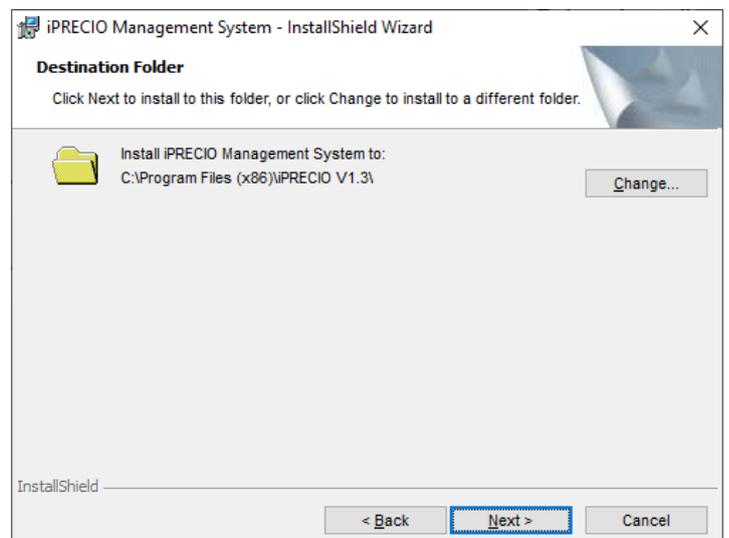


Click on <Install>

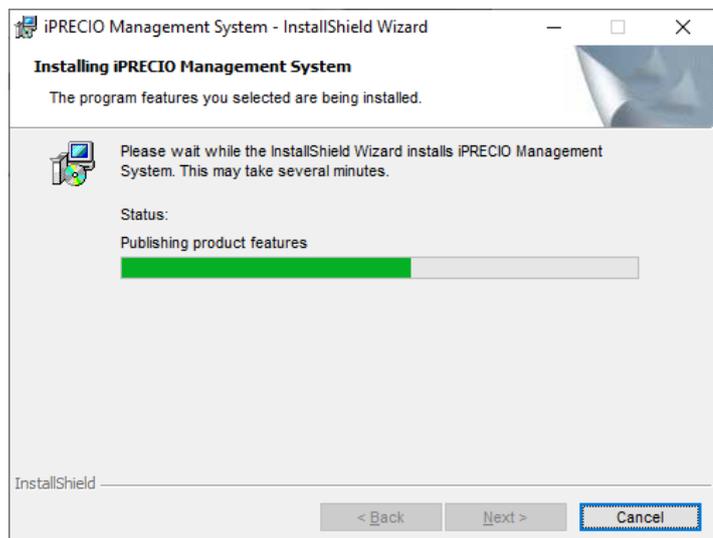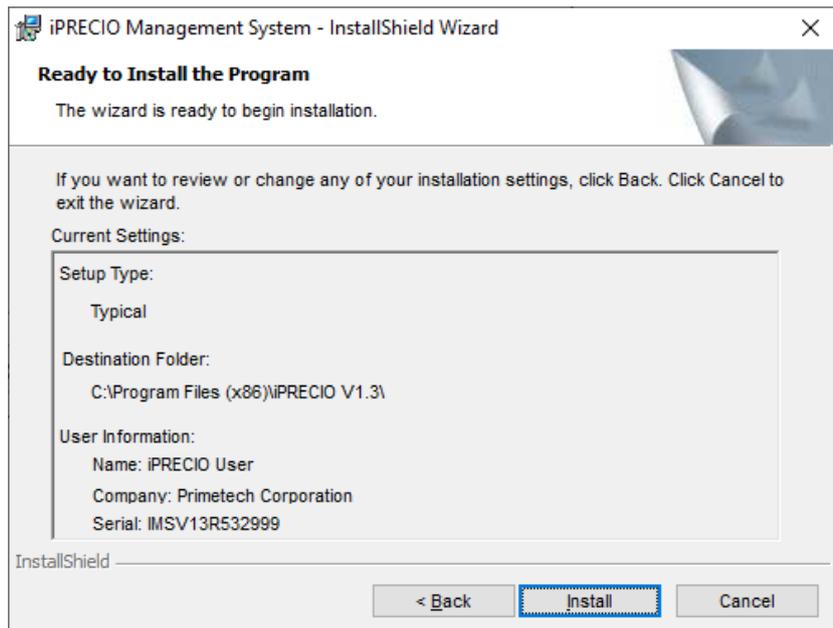

Click on <Finish> and close Setup Window as installation is complete.

If required setup shortcut to desktop.

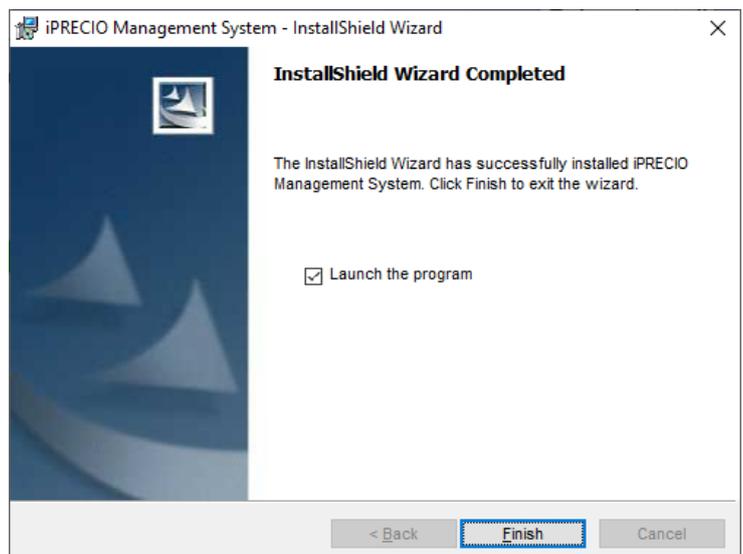


Run Application and See Page 16 of User Manual for software installation.

**Software Setup – Page 16 of User Manual**



The 'User Registration' dialog box contains the following elements:

- Title bar: 'User Registration' with a close button (X).
- Section Header: 'User Registration'.
- Form Fields:
  - Institution Name (\*)
  - Department Name (\*)
- Legend: (\*) Required field.
- Instructional Text: 'Director's Name, Login ID, and Password are required only if you would like to have controlled access to the software. Otherwise, please leave the fields blank.'
- Section Header: 'Input a Director's name, ID, and password'.
- Form Fields:
  - Name
  - Login ID (pre-filled with 'Admin')
  - New
  - Confirmation
- Buttons: 'Reg.' and 'Cancel'.

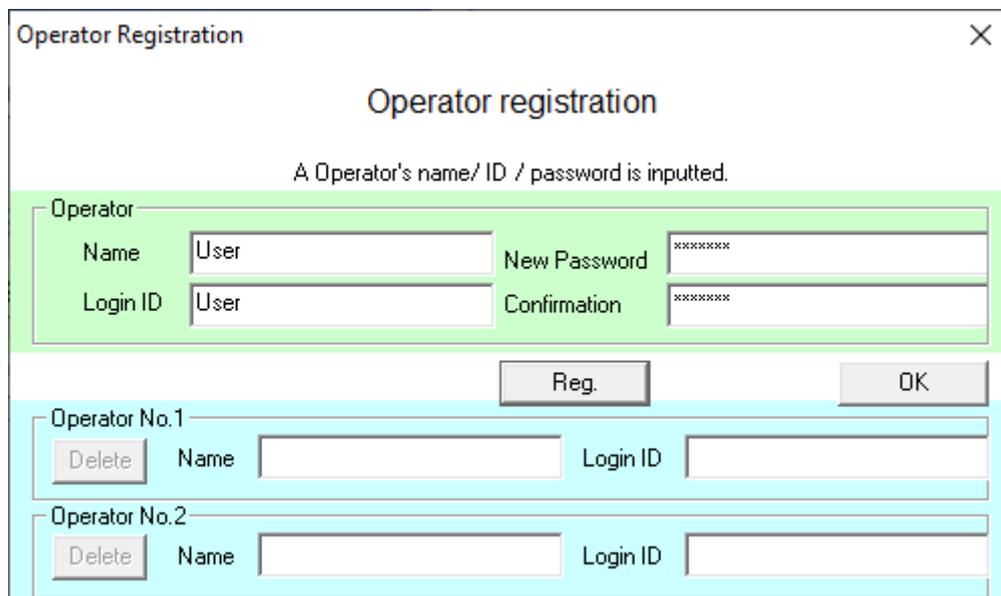

The 'Operator Registration' dialog box contains the following elements:

- Title bar: 'Operator Registration' with a close button (X).
- Section Header: 'Operator registration'.
- Instructional Text: 'A Operator's name/ ID / password is inputted.'
- Form Fields:
  - Operator:
    - Name: User
    - Login ID: User
    - New Password: [masked]
    - Confirmation: [masked]
- Buttons: 'Reg.' and 'OK'.
- Operator List:
  - Operator No.1: [Delete] Name [ ] Login ID [ ]
  - Operator No.2: [Delete] Name [ ] Login ID [ ]

Feb 21 2020