



## iPRECIO IMS-200 Application Software Version 1.3 Revision 532 Renew

- DISK1
- DISK2
- DISK3
- DISK4
- Documents for Returns
- Driver
- iPRECIO Workflow Examples
- Surgical Videos (links to download videos)
- Technical Notes
- User Manual - FAQ
- AUTORUN
- FirstStep
- Important Notice-Total Dead Volume
- Installation Instructions for iPRECIO install CD 13rev532 (Feb 2020)
- Readme (for Serial Number information)


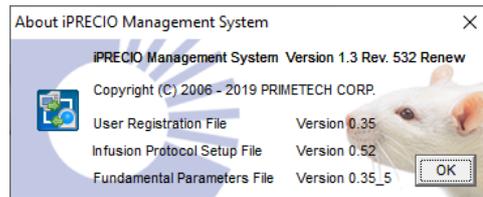


### From Install CD

1. Run <FirstStep> as Administrator (Administrator Rights Requested) for installation. Screen below <iPRECIO Management System Setup> will appear.
2. Plug UCD-200 device to PC – Required drivers will be installed (internet access required)
  - a. Driver Install Option shown below is no longer required as Windows 10 will automatically install the required drivers when UCD-200 device is plugged to PC USB port. **If the right drivers are not automatically installed, install manually using drivers on CD (Folder: Drivers 2020)**
  - b. Drivers are also on CD for manual installation for PCs without internet connection (Folder: Drivers 2020)
  - c. After completed, Run <Install> to start the installation process
3. Then follow instructions on the PC/Screen

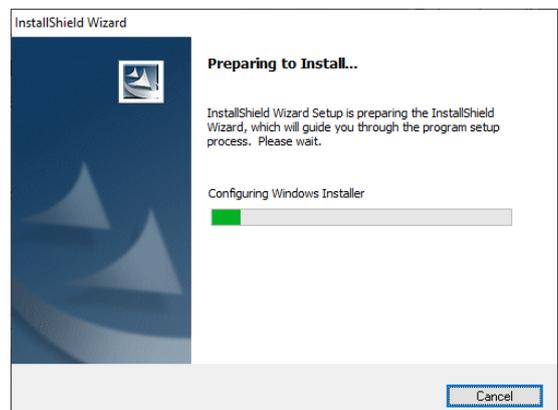


Install Shield Wizard appears, Click <Next> and then accept License Agreement and click <Next>

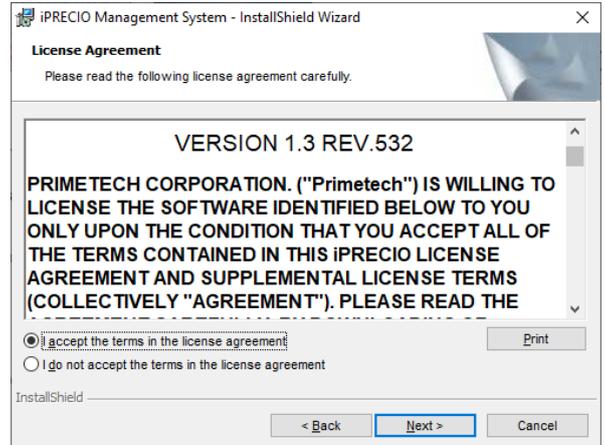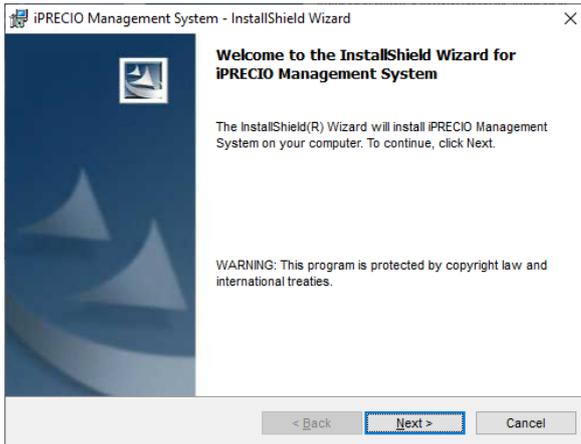


You will need to use the software key from the CD included in the package.

- The iPRECIO Software CD is labeled with serial number (S/N),
- IMSV13R532XXX where XXX is a unique number. Screenshot does not show XXX and only 532
- Refer to iPRECIO User Manual for full installation instructions. (User Manual Version is 1.3rev531)
- See page 14 for S/N details and page 13 for full installation guide. (User Manual Version is 1.3rev531)
- **Latest version is 1.3rev532**



Fill as required and click on <Next>

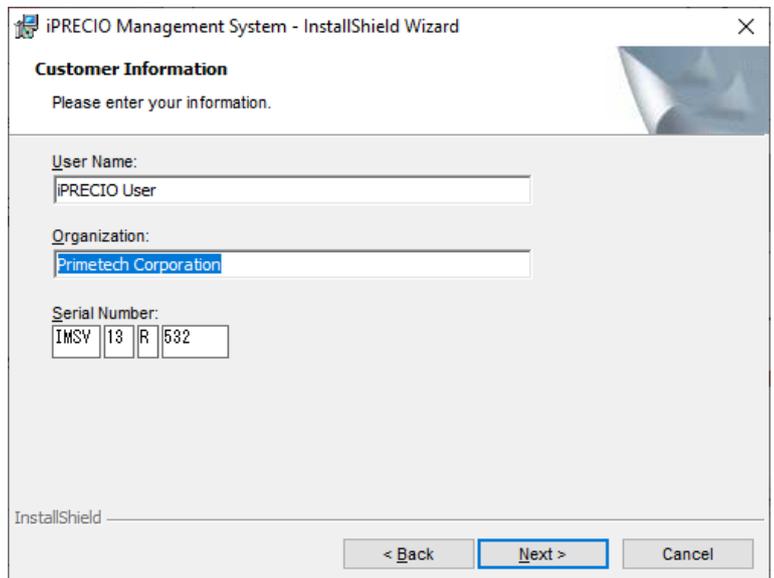

Confirm or change folder location and click on <Next> to proceed.

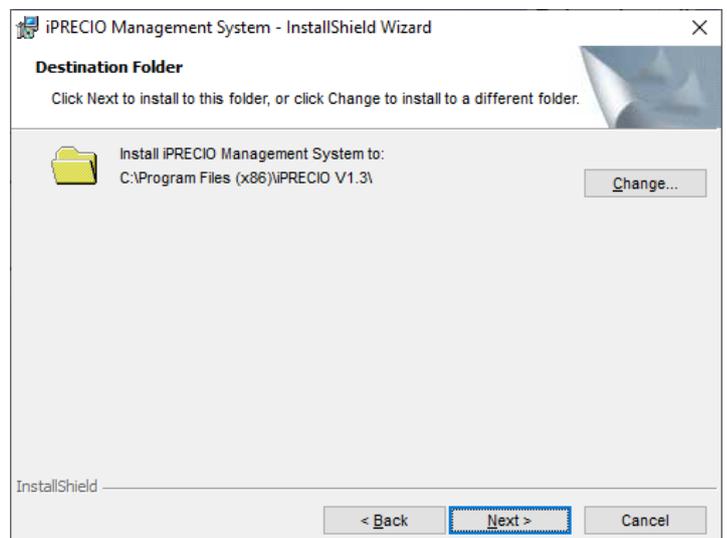


Click on **<Install>**

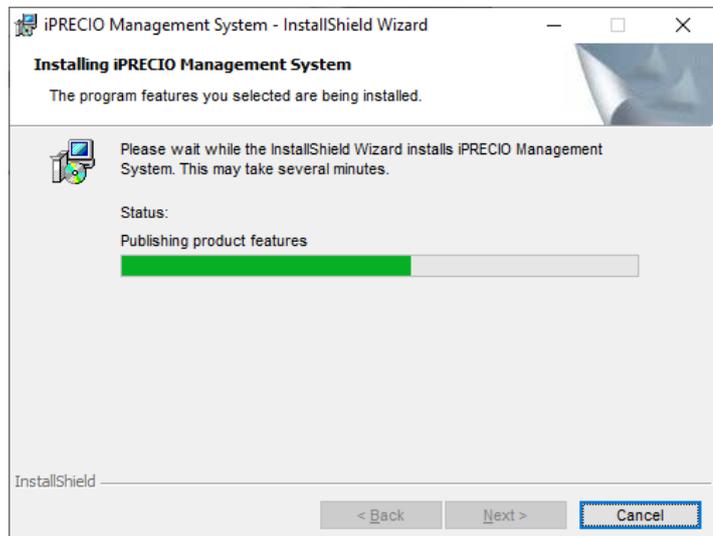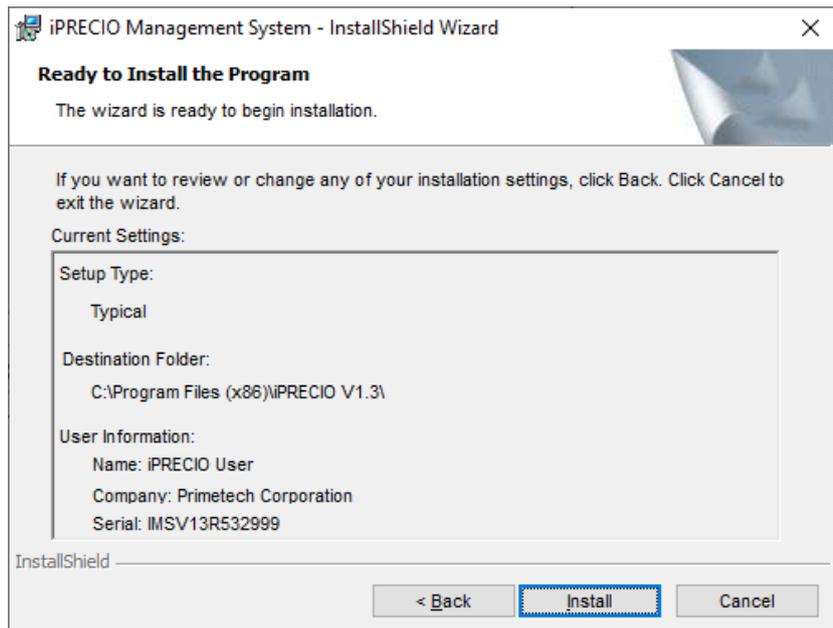


Click on **<Finish>** and close Setup Window as installation is complete.

If required setup shortcut to desktop.

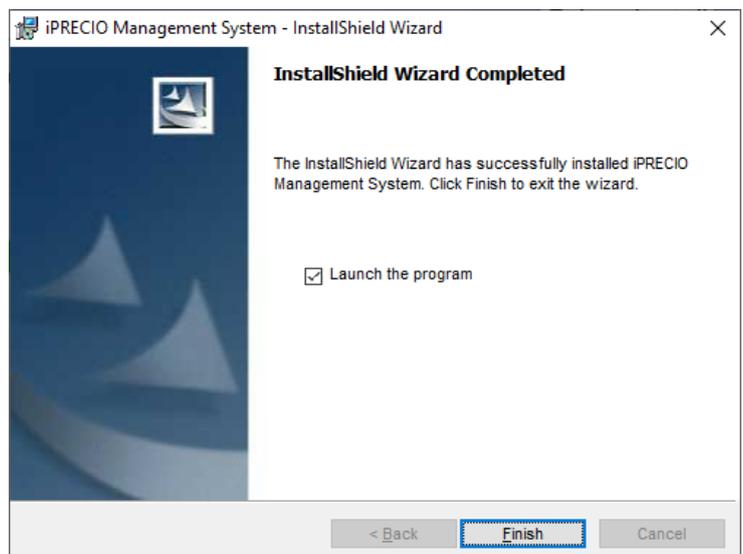


Run Application and See Page 16 of User Manual for software installation.

**Software Setup – Page 16 of User Manual**

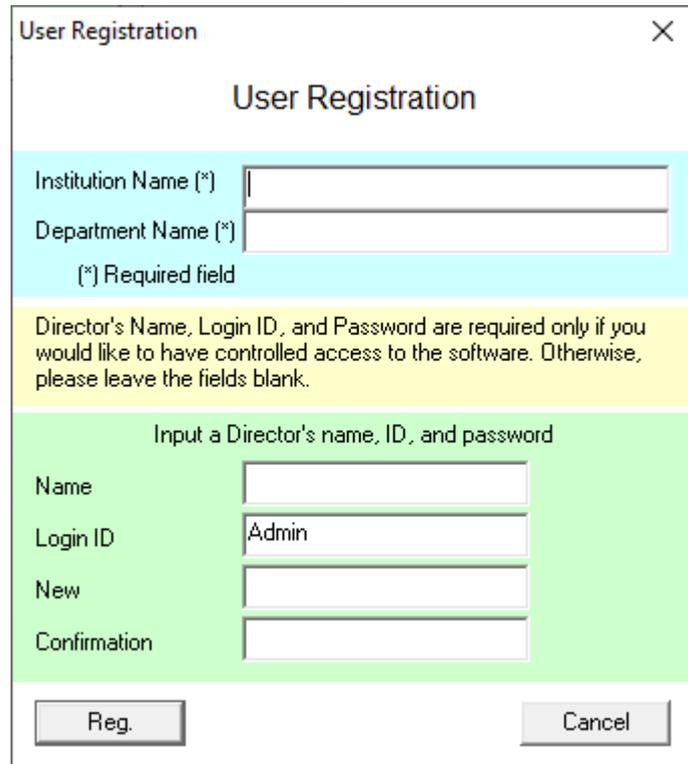



The 'User Registration' dialog box features a title bar with a close button. The main title is 'User Registration'. Below the title, there are two required text input fields: 'Institution Name (\*)' and 'Department Name (\*)'. A note below these fields states '(\*) Required field'. A yellow informational box contains the text: 'Director's Name, Login ID, and Password are required only if you would like to have controlled access to the software. Otherwise, please leave the fields blank.' Below this, a green section titled 'Input a Director's name, ID, and password' contains four input fields: 'Name', 'Login ID' (with 'Admin' pre-filled), 'New', and 'Confirmation'. At the bottom, there are 'Reg.' and 'Cancel' buttons.


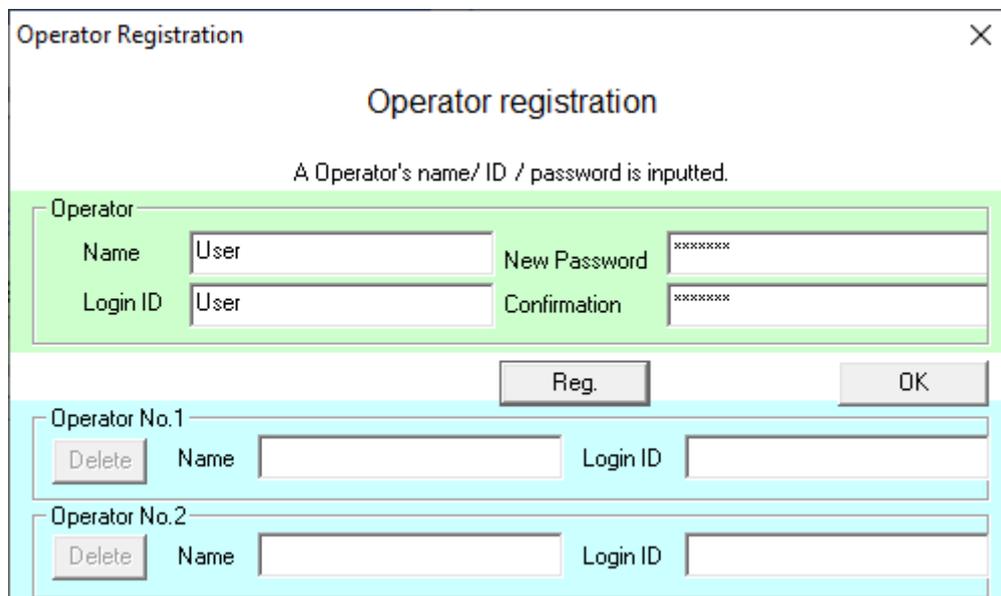



The 'Operator Registration' dialog box has a title bar with a close button. The main title is 'Operator registration'. Below the title, a message reads 'A Operator's name/ ID / password is inputted.' The form is divided into three sections. The top section, titled 'Operator', has a light green background and contains four input fields: 'Name' (with 'User' pre-filled), 'New Password' (with '\*\*\*\*\*' pre-filled), 'Login ID' (with 'User' pre-filled), and 'Confirmation' (with '\*\*\*\*\*' pre-filled). Below this section are 'Reg.' and 'OK' buttons. The bottom section, titled 'Operator No.1', has a light blue background and contains a 'Delete' button, a 'Name' input field, and a 'Login ID' input field. A second identical section for 'Operator No.2' is located below the first.

June 21 2021

**UCD-200 Manual Driver Install via provided drivers in folder driver2020.**  
**Use this procedure if automatic install from the internet does not work correctly /or has errors in Device Manager.**

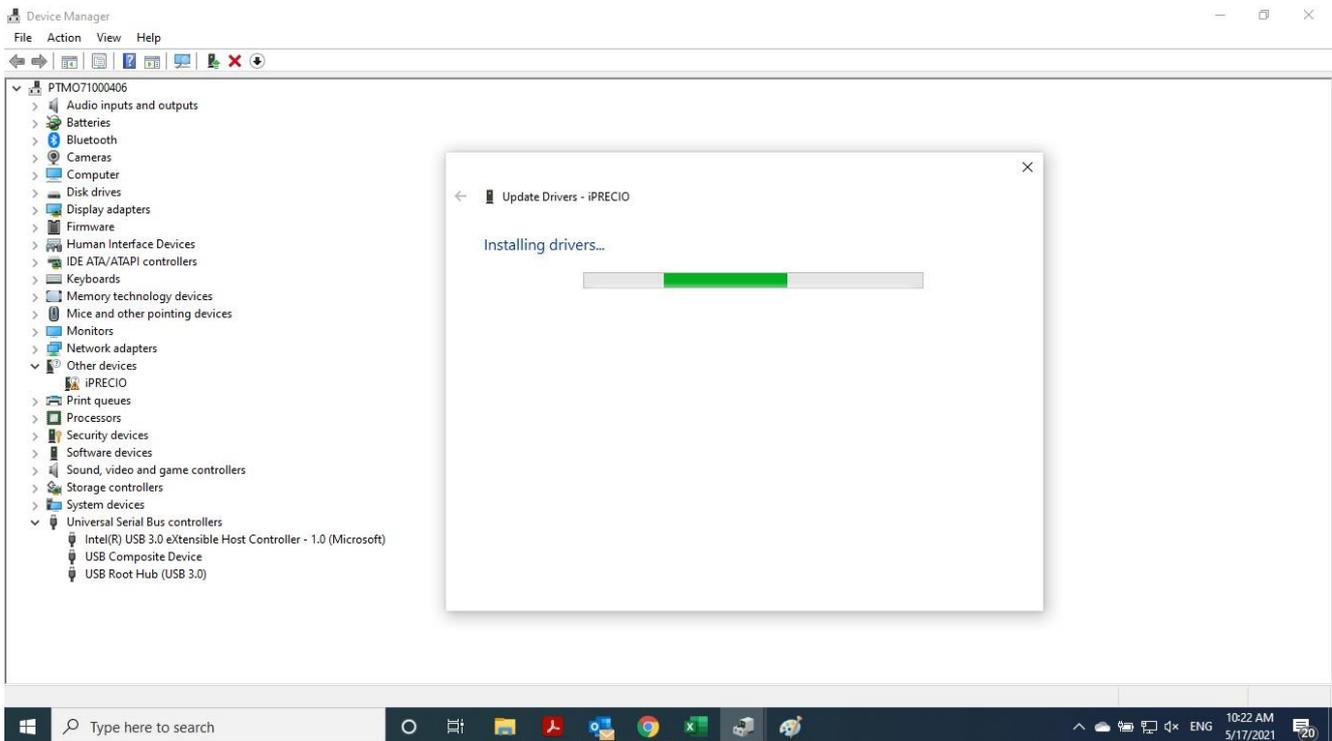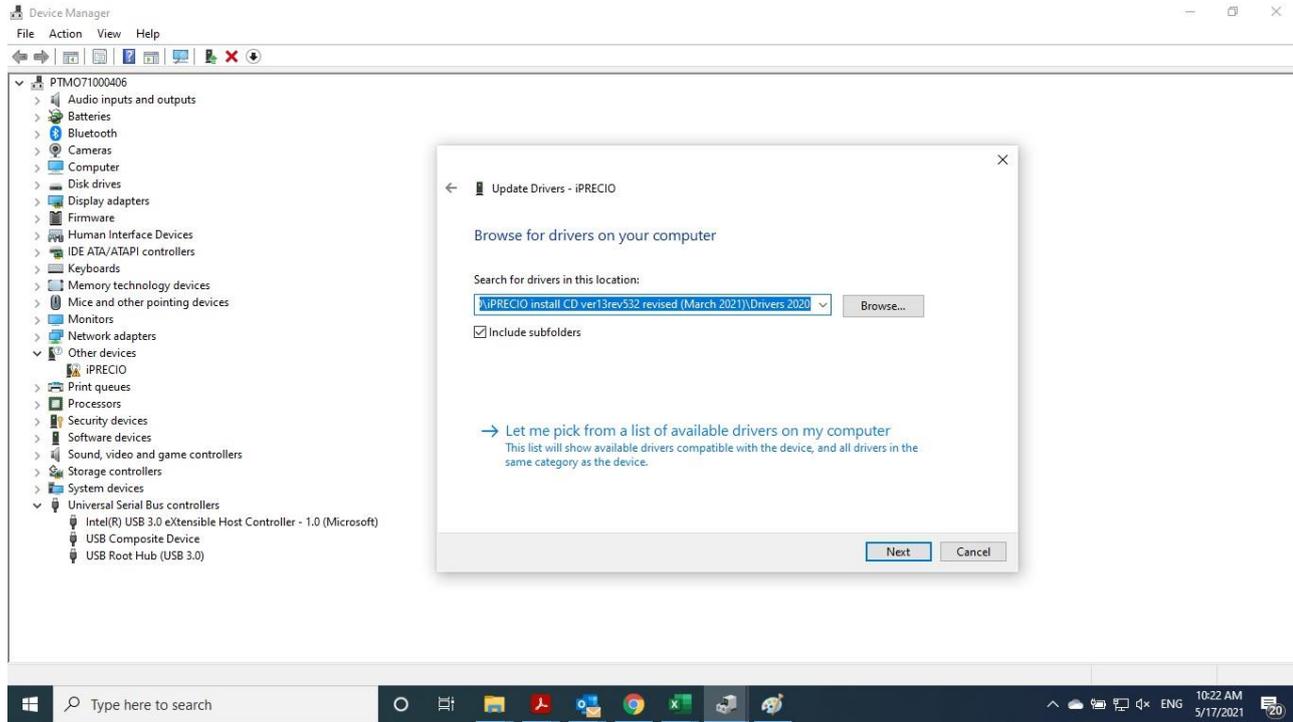

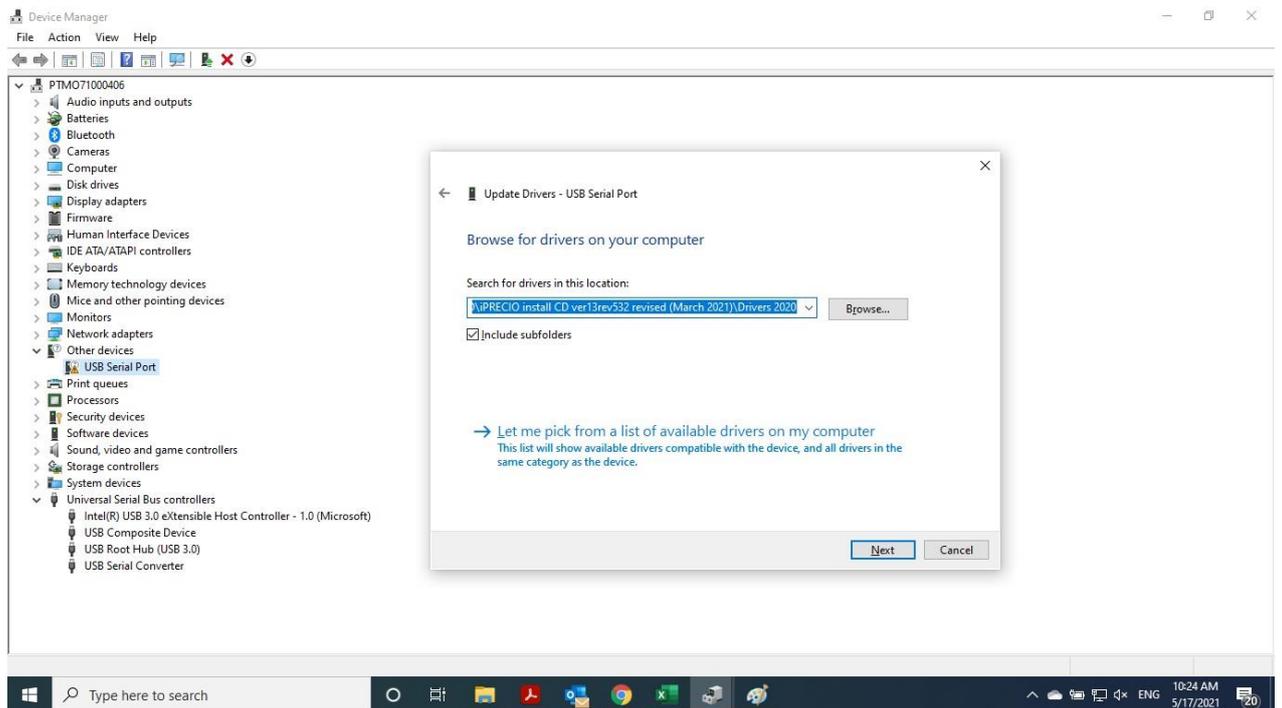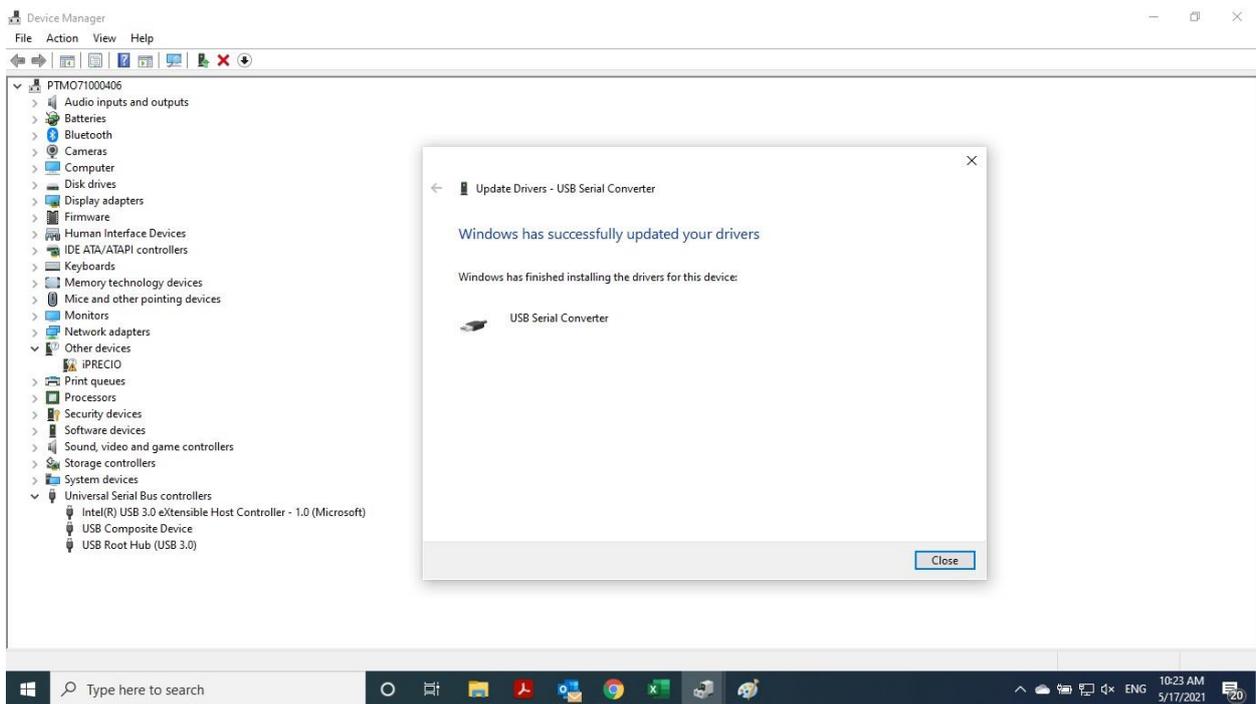

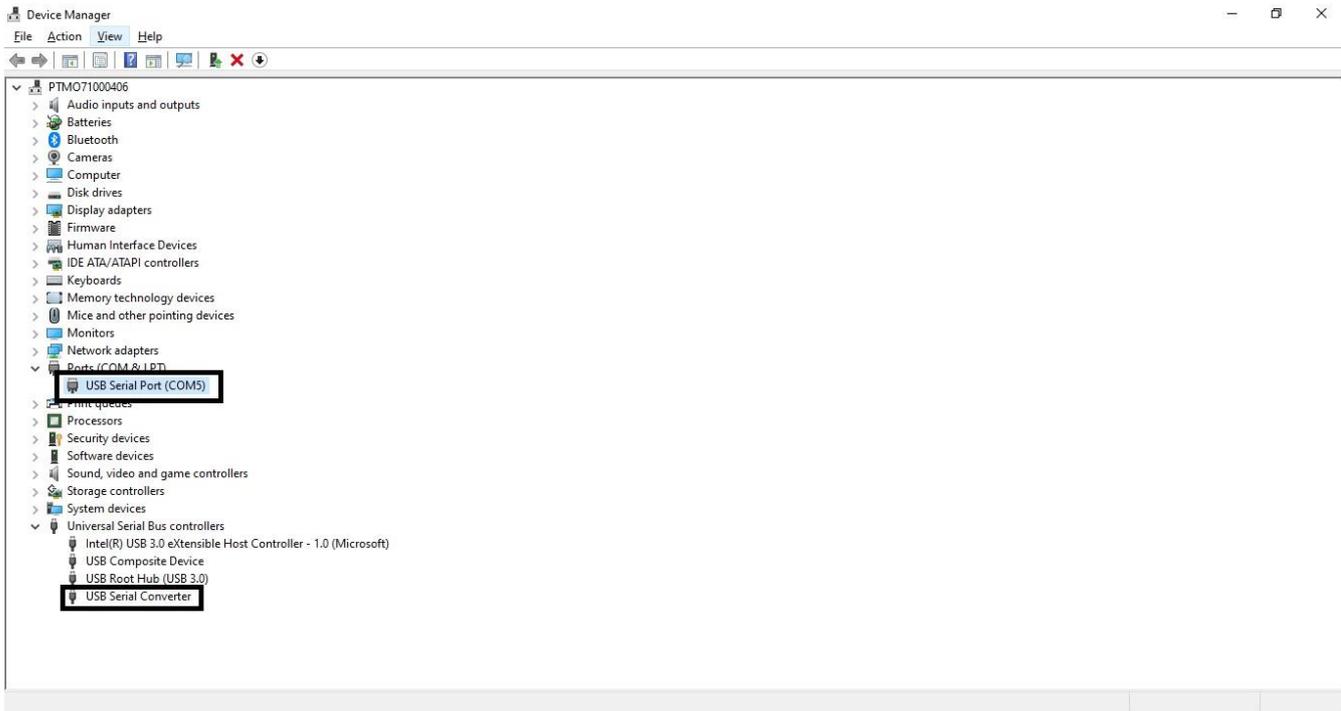

June 2021